

FRANKA RESEARCH 3

Hardware Manual

© Copyright 2025
Franka Robotics GmbH
Frei-Otto-Straße 20
80797 Munich
Germany

This document and extracts of it may only be reproduced or made available to third parties with the explicit authorization of Franka Robotics GmbH.

This Franka Research 3 hardware manual is designed to provide comprehensive guidance for both Franka Research 3 and Franka Production 3.

The content of this document has been carefully checked against compliance with the hardware and software described. However, discrepancies cannot entirely be ruled out, which is why we assume no liability for complete compliance.

In the interest of our customers, we reserve the right to make improvements and corrections to hardware, software, and documentation at any point in time without notice.

We are always grateful for your feedback at documentation@franka.de.

The English documentation is the ORIGINAL DOCUMENTATION. Other languages are translations of the original document.

Complementary documents to this hardware manual are:

- Datasheet Franka Research 3 with FR3 Arm v1 (Document number: R02212)
- Quick Guide for Installation FR3 (Document number: R02040)

Document name: Hardware Manual Franka Research 3

Document number: R02210

Hardware Version: Franka Research 3 with Arm v1

Change Log:

Content Iteration	Release Date	Release notes / Changes
1.0	November 2025	Release of Franka Research 3 Manual with Arm v1
1.1	November 2025	Removed X2 – Digital IOs description and added note in chapter Equipment Overview
1.2	December 2025	Updated information in chapter 4.7 "Installation of safety peripherals" and corrected LED color reference table in chapter 9.5



Get your manual and additional supporting material in English and other languages at www.franka.de/documents.

TABLE OF CONTENTS

1	ABOUT HARDWARE MANUAL.....	6
1.1	Software and hardware version.....	6
1.3	Applicable document.....	6
1.4	Before you begin	6
1.4.1	Intended audience and training requirements.....	6
2	RIGHTS OF USE AND PROPERTY RIGHTS.....	7
2.1	General.....	7
2.2	Identification.....	7
3	DECLARATION OF INCORPORATION AND CERTIFICATES.....	8
3.1	Declaration of Incorporation.....	8
3.2	Certificates.....	10
3.3	Further Statements	10
3.3.1	RoHS / REACH / WEEE / Battery Directive.....	10
3.3.2	China RoHS 2	11
3.4	Labeling on the Equipment.....	12
4	SAFETY	14
4.1	Safety Instructions and General Indications	14
4.2	Notice of Liability.....	15
4.3	Intended Use	15
4.4	Misuse.....	15
4.5	General Possible Dangers and Safety Measures when Working with Robots.....	16
4.6	Application Related Possible Dangers and Safety Measures.....	19
4.7	Installation of Safety Peripherals.....	21
4.8	Fail-safe locking system	22
4.9	Manually Moving the Arm	22
4.10	Safety Functions	25
4.11	Additional information for planning and initial installation of a robot system	28
5	EQUIPMENT OVERVIEW.....	32
5.1	The Arm	33
5.2	Control	39
6	SCOPE OF DELIVERY AND ADDITIONAL EQUIPMENT	40
6.1	Included in the box.....	40
6.2	Not included in the box.....	41
6.3	Available Spare Parts and Accessories	42
7	MOUNTING AND INSTALLATION	43
7.1	Unpacking the Equipment.....	44
7.2	Correct Installation Site.....	48
7.2.1	Maximum-and safeguarded space	48
7.2.2	Ambient conditions: Arm	49

7.2.3	Ambient conditions: Control	50
7.3	Preparing the Installation Site	51
7.3.1	Arm	52
7.3.2	Control	53
7.4	Mounting the Arm	54
7.5	Positioning the Control	55
7.6	Wiring and Electrical Installation	56
7.6.1	Connection diagram	57
7.6.2	Interfaces	57
7.6.3	Connecting the functional earth	60
7.6.4	Wiring	61
7.7	Mounting End Effectors	65
7.8	Practical Tips for Usage and Positioning of Franka Research 3	67
7.8.1	Energy consumption	67
7.8.2	ESD limits	67
7.8.3	Design of the workspace	68
7.8.4	Personal safety and ergonomics	69
8	USAGE	71
8.1	Switching On	71
8.2	Safety Relevant Tests of Franka Research 3	73
8.2.1	Self-test of the robotic system	73
8.2.2	Regular testing of safety functions	73
8.2.3	Testing the emergency stop	74
8.3	Switching Off and Restarting	75
9	ROBOT LED INDICATOR SYSTEM	77
9.1	Overview of the status indicators	77
9.2	LED Activation Behavior	77
9.3	Flashing Patterns	78
9.4	LED Priority Logic	78
9.5	LED Color Reference Table	79
10	MAINTENANCE AND DISPOSAL	81
10.1	Maintenance	81
10.2	Cleaning	81
10.3	Disposal	82
10.4	Mechanical exchange of Control	82
11	TROUBLESHOOTING	83
12	TECHNICAL DATA	83
12.1	Ambient conditions for delivery and transport	84
13	TRANSPORTATION AND HANDLING	84
13.1	Procedure	84
13.1.1	Step 1. Transportation pose of the Arm	85

13.1.2	Step 2. Handling and lifting	85
13.1.3	Step 3. Re-packing the Arm.....	87
14	APPENDIX	89
14.1	Stopping Times and Distances.....	89
14.2	Stop Category 1	90
14.2.1	Joint 1	90
14.2.2	Joint 2	92
14.2.3	Joint 3	94
14.2.4	Joint 4	96
14.3	Stop Category 2	98
14.3.1	Joint 1	98
14.3.2	Joint 2	100
14.3.3	Joint 3	102
14.3.4	Joint 4	104
14.4	Response times	106
14.5	Safe Position Accuracy	106
15	GLOSSARY	107
16	INDEX	112
17	TABLE OF FIGURES.....	114

1 ABOUT HARDWARE MANUAL

This hardware manual provides essential information to ensure the safe and correct use of the Franka robot. It includes detailed guidance on identifying the robot's components, performing general maintenance, understanding built-in safety functions, and following step-by-step instructions for initial setup and usage.

All personnel must read and fully understand this manual before beginning any work with the robot. Safe operation requires strict adherence to all safety guidelines and handling instructions provided in this manual.

1.1 Software and hardware version

This hardware manual applies to Franka Research 3 with Arm version v1
This version is compatible with system image version 5.4.0 to 5.8.

1.3 Applicable document

In addition to this manual, the following document also applies:

- **Document Number: R02216**
 - **Operating Manual:** Franka Research 3 with system image version 5.6.0
 - **Operating Manual:** Franka Research 3 with system image version 5.8
- **Product Manual Franka Research 3 (Document number: R02210)**
 - System Image v5.8.0
 - System Image v5.7.0
 - System Image v5.6.0
 - System Image v5.5.0
 - System Image v5.4.0

NOTICE

This Franka Research 3 with Arm v1 hardware manual configuration is derived from the Product Manual Franka Research 3 (Document number: R02210) and is compatible with System Image versions 5.4.0 through 5.8.0.

1.4 Before you begin

1.4.1 Intended audience and training requirements

This manual is intended for qualified technical personnel responsible for installation, operation, and maintenance of the Franka Research 3 system.

Users must:

- Be trained in handling industrial robots and familiar with applicable safety regulations (e.g. EN ISO 10218-2).
- Understand basic mechanical and electrical safety principles.
- Be authorized by their employer to perform the described tasks.

Untrained or unauthorized people are **not permitted** to install, operate, or service this product.

2 RIGHTS OF USE AND PROPERTY RIGHTS

2.1 General

Protected trademarks

This hardware manual refers to protected trademarks not explicitly designated in the continuing text. The absence of such indication may not imply that the corresponding product name is free of third-party rights. The following trademarks are protected trademarks:

Franka and Franka Robotics are registered trademarks.

Microsoft is a registered trademark, and Windows is an indication of the Microsoft Corporation in the United States and other countries.

GOOGLE, Mozilla, Firefox, CHROME, ITEM are registered trademarks.

Trademark rights

The responsible person is not granted any rights or claims to the trademark, logo, or trade names of Franka Robotics.

Use of open-source code

A complete list of all open-source licenses used by Franka Research 3 can be accessed via the Settings menu of the Franka UI.

2.2 Identification

Removal of identification

Copyright notices, serial numbers, and any other labeling that identifies the product or operating software may not be removed or modified.

3 DECLARATION OF INCORPORATION AND CERTIFICATES

3.1 Declaration of Incorporation

	Declaration of Incorporation according to directive 2006/42/EC on machinery (Annex II B) for partly completed machinery	
--	--	--

Description of the partly completed machinery:

Product identification: Franka Production 3 / Research 3 components: *Control, Arm*

Model/Type:

Control (#295341) in combination with *Arm FP3* (#290101) or *FR3* (#290102), each with running number starting from 2320001.

We declare that the product complies with the following essential safety and health requirements set out in Annex I of the Machinery Directive 2006/42/EC:

1.1.2; 1.1.3; 1.1.5; 1.1.6; 1.2.1; 1.2.2; 1.2.3; 1.2.4.1; 1.2.4.2; 1.2.4.3; 1.2.4.4; 1.2.5; 1.2.6; 1.3.1; 1.3.2; 1.3.3; 1.3.4; 1.3.6; 1.3.7; 1.3.8; 1.3.8.1; 1.3.8.2; 1.3.9; 1.4.1; 1.4.2.1; 1.4.2.2; 1.4.3; 1.5.1; 1.5.2; 1.5.3; 1.5.4; 1.5.5; 1.5.6; 1.5.8; 1.5.9; 1.5.10; 1.5.11; 1.5.13; 1.5.14; 1.6.1; 1.6.3; 1.6.4; 1.7.1; 1.7.1.1; 1.7.1.2; 1.7.2; 1.7.3; 1.7.4; 1.7.4.1; 1.7.4.2; 2; 2.2.1; 2.2.1.1; 4; 4.1.2.3; 4.2.1; 4.3.3; 4.4.2

In addition, the partly complete machinery is in conformity with the following EU Directives:

Directive 2014/35/EU relating to electrical equipment (LVD)

Directive 2014/30/EU relating to electromagnetic compatibility (EMC)

Directive 2011/65/EU relating to hazardous substances in electrical and electronic equipment (RoHS)

Directive 94/62/EC relating to packaging and packaging waste

We declare that the relevant technical documentation is compiled in accordance with part B of Annex VII.

Applied harmonized standards:

Electrical safety

Standard	Name
EN 60204-1:2018	Safety of machinery – Electrical equipment of machines
IEC 60204-1:2016	– Part 1: General requirements
EN 60664-1:2007	Insulation coordination for equipment within low-voltage systems –
IEC 60664-1:2007	Part 1: Principles, requirements and tests
EN 60664-4:2006	Insulation coordination for equipment within low-voltage systems –
IEC 60664-4:2005	Part 4: Consideration of high-frequency voltage stress
EN 60529:1991/A1:2000/A2:2013	Degrees of protection provided by enclosures (IP Code)
IEC 60529:1989/AMD1:1999/A2:2013	
EN 61010-1:2010/A1:2019/AC:2019-04	Safety requirements for electrical equipment for measurement,
IEC 61010-1:2010/AMD1:2016/COR1:2019	control, and laboratory use – Part 1: General requirements
EN IEC 61010-2-201:2018	Safety requirements for electrical equipment for measurement,
IEC 61010-2-201:2017	control, and laboratory use – Part 2-201: Particular requirements for control equipment
EN 61800-5-1:2007/A1:2017-04	Adjustable Speed Electrical Power Drive Systems
IEC 61800-5-1:2007	– Part 5-1: Safety requirements – Electrical, thermal and energy
/AMD1:2016	

Machinery safety

Standard	Name
EN ISO 10218-1:2011	Robots and robotic devices – Safety requirements for industrial robots
ISO 10218-1:2011	– Part 1: Robots
EN ISO 12100:2010	Safety of machinery – General principles for design
ISO 12100:2010	– Risk assessment and risk reduction

DECLARATION OF INCORPORATION AND CERTIFICATES

EN ISO 13849-1:2015	Safety of machinery – Safety-related parts of control systems
ISO 13849-1:2015	– Part 1: General principles for design
EN ISO 13849-2:2012	Safety of machinery – Safety-related parts of control systems
ISO 13849-2:2012	– Part 2: Validation
EN ISO 13850:2015	Safety of machinery – Emergency stop function
ISO 13850:2015	– Principles for design
EN ISO 14118:2018	Safety of machinery
ISO 14118:2017	– Prevention of unexpected start-up
EN 61310-1:2008	Safety of machinery – Indication, marking and actuation
IEC 61310-1:2007	– Part 1: Requirements for visual, acoustic and tactile signals
EN 61310-2:2008	Safety of machinery – Indication, marking and actuation
IEC 61310-2:2007	– Part 2: Requirements for marking

EMC

Standard	Name
EN IEC 61000-6-1:2019	Electrical compatibility (EMC) – Part 6-1: Generic standards
IEC 61000-6-1:2016	– Immunity standard for residential, commercial and light-industrial environments
EN IEC 61000-6-2:2019	Electrical compatibility (EMC) – Part 6-2: Generic standards
IEC 61000-6-2:2016	– Immunity standard for industrial environments
EN 61000-6-3:2007	Electrical compatibility (EMC) – Part 6-3: Generic standards
/A1:2011/AC:2012-08	– Emission standard for residential, commercial and light-industrial environments
IEC 61000-6-3:2020	
EN IEC 61000-6-4:2019	Electrical compatibility (EMC) – Part 6-4: Generic standards
IEC 61000-6-4:2018	– Emission standard for industrial environments
EN 61000-6-7:2015	Electrical compatibility (EMC) – Part 6-7: Generic standards
IEC 61000-6-7:2014	– Immunity requirements for equipment intended to perform functions in a safety-related system (functional safety) in industrial locations
EN 61326-3-1:2017	Electrical equipment for measurement, control and laboratory use – EMC
IEC 61326-3-1:2017	requirements – Part 3-1: Immunity requirements for safety-related systems and for equipment intended to perform safety-related functions (functional safety)
	– General industrial applications
CISPR 11:2015+AMD1:2016	Industrial, scientific and medical equipment
+AMD2:2019 CSV	– Radio-frequency disturbance characteristics – Limits and methods of measurement

We commit to transmit, in response to a reasoned request by the market surveillance authorities, relevant documents on the partly completed machinery. The industrial property rights remain unaffected!

Important note!

The partly completed machinery must not be put into service until the final machinery into which it is to be incorporated has been declared in conformity with the provisions of Directive 2006/42/EC on Machinery, where appropriate, and until the CE Declaration of Conformity according to Annex II A is issued.

Representative in EU, authorized to compile the relevant technical documentation:

Franka Robotics GmbH
Frei-Otto-Straße 20
80797 München
Germany

Manufacturer:

Franka Robotics GmbH
Frei-Otto-Straße 20
80797 München
Germany

Date, location

25.01.2024
Munich, Germany

Jin Zhengxun, CEO



Document No.: R02011

Page 2/2

3.2 Certificates

Certificates by TÜV SÜD RAIL and TÜV SÜD PS can be found under the following link:

www.franka.de/documents

3.3 Further Statements

3.3.1 RoHS / REACH / WEEE / Battery Directive

<p>Further Information</p> <p>Status: 01.04.2022 Franka Research 3</p>
<p>Restriction of Hazardous Substances (RoHS):</p> <p>The components <i>Control and Arm</i> do not fall within the scope of EU RoHS Directive 2011/65/EU but still meet the requirements of the restricted substances and maximum concentration values that are allowed in homogenous materials:</p> <ul style="list-style-type: none"> • Lead (0.1 %) • Mercury (0.1 %) • Cadmium (0.01 %) • Hexavalent chromium (0.1 %) • Polybrominated biphenyls (PBB) (0.1 %) • Polybrominated diphenyl ethers (PBDE) (0.1 %) <p>The following exceptions are also applied:</p> <p>6a: Lead as an alloying element in steel for machining purposes and in galvanized steel containing up to 0.35 % lead by weight</p> <p>6b: Lead as an alloying element in aluminum containing up to 0.4 % lead by weight</p> <p>6c: Copper alloy containing up to 4 % lead by weight</p> <p>7a: Lead in high melting temperature type solders (i.e., lead-based alloys containing 85 % by weight or more lead)</p> <p>7c-l: Electrical and electronic components containing lead in a glass or ceramic other than dielectric ceramic in capacitors, e.g., piezo-electronic devices, or in a glass or ceramic matrix compound</p>
<p>REACH:</p> <p>FRANKA ROBOTICS GmbH is a "downstream user" as defined in REACH. Our products are exclusively non-chemical products (manufactured items). In addition, no substances are released under normal conditions of use and conditions that can reasonably be predicted (Article 7, REACH).</p> <p>We confirm that our products do not contain more than 0.1 percent by mass of any of the listed substances on the published ECHA candidate list (SVHC) unless covered by RoHS exceptions (see above). Extensions published by the ECHA candidate list match our products, and if it is known that one of these newly added substances is contained in our products, we will inform you immediately.</p> <p>This confirmation was created based on currently available information from our suppliers.</p>
<p>WEEE Directive:</p> <p>The components <i>Control and Arm</i> are not subject to the WEEE Directive 2002/96/EC for collection, recycling, and recovery for electrical goods.</p>

DECLARATION OF INCORPORATION AND CERTIFICATES

Battery Directive:

The product Control contains a BIOS battery.

Disposal of batteries:

There is an obligation to return rechargeable and non-rechargeable batteries by Battery Directive 2006/66/EC; do not dispose of them with consumer waste. Dispose of them according to statutory orders and bring them to a recycler. Batteries will be recycled.

The signs below the crossed-out trashcan indicate the substances lead (Pb), cadmium (Cd), or mercury (Hg).



3.3.2 China RoHS 2



Part Name 部件名称	Toxic and Hazardous Substances and Elements / 有毒和有害的物质和元素					
	Lead 铅 (Pb)	Mercury 汞 (Hg)	Cadmium 镉 (Cd)	Hexavalent Chromium 六价铬 (Cr (VI))	Polybrominate d biphenyls 多溴联苯 (PBB)	Polybrominated diphenyl ethers 多溴二苯醚 (PBDE)
Control 控制器	X	O	O	O	O	O
Arm FP3 臂架 FP3	X	O	O	O	O	O
Robot Connecting Cable 连接线 2.5m / 5m /10m	O	O	O	O	O	O
Emergency Stop Device 紧急停止装置	O	O	O	O	O	O
External Enabling Device 外部支持设备	O	O	O	O	O	O
<p>This Table is prepared in accordance with the provisions of SJ/T 11364-2006. 本表根据SJ/T 11364-2006的规定编制。</p> <p>O: Indicates that said hazardous substance contained in all of the homogeneous materials for this part is below the limit requirement of GB/T 26572-2011. O: 表示该部分所有均质材料中的所述有害物质低于GB/T 26572-2011的限量要求。</p> <p>X: indicates that said hazardous substance contained in at least one of the homogenous materials used for this part is above the limit requirement of GB/T 26572-2011 X: 表示本试验用同质材料中至少有一种所含的有害物质高于GB/T 26572-2011的限量要求。</p> <p>(Enterprises may further provide in this box technical explanation for marking "X" based on their actual circumstances) (企业可根据自己的实际情况，在此框中进一步提供标示 "X" 的技术解释)</p>						

3.4 Labeling on the Equipment

Arm

Type label



Fig. 3.1: Type label

Emergency unlock label

Three Emergency Unlock Labels on the Arm point to the insertion points where the Emergency Unlocking Tool needs to be inserted to manually unlock the fail-safe locking system in case of an emergency.



Fig. 3.2: Emergency unlock label

WARNING

Falling heavy Arm when using Emergency Unlocking Tool

Risk of being trapped by the Arm when unlocking joints

- Do support the Arm before and during unlocking it.
- Do not place your head or other body parts between or underneath links of the Arm.
- Do not place body parts (especially hands, fingers) between links of the Arm, the end effector, or stationary objects.
- Do not use the Emergency Unlocking Tool while the Arm is powered.
- Only use the supplied Emergency Unlocking Tool.
- Store Emergency Unlocking Tool near the Arm.

DECLARATION OF INCORPORATION AND CERTIFICATES

Hot Surfaces label



Fig. 3.3: Hot Surface label

WARNING

Hot Surfaces and Guiding

At ambient temperatures above 30 °C the robot surface can become too hot to touch. Therefore, using the Assist feature in Execution mode is not permitted above 30 °C.

Functional earth label

The functional earth label points to the location where functional earth can be connected to the Arm.



Fig. 3.4: Functional earth label

Lifting position label

The lifting position label indicates the points at which the Arm may be lifted.



Fig. 3.5: Lifting position label

Control

Type label

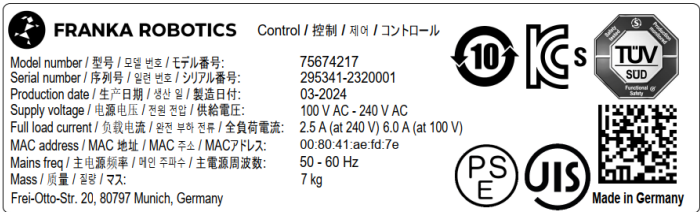


Fig. 3.6: Type label control

SAFETY

External Enabling Device

Type label



Fig. 3.7: Type label external device

Emergency Stop Device

Type label



Fig. 3.8: Type label emergency stop device

4 SAFETY

4.1 Safety Instructions and General Indications

Warning notices

Before installing, starting up, and operating the device, carefully read this manual and any additional documentation. Take NOTICE of the safety instructions as well as general indications.

Warning notices are posted as follows:

⚠ CAUTION

Type of hazard
Potential consequences of hazard
Evasive/avoidance actions to be taken

The following warning notices are used in this manual:

⚠ DANGER

DANGER indicates a hazardous situation that, if not avoided, will result in death or serious injury.

⚠ WARNING

WARNING indicates a hazardous situation that, if not avoided, could result in death or serious injury.

SAFETY

CAUTION

CAUTION indicates a hazardous situation that, if not avoided, could result in minor or moderate injury.

NOTICE

NOTICE indicates information considered important but not hazard related.

SAFETY- INSTRUCTION

SAFETY INSTRUCTION indicates processes that need to be strictly observed.

Indications



Indicates where further information can be obtained.

4.2 Notice of Liability

Franka Research 3 has been developed according to the relevant quality standards. A hazard and risk assessment according to EN ISO 12100 has been carried out through the course of development and is the basis for Franka Research 3 and this manual.

The present document includes assembly instructions for Franka Research 3 as a partly completed machinery. It contains descriptions of the conditions that must be met to correct incorporation in the final machinery not to compromise safety and health (e.g., Annex I of Machinery Directive 2006/42/EC).

4.3 Intended Use

Franka Research 3 is exclusively intended for use in research and development in academic and industrial environments. The system may only be used in sound technical conditions, for its intended purpose, and within the technical specifications and operating conditions, with awareness of safety and possible dangers.

The present Franka Research 3 system is intended exclusively for use as described in this manual.

For the normal and extended working conditions under which the robot should operate, see chapter Error! Reference source not found. "Technical Data" and Chapter 7.2 "Correct Installation Site".

4.4 Misuse

DANGER

Misuse of Franka Research 3

Danger to life and risk of injury, as well as risk of impairments, damage to the robot, and other material assets.

- Only use Franka Research 3 in sound technical condition.
- Only use Franka Research 3 in the ambience and operating conditions described in this document.

SAFETY

The misuse of Franka Research 3 voids the manufacturer's warranty and liability. Any usage different from the intended use is considered misuse and is not permissible.

Misuse is any use that deviates from the warnings, notices, and instructions in this manual and the Getting Started guide but not limited to, the following uses:

- Transport of people or animals
- Transport without packing pose and original packaging
- Use as climbing assistance
- Leaning against the Arm
- Use in potentially explosive areas
- Use below ground
- Use for the handling of radioactive objects
- Use outdoors
- Use as a medical product
- Use as a service Arm, e.g., for the care of the elderly
- Use in the vicinity of children
- Handling of liquids
- Use in any mounting position other than upright
- Use outside of the specified operating limits

Modifications to Franka Research 3 that Franka Robotics does not explicitly permit are not allowed and will lead to loss of warranty and liability claims. Not permitted modifications include but are not limited to the following:

- Any adaptation of the mechanical structure
- Varnishing
- Wrapping the robotic structure unless using FE certified equipment

Franka Robotics grants only the following modifications to Franka Research 3:

- Installation of an external cable guiding system (can impact the motion and control behavior of the system)
- Mounting equipment to the flange
- Covering screw holes

It is prohibited to open the Arm and the Control enclosure, and other equipment.

The robot may only be used in places where sufficient space and safe use can be ensured.

Franka Robotics is not liable for damage caused by mounted equipment or damage caused by misuse.

4.5 General Possible Dangers and Safety Measures when Working with Robots

Abstract of possible dangers

An extensive but not definitive list of dangers that generally may be presented by a robot system can be found under EN ISO 10218-1:2011 ANNEX A.

Special attention is drawn here to the following dangers that Franka Research 3 may present:

 **DANGER**

Electrical or fire hazards and dangerous fumes

SAFETY

Fire and fumes may cause breathing difficulties, eye irritation, lung damage, poisoning, and may lead to death.

- Do not use Franka Research 3 outside of its indicated specifications.

DANGER

Damaged wires or inadequate electrical installation

Risk of personal injury by electric shock, as well as material damage.

- Only use Franka Research 3 in sound technical condition.
- Only install the emergency stop system with qualified personnel.
- Check cables and electrical installations.

DANGER

Risk of smoldering fires

Too many devices connected to the power supply system may lead to an overload of the electrical installation and may result in smoldering fires, which may cause death or severe damage to persons.

- Connect the Franka Research 3 accordingly to avoid an overload of the electrical installation.
- Install the overload protection devices accordingly.

WARNING

Objects falling from end effectors due to a power supply cut off

Objects falling from the gripper can lead to injuries to hands, fingers, feet, and toes.

- Always wear personal protective equipment (e.g., safety shoes).
- Use the appropriate type of grippers to prevent objects from falling.

WARNING

Falling and unexpected motions of the robot, especially in earthquake-prone areas

Risk of severe injuries such as crushing, tearing of the skin, and puncturing.

- Level the platform.
- Only install the Arm on even, unmoving, and stable platforms. Accelerations and vibrations are not permissible.
- Do not install the Arm on hanging, tilted, or uneven platforms.
- Level the platform in an upright position.
- Ensure a correctly laid out and tight screw connection.
- Tighten the screws after 100 hours of operation with the correct tightening torque.
- If operated in an earthquake-prone area, take the respective hazard and risk assessments into consideration.

SAFETY

WARNING

Dangerous and uncontrolled motions of the Arm

Risk of severe injury, such as crushing, tearing of the skin, and puncturing from the Arm and end effectors.

- Ensure that the end effector and/or the gripped object mass and the Center of Mass (CoM) are parameterized correctly.
- Keep out of the workspace during operation.

CAUTION

Damaged cables, plugs, mechanical housing, or oil leaks

Contact with leaking oil may cause eye or skin irritation.

Electrical risks may lead to severe injury.

- Only use Franka Research 3 in sound technical condition.
- Do not use damaged cables, plugs, and mechanical housing for operation. In cases of doubt, contact Franka Robotics.

Sound condition

CAUTION

Leaking grease or oil through voids of structural elements of the robot

Irritation of skin and eyes.

- Stop operating the machine.
- Contact manufacturer.
- Use gloves.
- Seek medical attention after contact with eyes or skin.

Disruptions

NOTICE

The operator will be informed about possible malfunctions through Desk. Malfunctions need to be rectified before continuing the operation.

- To rectify the possible malfunctions, follow the instructions in the Franka UI. A reboot of the system may be required.

Overload of joints

NOTICE

An offline overload in all the robot's joints may lead to material damage to the robot.

- Perform a risk assessment with foreseeable misuse in mind.
- Adhere to the system's warning to take the robot out of operation if necessary.

4.6 Application Related Possible Dangers and Safety Measures

The following possibly safety-related aspects are to be considered when planning and designing the application and performing the hazard & risk assessment for completed machinery. The integrator is obligated to perform a risk analysis.

Functionalities and features of various safety levels

Franka Research 3 offers functionalities and features of various safety levels. All safety functions and their corresponding safety ratings are described in **section Safety Functions in chapter 4.10 "Safety Functions"**. All other functionalities described throughout this chapter are not classified as safety-rated according to EN ISO 13849-1. You may, therefore, not rely on the availability of these functionalities.

Please note that the integrator is obligated to perform a risk analysis.

Unexpected motion

CAUTION

Unexpected movement of Arm

The use of various applications, handled end effectors and surrounding objects, may cause crushing between arm segments and impact and collision.

- Ensure that the end effector and/or the object mass and the Center of Mass (CoM) are parameterized correctly.
- Keep out of the maximum workspace during operation.

Trapping within a completed machinery

WARNING

Risk of trapping body parts or person

Risk of severe injury, such as crushing, tearing of the skin, and puncturing from the Arm and end effectors.

- Do not place body parts between Arm segments.
- In case of acute mortal danger:
 1. Press the Emergency Stop Device to stop the robot from operating.
 2. Pull or push the Arm out of the dangerous position manually.

Refer to the chapter 4.9 "Manually Moving the Arm" for more information.

Wiring of end effectors

NOTICE

If an end effector is used with Franka Research 3, the maximum attachable payload is reduced by the weight of the end effector and its external wiring.

NOTICE

External wiring adds additional loads and torques to the Arm which may influence Franka Research 3's control performance.

SAFETY

Initiation of motion of externally wired end effectors, associated equipment

NOTICE

Because of different configurations, installed apps and services, Franka Research 3 is able to send protocols to possibly connect machines (incl. starting motion), externally wired end effectors, and other associated equipment. Please be aware of possible related risks of using external equipment.

Single Point of Control

Franka Research 3 can be controlled via a single Franka UI connection or field bus. Single Point of Control (SPoC) mechanisms ensure control by one source only. Field buses are covered by SPoC as well.

For more information about SPoC, please refer to Chapter 4.2: Single Point of Control (SPoC) in the respective Operating Manual corresponding to your system version (e.g., 5.6.0 or 5.8).

WARNING

Unexpected movement of Arm

Risk of severe injury, such as crushing, tearing of the skin, and puncturing from the Arm and end effectors.

- Ensure that the end effector and/or the object mass and the Center of Mass (CoM) are parameterized correctly.
- Keep out of the workspace during operation.
- In case of acute mortal danger:
 1. Press the Emergency Stop Device to stop the robot from operating.
 2. Pull or push the Arm out of the dangerous position manually.
- In case of non-life-threatening danger:
 1. Use the Emergency Unlocking Tool to move the Arm.

Surface temperature of the Arm (from base until axis 7, excluding flange)

WARNING

Hot Surfaces

Handling of the Arm's metal or plastic segments for a prolonged time after intense task execution may cause thermal burns.

- Do not touch the Arm's segments for more than 60 seconds after stopping Franka Research 3 after an intense task execution at maximum payload and extended temperature.

WARNING

Hot Surfaces and Guiding

At ambient temperatures above 30 °C the robot surface can become too hot to touch. Therefore, using the Assist feature in Execution mode is not permitted above 30 °C.

Effects depending on the task and environment.

The required safety instruction depends on risk assessment (regarding hot surfaces).

SAFETY

SAFETY- INSTRUCTION

At extended ambient temperatures between 25 °C and 45 °C (after intense task execution and once Franka Research 3 is set to “monitored stop” state), the integrator needs to implement measures and evaluate the risks of touching the Arm for a prolonged time (< 60 s) without being subjected to thermal burns (EN ISO 13732-1:2006). Measures include but are not limited to the following:

- Cool-down time for the robot.
- Switching off the robot for a certain amount of time.
- Notifying the operator.
- Marking the spots that are most likely to be hot.
- Prohibiting access to the robot.

SAFETY- INSTRUCTION

The integrator needs to implement measures for touching the surfaces of the Arm, the end effector, and the end effector's flange regarding possible heating which may lead to thermal burns (EN ISO 13732-1:2006). Measures include but are not limited to the following:

- Cool-down time for the robot.
- Switching off the robot for a certain amount of time.
- Notifying the operator.
- Marking the spots that are most likely to be hot.
- Prohibiting access to the robot.

4.7 Installation of Safety Peripherals

Emergency stop installation

The Emergency Stop Device needs to be installed according to the generally valid and accepted engineering standards, e.g., European standards EN 60204 and related.

The Emergency Stop Device provided by Franka Robotics needs to be connected to the X3.1 port. Other devices than the Emergency Stop Device provided by Franka Robotics may also be connected to the X3 port.

The devices connected to the emergency stop signal must follow EN 60947-5-5 or equivalent.

Store detached devices that no longer fulfill a safety function away from the device to prevent them from mistakenly being activated.

NOTICE

Place the connected Emergency Stop Device such that it is always reachable in case of an emergency, but that accidental use can be prevented.

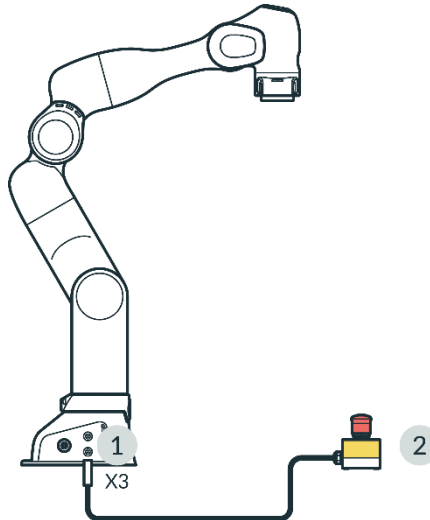


Fig. 4.1: Connection Emergency Stop Device

1	X3 - Safe inputs connector	2	Emergency Stop Device
---	----------------------------	---	-----------------------

Stopping time and distance

Stopping time (i.e., the time that passes between the request of an emergency stop and the complete stop of the Arm) and stopping distance (i.e., the distance covered by the Arm after activation of the emergency stop until it comes to a complete stop) were measured per EN ISO 10218-1, Annex B. Stopping time and distance can be found in the appendix of this document.

4.8 Fail-safe locking system

Fail-safe locking system

When the Arm is disconnected from the power supply, the locking bolts automatically lock all seven joints. The locking bolts mechanically lock any motion regarding the joints so that the Arm stays in position even when it is not supplied with power.

Due to the technology of these locking bolts, the exact position cannot be maintained when the power is cut off. The locking bolts lock with an audible click and the arm lowers a few centimeters. Especially in the joints, which are particularly influenced by gravity due to their alignment and position.

Unlocking the safety locking system

Each axis will move slightly as soon as the safety locking system unlocks.

4.9 Manually Moving the Arm

Moving the Arm without electrical power

CAUTION

Moving Arm

Risk of severe injury, such as crushing, tearing of the skin, and puncturing

- Always wear personal protective equipment (e.g., safety goggles).
- The integrator needs to perform a risk analysis for any attached end effectors.
- Do not stand in the maximum workspace during operation.

SAFETY

NOTICE

The integrator must assess the risk that a person may become trapped.

If a person is trapped by the Arm, follow one of the three release options below to free them even when the power supply is cut off.

- Use the Emergency Unlocking Tool at the corresponding openings of the joint to be unlocked to unlock the robot and move it by hand.
- Unscrew the Arm base from its mounting place.
- Manually move the Arm.

In case of non-acute danger and jamming of the Arm, use the Emergency Unlocking Tool.

WARNING

Falling heavy Arm when using Emergency Unlocking Tool

Risk of being trapped by the Arm when unlocking joints

- Do not support the Arm before and during unlocking it.
- Do not place your head or other body parts between or underneath links of the Arm.
- Do not place body parts (especially hands, fingers) between links of the Arm, the end effector, or stationary objects.
- Do not use the Emergency Unlocking Tool while the Arm is powered.
- Only use the supplied Emergency Unlocking Tool.
- Store Emergency Unlocking Tool near the Arm.

Action: Emergency unlock

SAFETY- INSTRUCTION

1. Press the Emergency Stop Device to stop the robot from operating.
2. Take the unlocking tool from the Pilot's base.
3. Hold the Arm's segments.
4. Insert the unlocking tool in the respective trapezoidal openings and unlock one or more joints one after another.

The openings are marked with the "Emergency Unlock" label.

The Arm segment can now be moved manually.

SAFETY

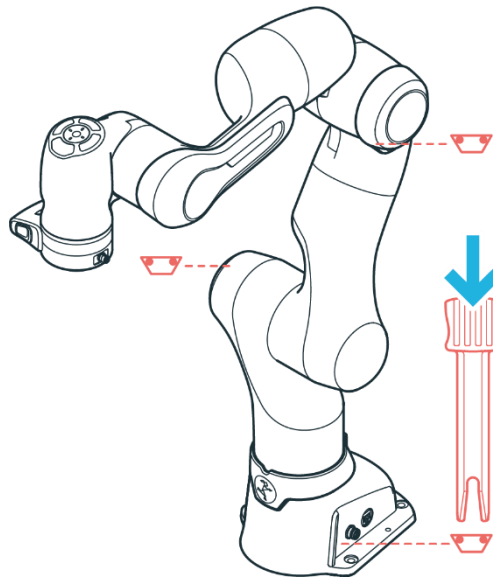


Fig. 4.2: Emergency unlock

NOTICE

Be aware that as soon as the unlocking tool is inserted, the Arm segment towards the wrist of the Arm can fall with gravity.

NOTICE

- The integrator needs to make sure that the unlocking tool is stored in a holder at the base of the robot.
- Do not remove the unlocking tool unless in an emergency.
- The unlocking tool must always be within reach.
- Only use the original unlocking tool.
- The unlocking tool should only be used in case of emergency.

Action: Pushing away manually

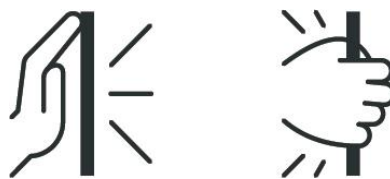


Fig. 4.3: Pushing away manually

SAFETY- INSTRUCTION

In case of acute mortal danger:

1. Press the emergency stop device to stop the robot from operating.
2. Pull or push the Arm out of the dangerous position manually.
3. Secure the robot arm from falling.

SAFETY

NOTICE

Manually pulling or pushing the Arm will damage the Arm as the joints will be overloaded.

- The Arm should only be manually pulled or pushed away in safety-critical situations.

4.10 Safety Functions

WARNING

Hot Surfaces and Guiding during Recovery

At ambient temperatures above 30 °C the robot surface can become too hot to touch. Therefore, in case of occurrence of a safety function violation which requires hand-guiding during recovery, the following must be observed:

- Recovery may be performed only by personnel specifically trained for this situation.
- Before recovery, surface temperatures must be evaluated to be within touchable limits. Cooldown times depend on previous operation and ambient temperatures.
- Wearing heat resistant safety gloves for this procedure is advised.

NOTICE

Franka Research 3 distinguishes between two kinds of safety functions: monitoring functions and stopping functions.

Monitoring functions guarantee that limits are not breached, e.g. speed (SLS-J), position (SLP-C).

Stopping functions are triggered upon a monitoring violation or a safety input. It is the Safety Operator's obligation to consider stopping times and stopping distances when configuring the limits.

NOTICE

Connecting external devices with a separate power supply may jeopardize the system's safety function if the electrical ratings are not complied with.

In addition, the voltages in the connected devices must either be SELV or suitably isolated from system-connected signals.



Safe inputs

Name	Description	Safety Rating	Stop Reaction
X3.1 - Emergency stop	The X3 connector in the robot base provides one safe input to connect an emergency stop.	PL d / Cat.3	Category 1 stop
X4 - External Enabling	The X4 connector at the robot base provides one safe input dedicated to a 3-position External Enabling Device.	PL d / Cat.3	Releasing or fully pressing the Enabling Button activates the SMSS safety function. The reaction in case of a violation of SMSS depends on the active safety scenario.
Enabling Button	A 3-position Enabling Button is provided near the flange on the Pilot-Grip of the robot.	PL d / Cat.3	Operating mode "Programming": Category 1 stop (see predefined scenario "Idle") Operating mode "Execution": Reaction depending on SMSS configuration in scenario "Work."

SAFETY

			During safety function violation or error recoveries, fully pressing or releasing the Enabling Button triggers a Category 1 stop.
X3.2 - Safe input 1 X3.3 - Safe input 2	The X3 connector on the robot base provides two additional safe inputs. The behavior of those two inputs can be configured in the safety setup.	PL d / Cat.3	Depends on the configuration in safety scenarios.


Monitoring functions

Name	Abbreviation	Description	Safety Rating	Recovery in case of violation NOTICE: The operator can recover all violations.
 When SLP-C is activated, the robot cannot be controlled by FCI!				
Safely Limited Cartesian Position	SLP-C	<p>Monitoring of the Cartesian position of certain points on the Arm. The position is checked against user-defined Cartesian area.</p> <p>The following points are monitored:</p> <ul style="list-style-type: none"> • Flange • Elbow • Wrist • Customer defined tool spheres <p>The monitoring can be configured to signal violation either when one or more points are inside the defined space or when one or more points are outside the defined limits.</p> <p>Parameterization:</p> <ul style="list-style-type: none"> • End Effector model (up to five spheres) <ul style="list-style-type: none"> ◦ Radius of each sphere ◦ Position of each sphere center relative to the flange <p>Hint: This is a general setting and will affect all safety functions using this tool model.</p> <ul style="list-style-type: none"> • Monitored Cartesian space (box) • Violation if inside/outside 	PL d / Cat.3	<p>Violated position or orientation limits will be shown in Desk of Franka UI.</p> <ul style="list-style-type: none"> • Unlock brakes of robot. • Guide robot out of Cartesian position limit. Franka UI shows if position limits are not violated anymore. • Finish recovery by pressing Confirm in Desk.
 When SLS-C is activated, the robot cannot be controlled by FCI!				
Safely Limited Cartesian Speed	SLS-C	<p>Monitoring of the Cartesian speed of certain points on the Arm structure. The following points are monitored:</p> <ul style="list-style-type: none"> • Flange • Elbow • Wrist • Centers of customer-defined tool spheres <p>Parameterization:</p> <ul style="list-style-type: none"> • Limit for Cartesian speed 	PL d / Cat.3	<p>A violation of the speed limit is shown in a dialog in the Franka UI.</p> <ul style="list-style-type: none"> • Confirm violation by pressing the button in Pop-up message. <p>No further recovery procedure is necessary.</p>
Safely Monitored Stand Still	SMSS	<p>Monitoring of the standstill in Cartesian space of certain points on the Arm structure. The following points are monitored:</p> <ul style="list-style-type: none"> • Flange • Elbow • Wrist • Centers of customer-defined tool spheres <p>The user cannot change the parameters of this safety function.</p>	PL d / Cat.3	<p>A violation dialog is shown in the Franka UI.</p> <ul style="list-style-type: none"> • Confirm violation by pressing the button. <p>No further recovery procedure is necessary.</p>

SAFETY

Safe End Effector Power Off	SEEPO	<p>Safely switch off the power provided to the end effector (48 V power line). The behavior of SEEPO can be configured in the safety setup. E.g., it can be configured that SEEPO switches off the power when an emergency stop is triggered.</p> <p>Parameterization:</p> <ul style="list-style-type: none"> General configuration, whether SEEPO is active or not Switch-off triggers of SEEPO 	PL b / Cat. b	The power of the end effector can be switched on again in Settings or the sidebar of Desk.
-----------------------------	-------	--	---------------	--

Internal monitoring functions (not parametrizable and configurable in safety rules)

Name	Abbreviation	Description	Safety Rating	Reaction	Recovery in case of violation
 When SLP-J is activated, the robot cannot be controlled by FCI!					
Safely Limited Position of Joint	SLP-J	<p>Monitoring of the position of each joint in joint space.</p> <p>This safety function is only used internally to protect the joint limits of the Arm, prevent self-collisions and local clamping. It is not available in user-defined scenarios.</p> <p>This is a limiting function.</p>	PL d / Cat.3	-	<p>A dialog within Franka UI informs the user about the violation and allows recovery.</p> <ul style="list-style-type: none"> Unlock the joint that shall be moved by clicking the unlock icon in the recovery dialog. To enable recovery motion, press the External Enabling Device. Move the joint by pressing the +/- buttons in the recovery dialog. <p>Hint: Joints in a violated state can only be moved away from violated limit. All other joints can be moved in both directions to move the robot into a more convenient pose.</p>
Safely Limited Speed of Joint	SLS-J	<p>Monitoring of the speed of a single joint in joint space.</p> <p>This internal safety function is used, e.g., to prevent fast motions during joint position recovery.</p>	PL d, Cat. 3	Cat. 1 Stop	<p>A violation dialog is shown in Franka UI.</p> <ol style="list-style-type: none"> Confirm violation by pressing the button. <p>No further recovery procedure is necessary.</p>
Safely Limited Distance	SLD	<p>SLD monitors a single joint to stay within a permitted position window.</p> <p>This internal safety function is used, e.g., to prevent excessive motion during the brake opening procedure.</p>	PL d, Cat. 3	Cat. 1 Stop	<p>A safety error dialog is shown in the Franka UI.</p> <ol style="list-style-type: none"> Acknowledge error by pressing the button. <p>No further recovery procedure is necessary.</p>

Stopping functions

Name	Description	Safety Rating
Category 0 stop	The Arm is stopped immediately by removing power from the motors and applying the brakes.	PL d / Cat.3
Category 1 stop	<p>The Arm is stopped in a controlled manner using the normal control of the motors until the standstill of each joint. The brakes are applied, and the power is removed from the Motors upon standstill.</p> <p>The Cartesian speed deceleration is monitored.</p>	PL d / Cat.3
Category 2 stop	<p>The Arm is stopped in a controlled manner using the normal control of the motors until the standstill of each joint. Upon standstill, the standstill is safely monitored.</p> <p>The Cartesian speed deceleration is monitored.</p>	PL d / Cat.3

SAFETY

Safe outputs

Name	Description	Safety Rating
Safe End Effector Power Off	Switch the power provided to the end effector (48 V power line) off.	PL b / Cat.b

Further safety ratings

The 3-stage Enabling Button near the robot's flange is realized in compliance with IEC 60204-1:2016 and IEC 60947-5-8:2006.

The 3-stage External Enabling Device provided by Franka Robotics is realized in compliance with IEC 60204-1:2016 and IEC 60947-5-8.

The Emergency Stop provided by Franka Robotics complies with IEC 60204-1:2016 and EN ISO 13850:2015.

4.10.1 Other safety-related recoveries (in case of safety errors)

Joint position error recovery

Only Safety Operators can recover joint position errors.

A dialog in Franka UI informs the user about the error and allows recovery.

Detailed instructions on how to correct the error can be found in Chapter 8 Troubleshooting in the in the respective Operating Manual corresponding to your system version (e.g., 5.6.0 or 5.8).

Safe input error recovery

Safe input errors are recoverable by confirming the respective dialog in Franka UI if confirmation for those inputs is configured in Watchman.

Other safety errors

Other safety errors are typically not recoverable. Please try to restart the system to recover from such errors. If the error persists, contact your vendor or Franka Robotics.

General info for all cases

- In case of a safety violation, the robot does not allow motion until the recovery has been accomplished.
- In case of a safety violation, the base flashes slowly in red.
- If necessary, Franka UI will show a recovery wizard to perform the recovery procedure.
- Only the safety operator can recover joint position errors.
- The operator can perform all other recoveries.

Further measures for possible troubleshooting are described in chapter in Chapter 8 Troubleshooting in the in the respective Operating Manual corresponding to your system version (e.g., 5.6.0 or 5.8).

4.11 Additional information for planning and initial installation of a robot system

According to the EN ISO 10218-2 and EN ISO 8373 standards, a robot system is understood to be a robot that is set up as a complete system with peripheral devices such as robot tools, workpieces, conveyor technology and all devices and protective equipment involved. Due to the motion of robots and the integrated applications, a robot system represents a potential hazard for those people involved with activities of operation, assembly or maintenance. The job of both the manufacturer and the installer of a robot system is to analyze and evaluate these hazards and to ensure appropriate protective measures.

This specification is based on laws, regulations and guidelines that are country-specific and thus depend on the respective location (place of the operation) of the robot.

SAFETY

In the European Economic Area (EEA), overarching regulations apply, which may be supplemented by respective country-specific legislation, industry-specific regulations and internal company regulations.

When planning a robot system, it is therefore necessary to find out about the regulations at the installation site and to take these into account accordingly.

This type of industry can also lead to different specifications. For example, will the robot system be used in industrial Research or in research?

As mentioned above, the location of the robot system determines which various rules, regulations and laws must be observed. In the European Economic Area (EEA), the Machinery Directive and harmonized European standards apply across all countries. Furthermore, local legislation such as the Product Safety Act, the Product Liability Act and the Ordinance on Industrial Safety and Health in Germany must be considered.

The most important rules and regulations for the build-up of a robot system are listed below.

Standards / Directives	Description
RL 2006/42/EG	Maschinenrichtlinie des Europäischen Parlaments und des Europäischen Rates
ISO 12100	Safety of machinery – General principles for design – Risk assessment and risk reduction
ISO 10218-2	Robots and robotic devices – Safety requirements for industrial robots – Part 2: Robot systems and integration
ISO/TS 15066	Robots and robotic devices – Collaborative robots
ISO 13854	Safety of machinery – Minimum gaps to avoid crushing of parts of the human body
ISO 13855	Safety of machinery – Positioning of safeguards with respect to the approach speeds of parts of the human body
ISO 13850	Safety of machinery – Emergency stop function – Principles for design
ISO 11161	Safety of machinery – Integrated manufacturing systems – Basic requirements
IEC 60204-1	Safety of machinery - Electrical equipment of machines - Part 1: General requirements
ISO 13849-1	Safety of machinery – Safety-related parts of control systems – Part 1: General principles for design
ISO 13849-2	Safety of machinery – Safety-related parts of control systems – Part 2: Validation
ISO 13482	Robots and robotic devices – Safety requirements for personal care robots

Basically, the same objective always applies: =>reducing the risk of injury to people.

Thus, the statement can be made that no robot system may be operated without appropriate protective measures.

Protective measures can be, for example:

- Safety protective device
- Electro-sensitive protective equipment
- Fencing and/or physical barriers
- Marked areas
- Signs

SAFETY

- Emergency stop buttons
- Indicating elements
- Control system safety devices
- Robot internal safety functionalities (*For the details please refer to section 4.10 in this Manual*)

Due to the different possible applications of a robot, Franka Robotics cannot provide a uniform guideline for determining the required protective devices when integrating a robot system. The responsibility for the safe realization of the robot system lies with the integrator / operator.

A very good and detailed guide to the design of robot systems is also described on the DGUV (German Social Accident Insurance) website => [DGUV Information 209-074](#).

The following explanations describe measures that are an essential part of planning a robot system. These measures must then be supplemented by the necessary details from the regulations and specifications of the application.

I. Analysis:

The analysis includes the description of the automation solution and at the same time a narrowing down of the function of the planned system. This part is called the intended use. The non-intended use, i.e. the conditions and activities that must not be performed with this system, must also be documented. A detailed description of the task is part of the subsequent risk assessment and simplifies the risk determination for the respective hazards in the further course of planning.

Creating a concept layout facilitates the overview of the planned system. This should show all components including the associated peripherals.

The next step is the analysis of the sources of the potential danger of the robot system. It brings clarity to the subsequent risk assessment if it is listed which component poses which hazard. If necessary, this also allows the effects of hazards to be grouped together.

All components of the robot system to be integrated are to be considered as possible sources of danger. In addition to the robot, this includes all components to be integrated, such as tools, devices, conveyor systems, control cabinets and protective devices, but also the hazards that can result from the combination of the components with each other.

II. Risk assessment:

The risk assessment is used to analyze and evaluate the sources of danger and the level of the resulting risks for personal injury, as well as to determine the necessary measures for risk reduction.

The procedure of the required risk assessment according to the Machinery Directive is described in the standard DIN EN ISO 12100. There are various tables and tools that allow a structured implementation (see reference above these listings). The basic structure of a risk assessment consists of the following elements:

- Data on the planned system (machine designation, serial number, etc.)
- Limits of the robot system
- Determination of the required regulations and standards
- Layout of the robot system
- Marking of the sources of danger in the layout
- Assessment of the sources of danger with regard to the respective activities and operating modes. For this purpose, different evaluation procedures can be applied. The procedure is explained in detail in the standard.

The procedure for determining the risk reduction measures is also described in the standard and in the guidelines and literature references.

In principle, the following prioritization of measures applies to risk reduction:

- Avoidance of the hazard
- Reduction through inherently safe design

SAFETY

- Reduction by mechanical protective devices
- Reduction through control-related protective devices
- Reduction through organizational measures

When determining the measures, the specifications from the harmonized standards should always be met. This reduces the burden of proof due to the presumption of conformity according to the standard.

III. Layout:

In the final layout of the robot system, all protective measures should be drawn to scale. There should be a clear assignment to the protective measures identified in the risk assessment.

IV. Realization phase:

Set-up of the system and implementation of the defined protective measures.

V. Verification:

After the system has been set up, including all protective measures, a verification of the protective measures must be performed in accordance with the respective standards. For example, the testing of the control technology protective measures is referred to as "functional safety verification" in standard 13849-2, and the requirements for validations are regulated therein.

This protocol of verification is an essential component for the acceptance of a robot system.

VI. Acceptance:

The final acceptance of a robot system includes a detailed logging of all individual steps mentioned above. In industrial areas of application, a declaration of conformity (CE) is required by the distributor in accordance with the Machinery Directive. A declaration of conformity (CE) is also required if a robot system is set up for "own use" in in-house Research. For robot systems in research and in the laboratory, it is also necessary to design the operation of robots to be safe for people and to implement appropriate protective measures. The Machinery Directive defines robot systems for research purposes as systems designed for a specific research purpose and built for temporary use only. Thus, the decisive factor is whether the system has temporary use (e.g., a one-time experiment that will be dismantled afterwards - no CE - or permanent use as equipment in the laboratory - CE required).

The diagram illustrates the assembly and connection of the robot system, numbered 1 through 12:

1. The robot arm is shown with joints labeled X1, X2, X3, X4, X5, and X6. Red dashed lines indicate the assembly path.
2. The robot arm is shown with the 'Left' and 'Right' labels.
3. The robot arm is shown with the 'Left' and 'Right' labels.
4. The robot arm is shown with the 'Left' and 'Right' labels.
5. The robot arm is shown with the 'Left' and 'Right' labels.
6. The robot arm is shown with the 'Left' and 'Right' labels.
7. The robot arm is shown with the 'Left' and 'Right' labels.
8. The robot arm is shown with the 'Left' and 'Right' labels.
9. The robot arm is shown with the 'Left' and 'Right' labels.
10. The robot arm is shown with the 'Left' and 'Right' labels.
11. The robot arm is shown with the 'Left' and 'Right' labels.
12. The robot arm is shown with the 'Left' and 'Right' labels.

1	Arm	7	Power cable
2	Emergency Stop Device	8	Mains power outlet
3	Connecting cable	9	Interface device (not included) with Franka UI
4	Control	10	External Enabling Device
5	Ethernet (Network)	11	Emergency Unlocking Tool
6	Power switch	12	Connection to functional earth

5.1 The Arm

The Arm has the following components:

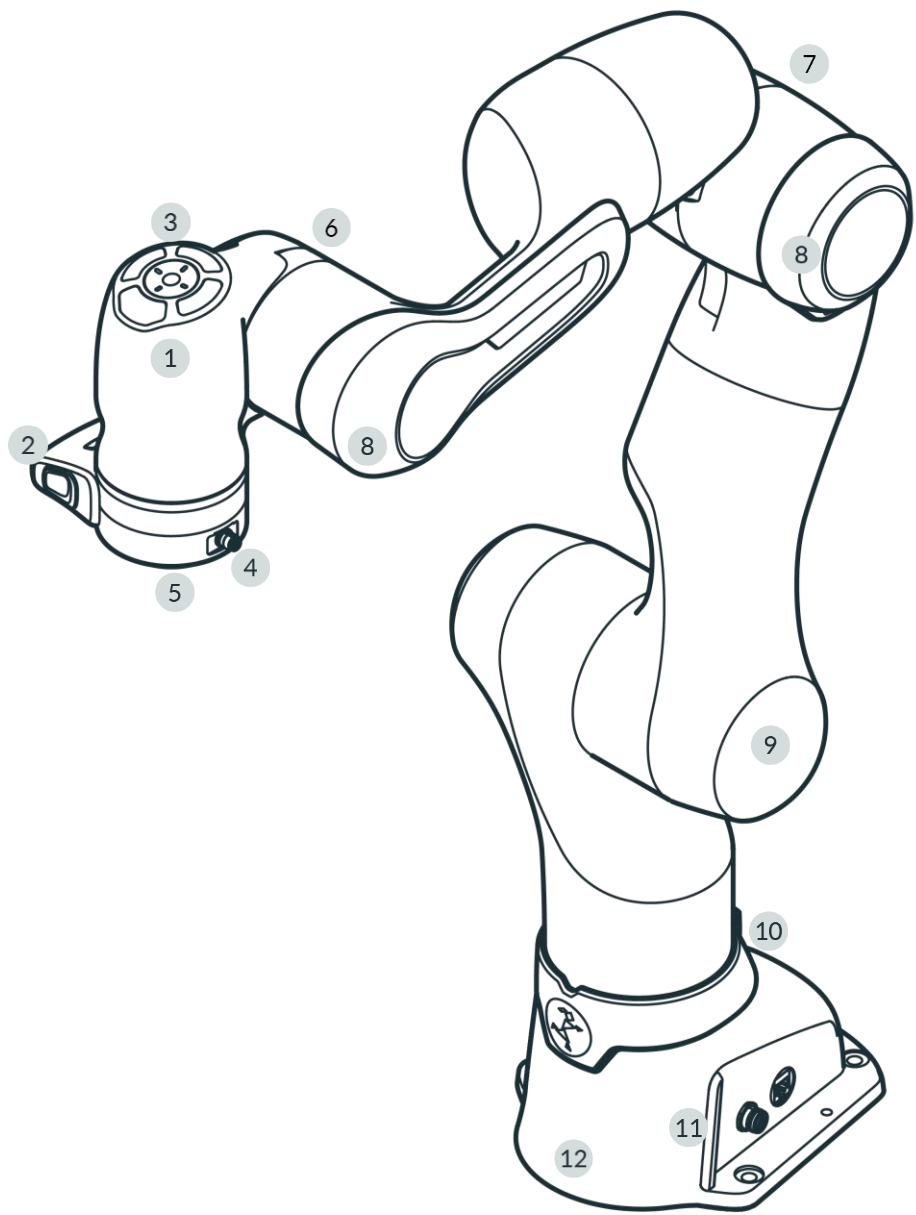


Fig. 5.2: Overview Arm

1	Pilot	7	Elbow
2	Pilot-Grip	8	Soft protector
3	Pilot-Disc	9	Shoulder
4	X6 - End effector connector	10	Holder for Emergency Unlocking Tool
5	Flange for end effector	11	Status light
6	Wrist	12	Base

EQUIPMENT OVERVIEW

Pilot - top view

Parts of Desk and the integrated end effectors can be operated directly from the robotic Arm via the Pilot-Disc.

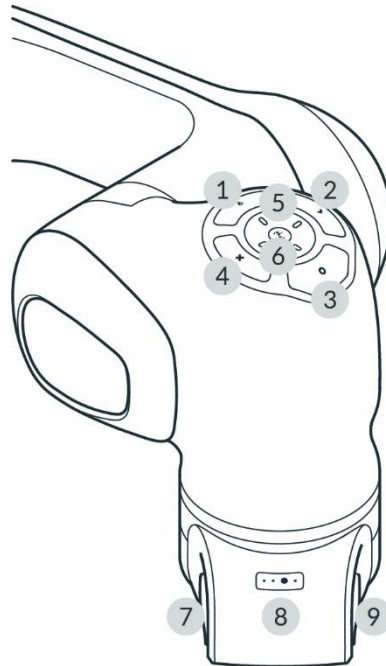


Fig 5.3: Pilot

1	Pilot-Mode button	6	Status light
2	Confirm button	7	Enabling Button
3	Teach button	8	Guiding-Mode Button
4	Delete button	9	Guiding Button
5	Arrow keys		

Pilot

The Pilot is the user interface integrated directly into the Arm for Guiding the robot and the easy interaction with end effectors and Desk. The Pilot consists of Pilot-Disc (1-6) and Pilot-Grip (7-9).

For more information on Desk, please refer to Chapter 5 Desk in the respective Operating Manual corresponding to your system version (e.g., 5.6.0 or 5.8).

Pilot-Disc (1-6)

The Pilot-Disc is located on top of the Pilot and is used for interacting with the robot system. You can switch between the control of Arm or end effector by pressing the Pilot-Mode button (1) on the Pilot-Disc. Select individual Apps, parameterize them or enter poses by manually guiding the Arm to the desired pose and pressing the Teach button (3).

Pilot-Grip (7-9)

The Pilot-Grip is located near the robot's tip as part of the robot structure. The Pilot-Grip provides a Guiding Button, an Enabling Button, and a Guiding-Mode Button.

Pilot-Mode button (1)

By pressing the Pilot-Mode button (1), the user switches between using the arrow keys of the Pilot-Disc to navigate Desk or to control the integrated end effectors (e.g., Franka Hand).

Confirm button (2)

If the Confirm button is illuminated, all changes made in the context menu will be confirmed and skipped to the next section. The Confirm button saves any selections made.

EQUIPMENT OVERVIEW

Teach button (3)

Save an arm pose or an end effector pose by moving the arm or end effector to the desired configuration and pressing the Teach button.

Delete button (4)

If illuminated, pressing the Delete button deletes a selected pose or section.

Arrow keys (5)

The Arrow keys can either navigate Desk or control integrated end effectors depending on the Pilot-Mode. In end effector mode, the key assignment depends on the active end effector.

Enabling Button (7)

The Enabling Button is located on the left side of the Pilot-Grip and enables robot motions if pressed to the middle position. To move the robot, half-press the Enabling Button while pressing the Guiding Button simultaneously. The Enabling Button is safety rated per the requirements in EN ISO 10218-1. The three positions of the Enabling Button either stop, re-enable, or move the robot. To immediately stop the robot, release or fully press the Enabling Button. After stopping, release it first completely and then press it again to its middle position to re-enable the robot.



Fig. 5.4: switching status Enabling Button

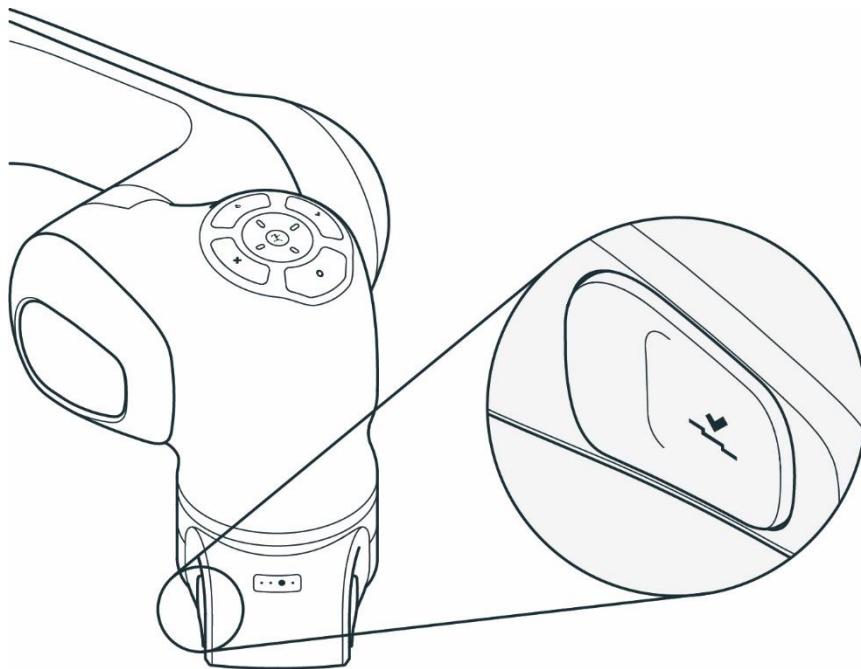


Fig. 5.5: Enabling Button

EQUIPMENT OVERVIEW

Guiding-Mode Button (8)

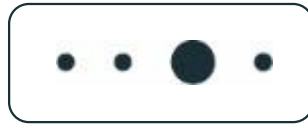


Fig. 5.6: Guiding Mode Button

The Guiding-Mode Button is located on the top of the Pilot-Grip and allows the user to change between different Guiding modalities by pressing the Guiding-Mode Button. The possible Guiding-Modes are translation only, rotation only, free moves, and user-defined movement.

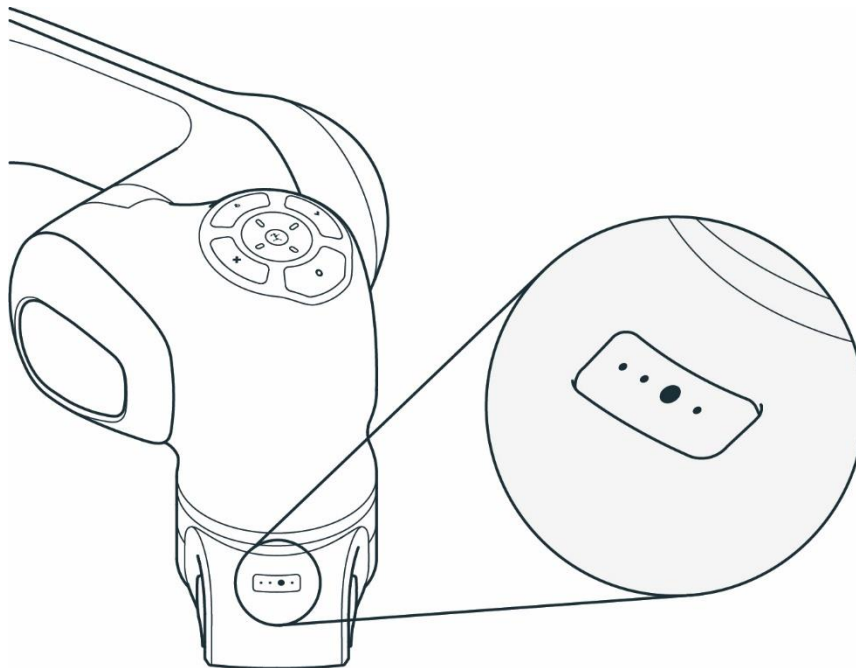


Fig. 5.7: Guiding Mode Button

Guiding Button (9)



Fig. 5.8: Guiding Button

The Guiding Button is located on the right of the Pilot-Grip. Press the Guiding Button while half-pressing the Enabling Button (7) simultaneously to move the robot.

EQUIPMENT OVERVIEW

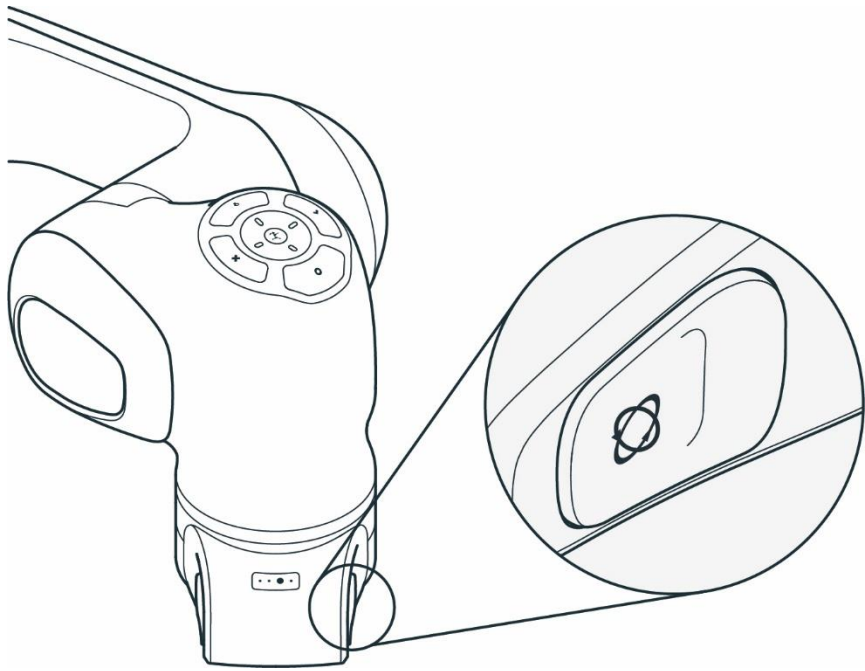


Fig. 5.9: Guiding Button

Base of the Arm

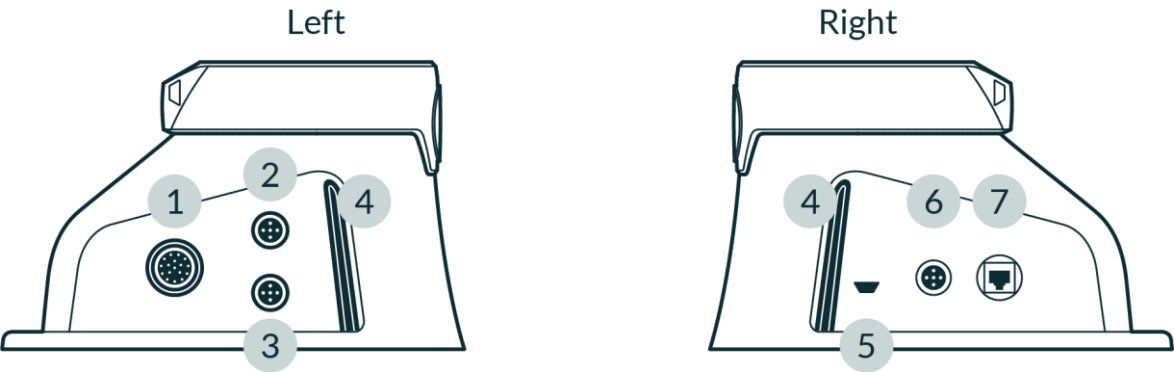


Fig. 5.10: Connection ports at robot base

1	X1 - Connection to Control connector	5	Insertion points for Emergency Unlocking Tool
2	X2 - Reserved	6	X4 - External enabling connector
3	X3 - Safe inputs connector	7	X5 - Robot network connector
4	Status light		

The base of the robot has several ports to connect and support various equipment:

- X1 - Connection to Control
- X2 - Digital IO's

NOTICE

For internal use only. This feature is not supported for customer use. No software implementation is available to enable functionality.

EQUIPMENT OVERVIEW

- X3 - Safety Signals
 - X3.1 - Emergency stop: A category 1 stop command will be triggered to stop the robot and optionally cut the power to the end effector by pressing the Emergency Stop Device. This behavior can be configured in Watchman.
 - X3.2, X3.3 - Safe inputs: The X3 port allows two additional safe inputs. Their behavior can be configured in Watchman, *please refer to Chapter 5.3 in the respective Operating Manual corresponding to your system version (e.g., 5.6.0 or 5.8).*
 - X4 - External Enabling

The 3-position External Enabling Device enable "Test & Jog" when the system is in Programming mode. It allows Franka Research 3 to move. Programs can be started via Desk.

- X5 - Robot network

The interface device running the browser-based Franka UI can be connected to the X5 Ethernet port.

NOTICE

A custom connector with the corresponding safety signals must be assembled to utilize the configurable Safe inputs (X3.2, X3.3). When doing so, the existing Emergency stop can no longer be used. Therefore, the Emergency stop functionality must be integrated for channel X3.1 in the custom connector.

NOTICE

By default, the Safe input channels X3.2 and X3.3 are assigned to the SMSS safety function in the "Work" scenario in Watchman. Suppose no external safety devices are connected to X3.2 and X3.3 (only Franka Robotics emergency stop is connected to X3.1). Those inputs will be considered "activated," meaning it will not be possible to move the robot with the default "Work" scenario rules. Depending on the application-specific risk & hazard analysis, the default rules may be changed to allow robot movements regardless of X3.2 and X3.3.

End effector flange

End effectors, such as Franka Hand, can be connected via the end effector flange. The end effector flange has been developed according to the relevant quality standards of DIN ISO 9409-1-A50.

For more information, see chapter 7.7 Mounting End Effectors in this Manual.

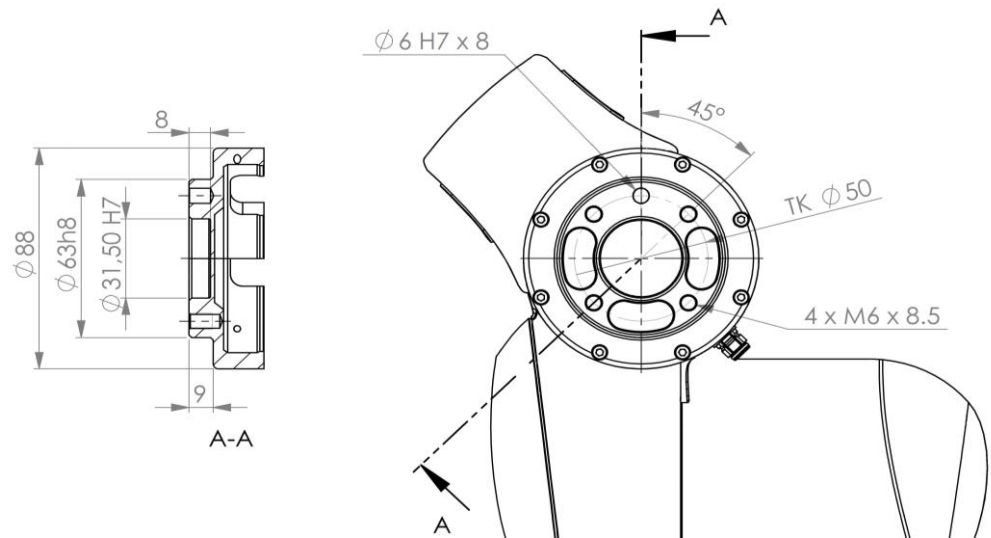


Fig. 5.11: End effector flange

NOTICE

Franka Hand is not part of the certified machinery.

5.2 Control

NOTICE

The operation of the Arm is only permitted with the Control provided by Franka Robotics.

The Control is the main control unit and is part of Franka Research 3. The main control unit realizes monitoring and control of the mechanical structure of the robot.

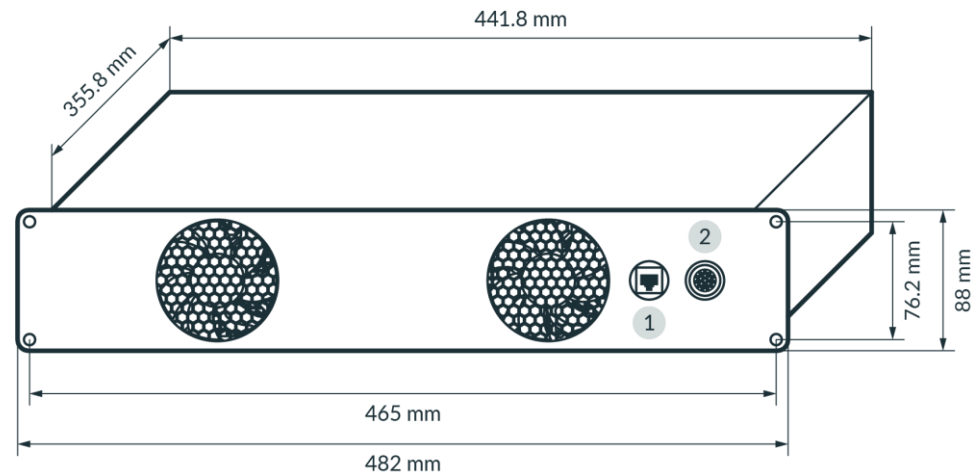


Fig. 5.12: Dimensions and connection ports of control

1	C2 - Shop floor network connector	2	C1 - Connection to Arm connector
---	-----------------------------------	---	----------------------------------

SCOPE OF DELIVERY AND ADDITIONAL EQUIPMENT



Fig. 5.:13

1	C3 – Power connector	2	Power switch
---	----------------------	---	--------------

Installation

The Control fits in a 2RU 19" rack.

6 SCOPE OF DELIVERY AND ADDITIONAL EQUIPMENT

6.1 Included in the box

Arm

- 1x Arm
- 1x Emergency Unlocking Tool
- 4x screw (ISO 4762, M8x20, ST 10.9 A2K)
- 4x washer (ISO 7089, M8, ST HV300 A2K)
- 1x screw (ISO 4762, M5x8, ST 8.8 A2K)
- 1x tooth washer (DIN 6797-A, M5, ST A2K)
- 1x Quick Guide for Installation FR3 (Document number: R02040)



Fig. 6.1: scope of delivery Arm

Control

- 1x Control
- 1x country-specific power cable

SCOPE OF DELIVERY AND ADDITIONAL EQUIPMENT



Fig. 6.2: scope of delivery Control

Devices

- 1x External Enabling Device
- 1x Emergency Stop Device

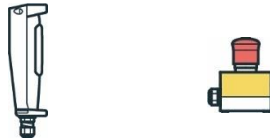


Fig. 6.3: Scope of delivery devices

Accessories

- 1x Connecting cable



Fig. 6.4: Scope of delivery Connecting cable

6.2 Not included in the box

For additional accessories, e.g., Cobot pump, visit <https://franka.world/>.

The following equipment is not included:

- Interface device
 - Tablet/Notebook/PC
The interface device should be equipped with a browser (Chrome, Chromium, or Firefox), an Ethernet port, and ideally with touch functionality.
- Material
 - Ethernet cable with RJ 45 connector for connecting the interface device to the Arm
 - Ethernet cable with RJ 45 connector for optional connection of the Control to the company network or PC workstation
 - Mounting accessories (recommended by Franka Robotics): 2x 6 mm h8 pins for precise mounting of the Arm, if applicable
 - Baseplate to mount the Arm (depending on the baseplate, different screws and washers may be needed, **see table in chapter 7.4 Mounting the Arm in this Manual.**)
 - Functional earth cable with eye
- Tools
 - Hex key for mounting the Arm on the baseplate
 - Screwdriver for connecting the functional earth cable
 - Level for ensuring the horizontal installation of the Arm
 - Torque spanners to tighten screws with 30 Nm

6.3 Available Spare Parts and Accessories

The spare parts for Franka Research 3 include but are not limited to:

- Arm
- Control incl. country-specific power cable
- External Enabling Device
- Emergency Stop Device
- Connecting cable (2.5 m, 5 m, or 10 m)
- Emergency Unlocking Tool
- Franka Hand (Not part of the certified machinery)
- Cobot pump (Not part of the certified machinery)

7 MOUNTING AND INSTALLATION

⚠ WARNING

Heavy equipment

Due to the dead weight and geometric design, lifting and handling the equipment may cause back injury, and, if it falls, serious injury to fingers, hands, toes, and feet.

- Always wear personal protective equipment (e.g., safety shoes) when transporting, mounting, or demounting the equipment.
- Always lift the equipment with the help of a second person.
- The equipment must be placed on even surfaces to prevent it from tilting or sliding.
- Follow the company regulations on lifting loads and personal protective equipment.

NOTICE

Before mounting and installing the system, please read Chapter 4 "SAFETY".

NOTICE

Arm does not stand stable without being screwed at the base.

Overview of interfaces provided by Control and Arm

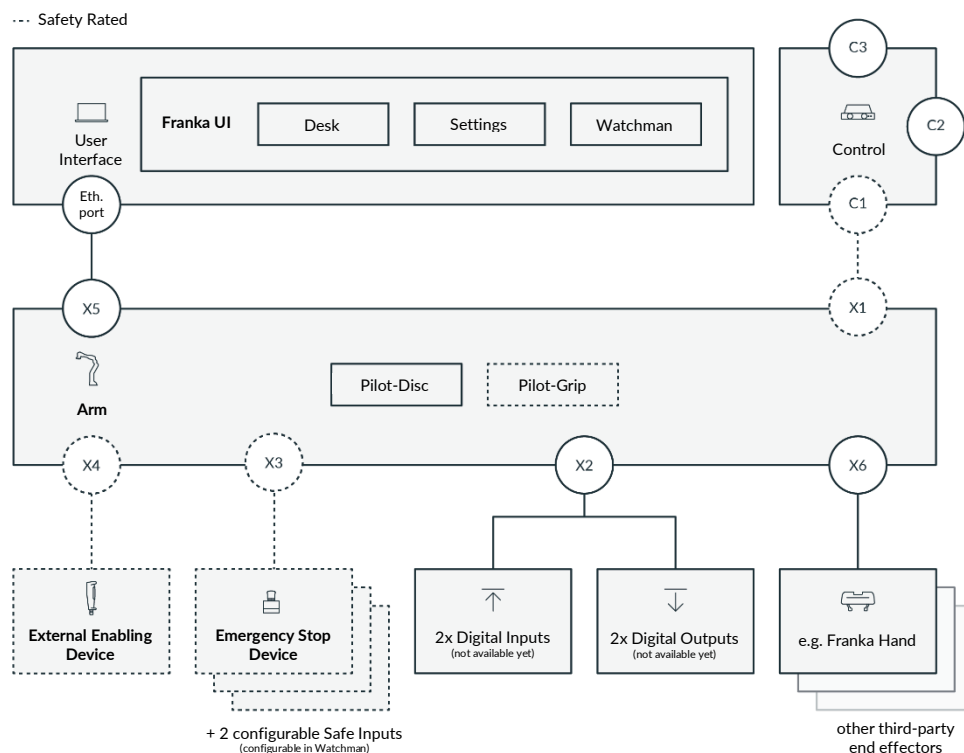


Fig. 7.1: Overview of interfaces

7.1 Unpacking the Equipment

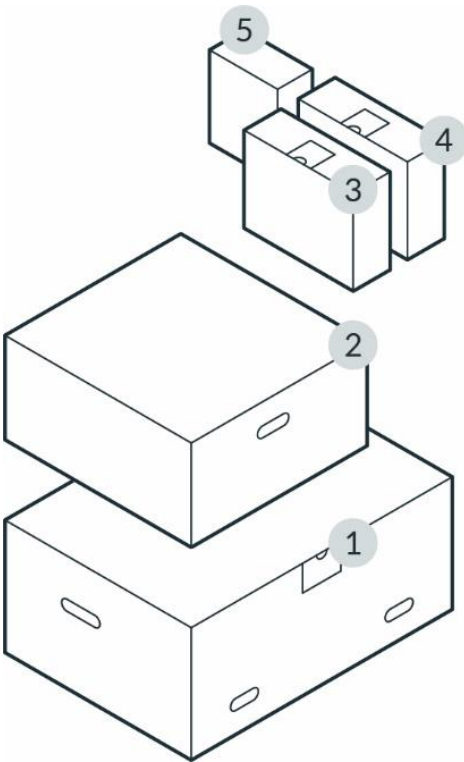


Fig. 7.2: Packing

1	Arm	4	Emergency Stop Device & External Enabling Device
2	Control	5	Optional (e.g. Franka Hand)
3	Connecting cable		

NOTICE

Always keep the original packaging in case of relocating the robot.

Unboxing

Procedure

1. Take off the top lid of the outer box.

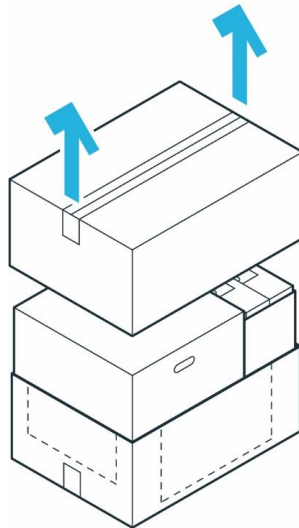


Fig. 7.3: Unpack main carton

2. Lift the top inner boxes and put them aside.

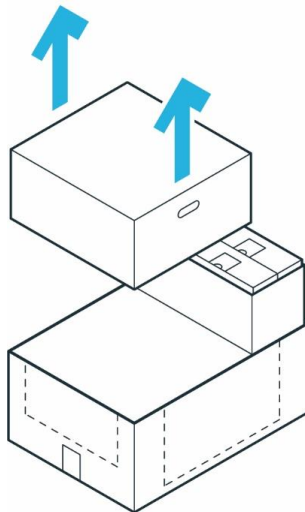


Fig. 7.4: Removal of individual cartons

3. Pull the outer box apart to access the bottom inner box.

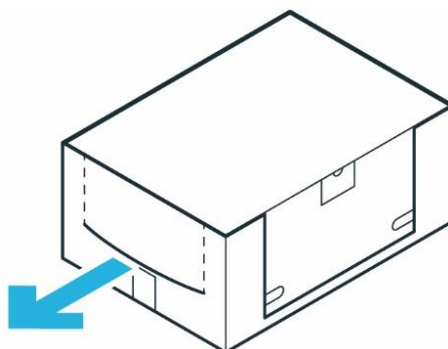


Fig. 7.5: Inner cartons

Unpacking the Arm

Procedure

MOUNTING AND INSTALLATION

1. Carefully open the box by removing the sealing adhesive strips on top of the cardboard box.
2. Open the foil coating.
3. Remove the top protective layer.

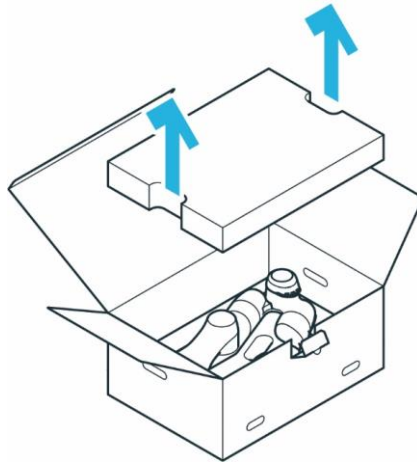


Fig. 7.6: Open box from Arm

4. Remove the middle protective layer.

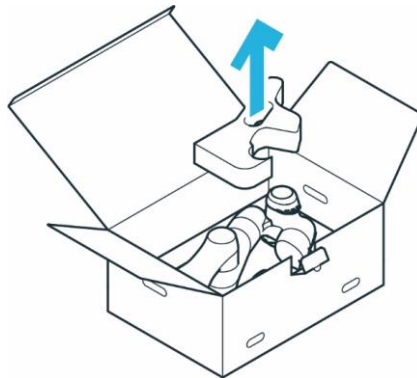


Fig. 7.7: Unpack arm

5. Grasp the Arm at the indicated lifting positions carefully, lift it out of the bottom protective layer, and set it aside.

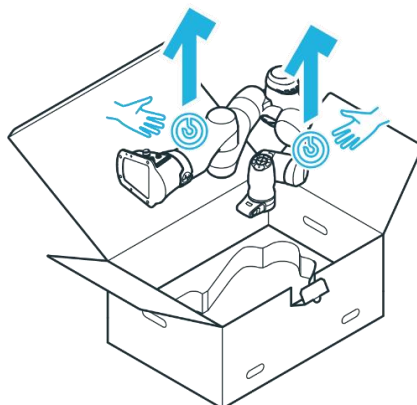


Fig. 7.8: Lift out arm

MOUNTING AND INSTALLATION

Unpacking the Control

Procedure

1. Carefully open the box by removing the sealing adhesive strips on top of the cardboard box.
2. Open the foil coating.
3. Remove the power cable and the top lid.

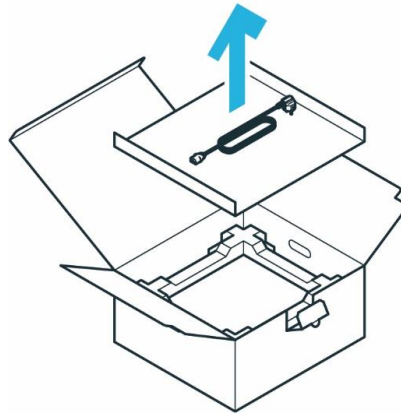


Fig. 7.9: Open box of control

4. Remove the top protective layer.

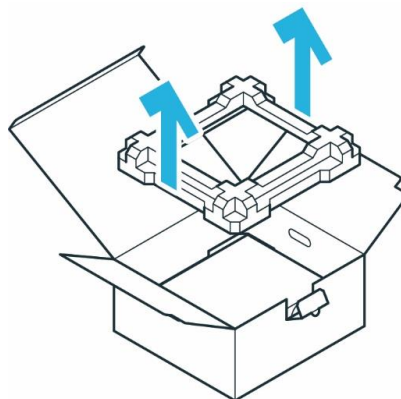


Fig. 7.10: Removal packaging

5. Grasp the Control at the indicated lifting positions, carefully lift it out of the bottom protective layer, and set it aside.

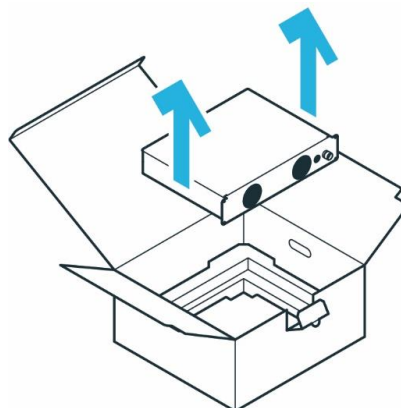


Fig. 7.11: Lift out control

7.2 Correct Installation Site

7.2.1 Maximum-and safeguarded space

Classification of spaces

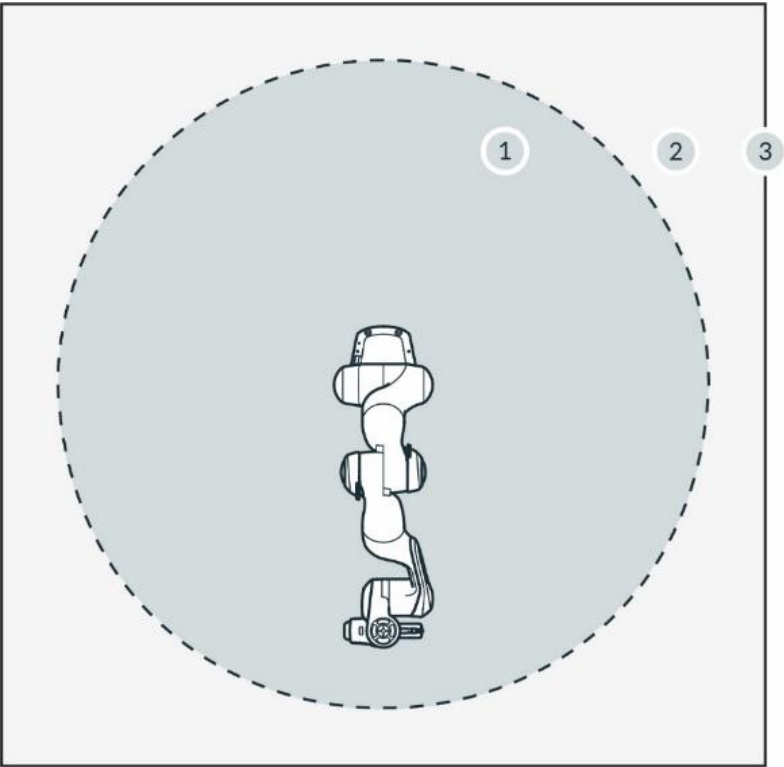


Fig. 7.12: Classification of spaces

1	Maximum workspace	3	Perimeter safeguarding
2	Safeguarded space		

- Maximum workspace
Space that can be swept by the robot's moving parts plus the space which the end effector and the workpiece can sweep.
- Safeguarded space
Defined by the perimeter safeguarding (see graphic)

NOTICE

Utilize the provided safety functions to mitigate any risks of accidental collisions from an unexpectedly moving Arm. Monitoring functions are only triggered upon violation. The safety integrator must consider stopping times, distances, and tolerances.

NOTICE

Application-specific risk analysis may define a hazardous zone larger than the maximum workspace in some applications.

MOUNTING AND INSTALLATION

7.2.2 Ambient conditions: Arm

Permissible conditions at the site of installation

Ambient temperature

- +15 °C to +25 °C (normal)
- +5 °C to +45 °C (extended)
- IP 20
- Normal operations (no-derating): +15°C to +25°C, 60% not condensing humidity
- Extended operations (no de-rating of safety system, de-rating of performance possible): +5°C to +45°C, 90% not condensing humidity
- Storage and transport: -10°C to +60°C

Relative air humidity

- 20 % to 80 %, non-condensing

Site of installation

- Indoors, in enclosed buildings
- Not exposed to direct sunlight
- No vibrations, no accelerating fundamentals
- Magnetic fields are only permissible in the stated specification range.

Please refer to 4.2 Notice of liability in this Manual.

Orientation of installation

- Arm may only be installed vertically (base horizontal to the earth's surface, no hanging Arm)

Ambient medium

- Air
- Free from flammable substances (dust, gas, liquid)
- Free from aggressive media
- Free from corrosive substances
- Free from flying objects
- Free from spraying liquids
- Free from pressurized air streams

Pollution degree

- Degree 2 (according to EN 60664)
- Only dry, non-conductive pollution occurs; occasionally, temporary conductivity caused by condensation may occur

Setup altitude

- ≤ 2,000 m above sea level

Electromagnetic compatibility

- Environmental conditions must adhere to general industrial equipment according to EN 61000-6-4 as the system is designed for respective emission tolerance according to EN 61000-6-2.

NOTICE

In order not to endanger the safety functionality of the system, ensure pollution degree 2 according to EN 60664.

MOUNTING AND INSTALLATION

Adequate ventilation

NOTICE

The heat produced by power electronic components and modules inside the Arm is dissipated via the Arm's surface.

- Install the Arm in an adequately ventilated place.
- Do not expose the Arm to direct sunlight.
- Do not repaint, paste up, or wrap up the Arm.

Ergonomic consideration

NOTICE

To avoid overheating, the system will stop working when it exceeds the extended temperature range. The user will be informed through Franka UI.

Follow further instructions in Franka UI.

NOTICE

To avoid overheating of the motors, the system will stop working if the internal sensors detect excess temperatures in the windings. The user will be informed through Franka UI.

Follow further instructions in Franka UI.

NOTICE

Install the Arm in an ergonomic teaching position.

7.2.3 Ambient conditions: Control

Permissible conditions at site of installation

Ambient temperature

- +15 °C to +25 °C (normal)
- +5 °C to +45 °C (extended)

Relative air humidity

- 20 % to 80 %, non-condensing

Site of installation

- Indoors, in enclosed buildings
- Not exposed to direct sunlight
- No vibrations
- Magnetic fields are only permissible in the stated specification range.

Please refer to 4.2 Notice of liability in this Manual.

- Enclosure needs a minimum degree of protection corresponding to IP4X or IPXXD if located in places open to all people.

Orientation of installation

- Device may only be installed horizontally on the earth's surface
- Mounting in angle brackets, e.g., under tables

MOUNTING AND INSTALLATION

- Mounting in control cabinets (2U, 4HP)

Power supply

- To ensure stability and safety of the system, provide a stable power supply that maintains enough power to the control to shut down in a controlled fashion, when the power supply is shut off.

Ambient medium

- Air
- Free from flammable substances (dust, gas, liquid)
- Free from aggressive media
- Free from corrosive substances
- Free from flying objects
- Free from spraying liquids
- Free from pressurized air streams

Pollution degree

- Degree 2 (according to EN 60664)
- Only dry, non-conductive pollution occurs; occasionally, temporary conductivity caused by condensation may occur

Setup altitude:

- ≤ 2,000 m above sea level

NOTICE

If it is not open to all people, only pollution degree 2 is relevant and needs to be ensured.

NOTICE

In order not to endanger the safety functionality of the system, ensure pollution degree 2 according to EN 60664.

The enclosure mentioned above is not suitable to protect us from higher pollution degrees. Here, an even higher IP is needed.

7.3 Preparing the Installation Site

Correct installation site

Before installation, prepare the installation site. *Please refer to Chapter 7.4 Correct installation site in this Manual.*

WARNING

Malfunctions and unexpected motions due to improper installation

Risk of severe injury, such as crushing fingers, hands, upper body, head.

- Only switch on the robot when the Arm is properly installed on the platform.
- Only install the Arm on even, unmoving, and stable platforms. Accelerations and vibrations induced by the platform are not permissible.
- Do not install the Arm hanging or on tilted or uneven platforms.
- Level the platform and install the robot in an upright position.

MOUNTING AND INSTALLATION

- Tighten the screws after 100 hours of operation with the correct tightening torque.

7.3.1 Arm

Derating

When operating Franka Research 3 within the extended temperature range, the user may need to reduce the dynamic parameters (acceleration, maximum speed, etc.) to avoid overheating the system and its components. Otherwise, Franka Research 3 stops its operation.

Stable platform

The Arm is equipped with highly sensitive sensor technology and fine-tuned control algorithms. The control algorithm requires installation on a stable, leveled, non-moving, and non-vibrating platform in an upright position. The maximum permitted tilt angle is 0.1°.

The following maximum forces must be supported during static and dynamic operation from mounting base:

- tilting torque: 280 Nm
- torque around axis: 190 Nm
- horizontally force: 300 N
- vertical force: 410 N

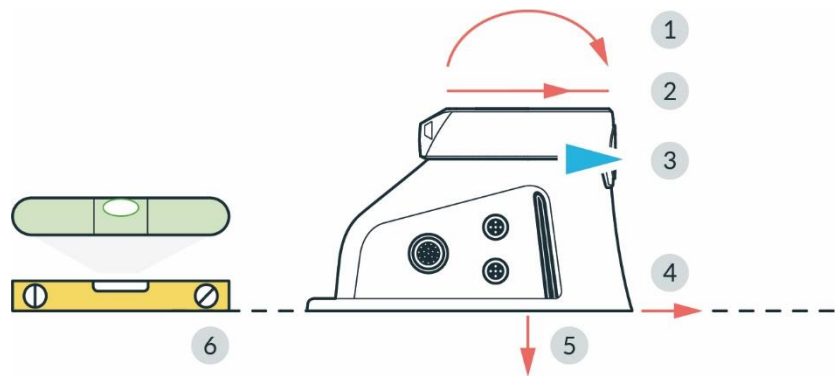


Fig. 7.13: Preparing the baseplate

1	Tilting torque	4	Horizontal force
2	Torque around axis	5	Vertical force
3	Front	6	Leveled surface

Preparing the baseplate

Required material

- Detailed mounting layout for baseplate

Procedure

- Use the technical drawing for positioning the holes.

NOTICE

The position of the Arm in the technical drawing and align it to the baseplate accordingly.

The hole spacing is designed to be compatible with flexible assembly parts by ITEM. Two holes (Ø 6 mm H7) for dowel pins in the mounting flange allow for accurate, repeatable assembly of the Arm using 2 x Ø6 h8 pins (see table in chapter 7.4 “Mounting the Arm in this Manual”).

MOUNTING AND INSTALLATION

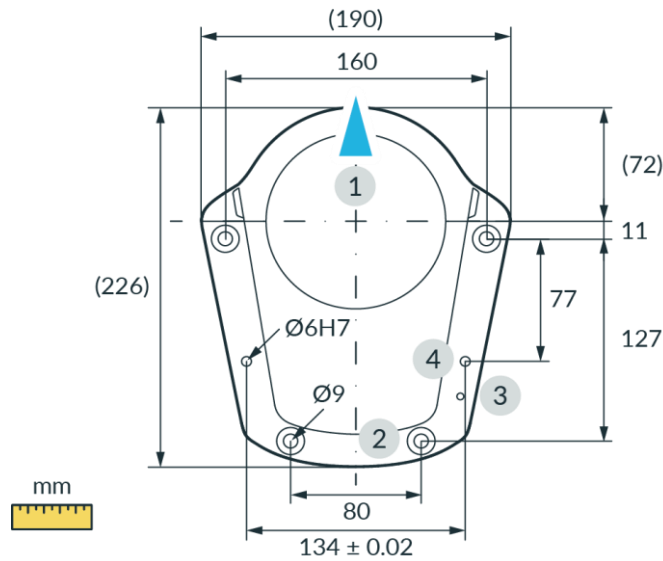


Fig. 7.14: Drilling template

1	Front	3	Thread for functional earth M5
2	Holes for M8 screws	4	Holes for alignment pins Ø6H7

7.3.2 Control

Installation site

Place the Control horizontally in its intended location.

Alternative:

Install the Control in a rack designed for 19-inch equipment.

For further information, please refer to chapter 7.2 Correct installation site in this Manual,

NOTICE

The power supply connection must be established via suitable equipment, e.g., by using the supplied country-specific cable.

Make sure that the main supply and main switch are easily accessible.

Adequate ventilation

NOTICE

The heat produced by power electronic components and modules inside the Control is dissipated through an internal ventilation system.

- Install the Control in an adequately ventilated place.
- Do not expose the Control to direct sunlight.
- Place the Control at a sufficient distance between the front/back fans and covering components (40 mm on both sides).
- Make sure that the Control's fans are not covered with dirt.

⚠ WARNING**Heavy equipment**

Due to the dead weight and partly due to the geometric design, lifting and handling the equipment may cause back injury and, if it falls, serious injury to fingers, hands, toes, and feet.

- Always wear personal protective equipment (e.g., safety shoes), when transporting, mounting, or demounting the equipment.
- The equipment must be placed on even surfaces to prevent it from tilting or sliding.
- Follow the company regulations on lifting loads and personal protective equipment.

7.4 Mounting the Arm

The Arm must be connected securely to the baseplate with four suitable-sized screws. For this purpose, four drill holes with a diameter of 9 mm are provided in the base flange of the Arm.

Use only the designated lifting points to lift the Arm.

Required tools and material

- Washers and screws depend on the surface the robot is mounted on. Please check the table below for details.
- 1x cylindrical head screw with hexagon socket M5x8 (strength class 8.8 A2K)
- 1x tooth washer M5 (strength class A2K)
- Torque spanners to tighten screws with 30 Nm

	Robot on aluminum table	Robot on steel table	Robot on ITEM aluminum profiles
Screws	ISO 4762 - M8x25 - 10.9	ISO 4762 - M8x20 - 10.9 (included in the scope of delivery)	
Washers	ISO 7089-8,4-HV300 Washers (included in the scope of delivery)		
Minimal thread length	16 mm	11 mm	Line 8 construction profiles
Tightening torque	30 Nm		
Other			Only use ITEM 0.0.420.83 Heavy Duty T-Slot M8 Nuts.

NOTICE**Material damage to the Arm**

Moving the Arm by force in a locked state will lead to a momentary slipping of internal parts which causes loss of calibration and damage to the Arm.

- Handle, lift, and transport the Arm only at the points indicated in this manual to avoid overstressing the joints of the Arm.
- The Arm is to be handled gently even when set up and switched on or off.

NOTICE

Make sure that the maximum forces and torques are supported during static and dynamic operation. For more information,

please refer to chapter 7.5 Preparing the installation site in this Manual.

MOUNTING AND INSTALLATION

Precondition

- Two people are required to mount the Arm.
- Prepared baseplate.

Please refer to chapter 7.5 Preparing the installation site section 'Preparing the Baseplate' in this Manual

Procedure

1. Lift the Arm.
2. Carry the Arm to its designated position.
3. Align the Arm according to the pre-fixed holes on the baseplate.
4. Person 1: Hold the Arm.
Person 2: Use the four screws to mount it to the baseplate with a tightening torque of 30 Nm.

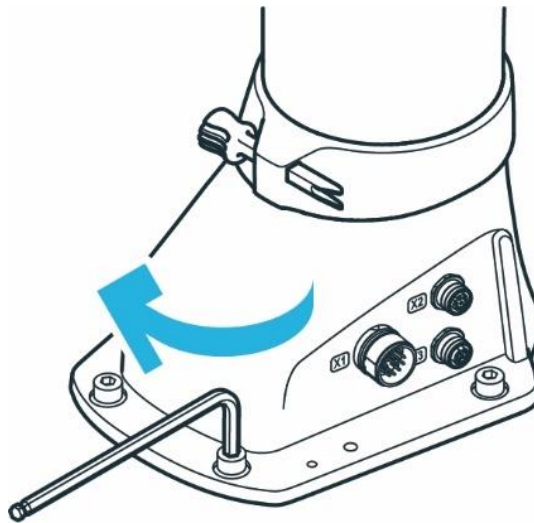


Fig. 7.15: Mounting of the arm

5. Connect the functional earth to the base of the Arm.

The Arm is successfully mounted to the baseplate.

NOTICE

The Arm should not be powered until the correct mounting is validated again.

7.5 Positioning the Control

WARNING

Heavy equipment

Due to the dead weight and partly due to the geometric design, lifting and handling the equipment may cause back injury and, if it falls, serious injury to fingers, hands, toes, and feet.

- Always wear personal protective equipment (e.g., safety shoes), when transporting, mounting, or demounting the equipment.
- The Control must be placed on even surfaces to prevent it from tilting or sliding.
- Follow the company regulations on lifting loads and personal protective equipment.

NOTICE

Material damage to Arm and Control

Moving the Arm by force in a locked state will lead to a momentary slipping of internal parts which causes loss of calibration and damage to the Arm.

- Avoid shocks.
- Carefully set down the devices.
- Always store and transport the devices in their original packaging, even inside buildings.

Positioning

Procedure

1. Person 1: Grasp the Control at the indicated lifting positions.
2. Person 2: Remove the foam packaging from the Control.
3. Place the Control horizontally in its designated position and ensure that proper ventilation is provided.

Alternative option:

Attach the Control in a rack designed for 19-inch units. **Please refer to chapter 7.4 Correct installation site in this Manual**

7.6 Wiring and Electrical Installation

Sound condition

DANGER

Damaged wires or inadequate electrical installation

Risk of personal injury by electric shock, as well as material damage

- Only use Franka Research 3 in sound technical condition.
- Only install the emergency stop and safety peripherals system with qualified personnel.
- Check cables and electrical installations.

CAUTION

Exposed wires and cables

Operators may trip and fall due to exposed wires and cables in the maximum workspace. Therefore:

- Always lay cables safely.

NOTICE

Only devices with galvanic isolation up to 60 V in the Ethernet port may be connected to the system.

NOTICE

Do not swap or unplug the connected Arm while the Control is powered on.

7.6.1 Connection diagram

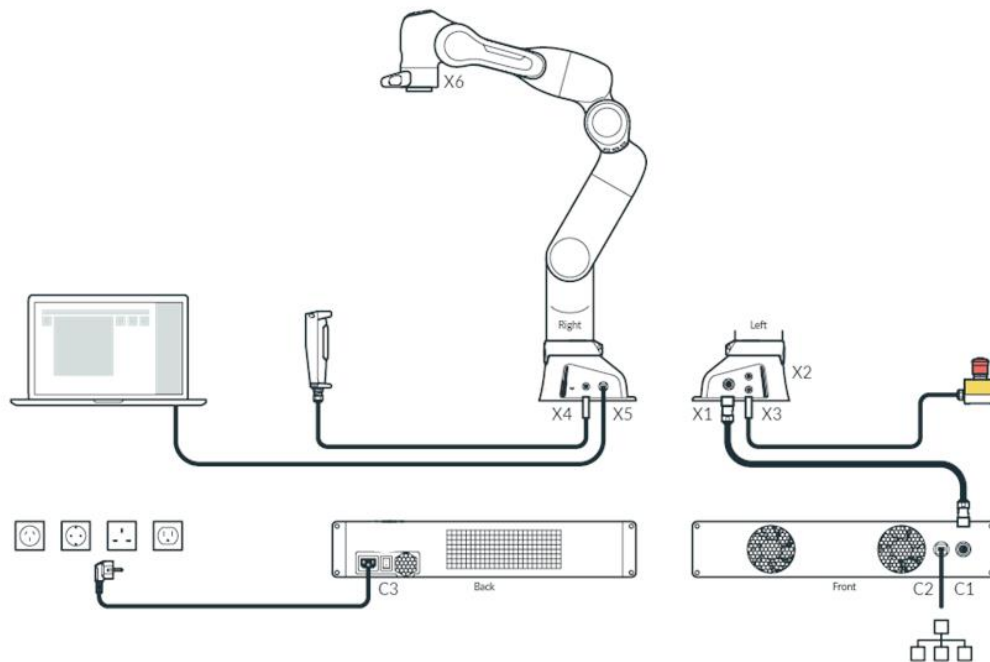


Fig. 7.16: Overview connection diagram

7.6.2 Interfaces

X2 - Digital IOs

The Interface X2 is located at the robot's base. This provides a hardware interface for non-safe digital inputs and outputs. This interface is not enabled in all - especially older - software versions.

The inputs and outputs and the 24 V power supply provided on X2 are galvanically isolated from the robot system and other interfaces at the robot.

The connector is a female 8-pin M12 connector with A coding.

The 24V power supply is driven by robot. Feeding power into these pins is neither necessary nor permitted. The maximum total output current for the 24 V pins and the digital outputs is 500 mA.

Input characteristics

- 24 V input according to IEC 61131-2 type 3
- 15 kHz low-pass input filter
- Sampling frequency 1 kHz

Output characteristics

- Output high level 24 V DC; Output current per output max. 250 mA

MOUNTING AND INSTALLATION

- Update frequency 1 kHz

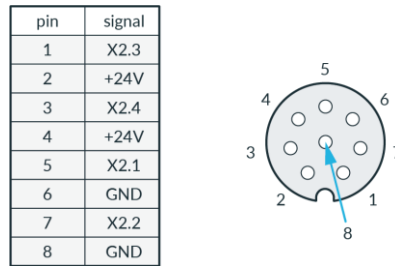


Fig. 7.17: X2 interface

X3 - Safe inputs

Safe inputs (X3.n) always have two separate channels, designated A and B. Each channel is implemented via its p- and n-pins which must be connected via a floating switch. During regular operation, both channels must be in the same state (open/closed) and not be connected; any other state will trigger a failure in the safety system.

The safe inputs are galvanically isolated from the robot system and other interfaces at the robot, although all safe inputs share a common electrical domain, regardless of their interface connector.

Interface X3 is located at the base of the robot arm and carries three safe input signals. X3.1 provides the emergency stop integration of the robot, X3.2 and X3.3 provide two freely configurable safety inputs. The connector is a female 12-pin M12 connector with A coding.

Safe Inputs electrical domain characteristics for X3:

- signal voltage 24 V; signal current 30mA

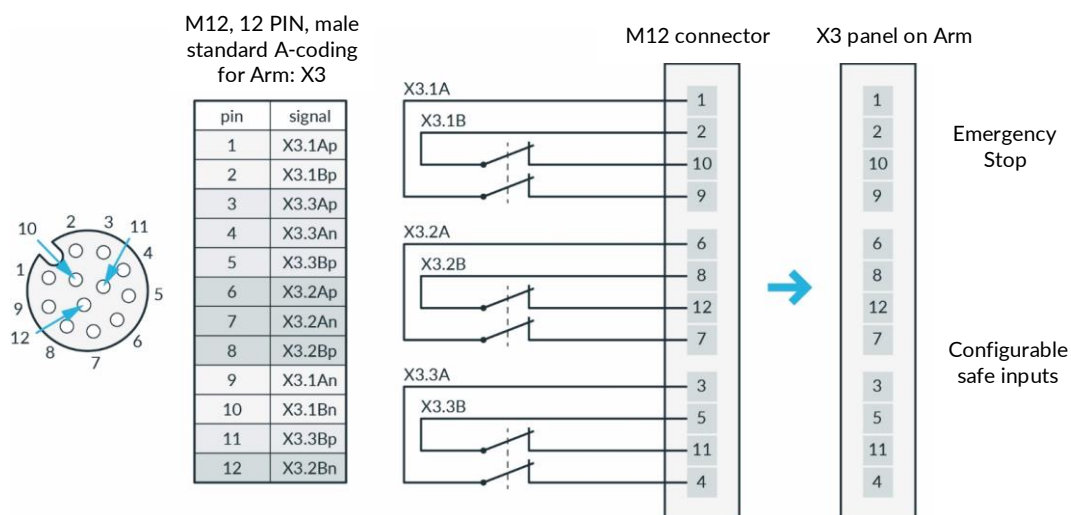


Fig. 7.18: X3 - Safe Inputs

X4 - External enabling

Interface X4 is located at the base of the robot arm and carries one safe input signal. The connector is a female 4-pin M12 connector with A coding. This connector is intended for the temporary connection of the External Enabling Device during operating phases when this is required.

If possible, always use the external enabling device supplied.

Notice

If a separate external enabling device is to be used, the external enabling device must comply with IEC 60204 1 and DIN EN 60947 5 8.

MOUNTING AND INSTALLATION

X5 - Robot network

The X5 connector is located on the robot base and provides the internal robot network via an Ethernet socket. The robot network has an integrated DHCP server. The operating device can be connected to X5. By entering the URL robot.franka.de, you can access the Franka UI web interface of the robot. The IP address of the X5 interface can be configured in the settings.

The default setting has the stored IP address 192.168.0/24. The robot can then be reached under the IP address 192.168.0.1. The DHCP server assigns addresses in the range 100 to 150 to the clients, i.e. with default settings 192.168.0.100 to 192.168.0.150.

X6 - End effector

Interface X6 is located at the wrist of the robot arm and carries signals from the robot connecting to the end effector. The connector is a Binder 8-pin female Snap-in IP67 Series 620.

pin	signal
1	48V
2	CAN_H
3	CAN_L
4	reserved
5	reserved
6	reserved
7	reserved
8	GND



Fig. 7.19: X6 interface

- Nominal voltage 48 ± 3 VDC.
- Nominals hold current 0.5 A at 25°C.
- Maximal capacitive load 220 μ F.

The end-effector interface does not transmit any safety-related information. It provides neither discrete signals nor protocol-based safe data transfer. When SEEPO (Safe Electrical Power Off) is active, no 48 V power supply is available on this interface. Additionally, the end-effector supply does not include an earthing connection.

If additional devices are connected, please test whether the intended function works as expected.

C2 Network connection

The C2 interface is located on the front of the control unit. It provides an Ethernet connection via which the control unit can be connected to a system/company network and to the Internet.

The network connection can be configured in the settings. The DHCP client is activated for this interface in the default settings. It is also possible to manually set the network connection for integration into an existing network. Please note that the robot network and the company network must not have an identical IP address range.

7.6.3 Connecting the functional earth

NOTICE

It is required to connect the functional earth to meet the stated EMC levels.

Required material

- M5 screw thread
- 1x tooth washer M5
- Functional earth cable

We recommend using a minimum 1.5 mm² section Cu cable, with a maximum length of 5 m.

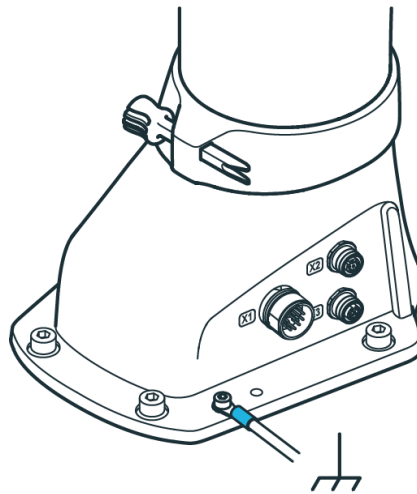


Fig. 7.20: Connection with the functional earth

Procedure

1. Connect the functional earth to the M5 thread of the Arm's base at the indicated position on one side and to a nearby, well-grounded part (e.g., a solid metal grounding bar) on the other side.
2. Position the tooth washer M5 at the indicated position at the Arm's base for the functional earth.
3. Secure the cable lug of the functional earth cable with the M5 screw.
4. Connect the other side of the cable to a nearby, well-grounded part (e.g., a solid metal grounding bar).

NOTICE

The electrical safety of the system does not depend on a functional earth connection. A functional earth connection is not suited to propagate protective bonding to attached equipment like end effectors. All equipment in the vicinity of the robot must be installed in accordance with their respective electrical requirements, including protective bonding, if applicable.

7.6.4 Wiring

NOTICE

The Arm connection cable, emergency stop cable, External Enabling Device cable, and user specific cables shall not be extraordinarily subjected to the following:

- Mechanical handling and dragging across rough surfaces (abrasion)
- Operation without guides (kinking)
- Guide rollers and forced guiding, being wound and re-wound on cable drums (stress)
- High tensile stress, small radio, bending into another plane and/or frequent duty cycles

Connecting Arm to Control

Required material:

- Connecting cable

NOTICE

Only connecting cables provided by Franka Robotics shall be used for electrical connection between Arm and Control.

Procedure

1. Carefully place the connector port (female site) onto the connector X1 and ensure that the triangular marking points upward.

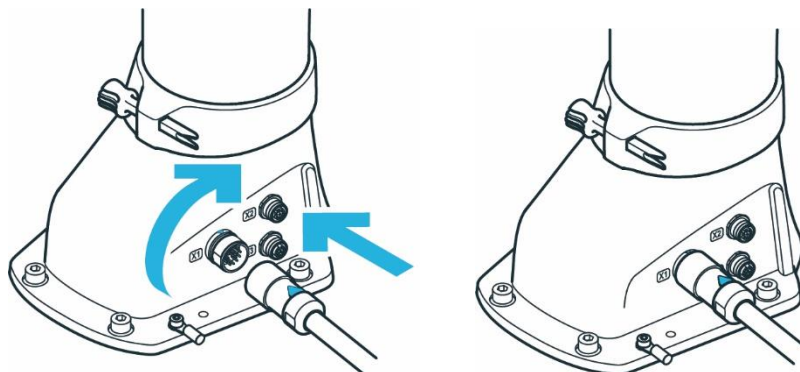


Fig. 7.21: Connection of connecting cable the arm

2. The plug itself is pulled into the connector port by turning the movable front part of the connector.
3. Turn hand-tight and test the correct fit by slightly pulling the plug.
4. Apply the same principle to connect the other end of the connecting cable (male site) with the connector C1 on the front of the Control.

MOUNTING AND INSTALLATION

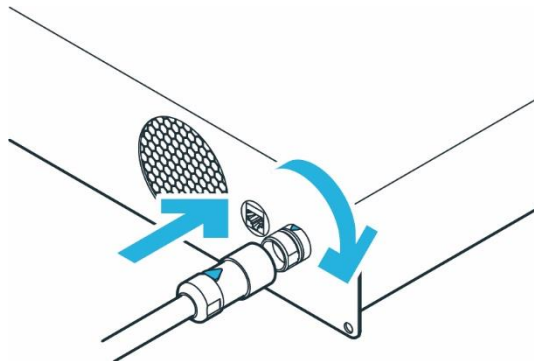


Fig. 7.22: connection of connecting cable to the control

Connecting External Enabling Device

Required material:

- External Enabling Device supplied

Procedure

1. Make sure that the guide pin is pointing in the right direction.
2. External Enabling Device to the X4 connector.

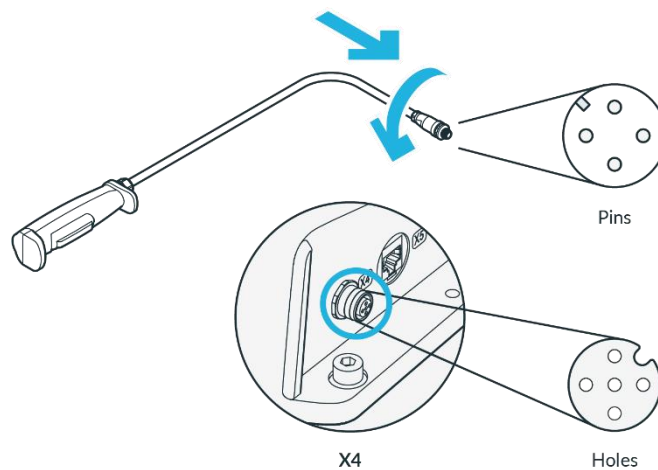


Fig. 7.23: Connection of external Enabling Device

3. The plug itself is pulled into the connector port by turning the movable front part of the connector.
4. Turn hand tight.

Connecting your operating device (for operation via Franka UI)

Required material:

Interface device **Please refer to Chapter 2.1.1 Connecting a user interface device in the respective Operating Manual corresponding to your system version (e.g., 5.6.0 or 5.8).**

- Ethernet cable with RJ 45 connector (not included)

Procedure

- Connect your interface device and the connector X5 at the Arm base with the Ethernet cable.

MOUNTING AND INSTALLATION

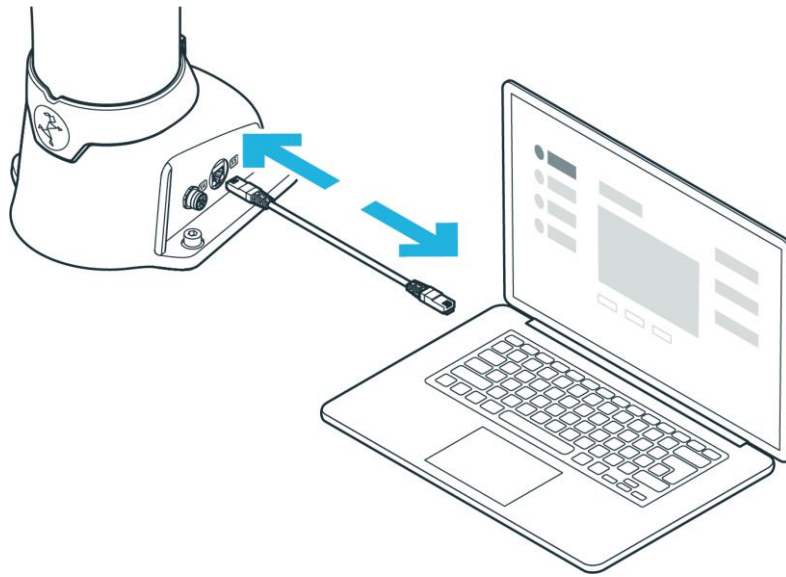


Fig. 7.24: Connection of operating device

Connecting the Control with the power supply

NOTICE

Make sure to connect the mandatory locking mechanism to the C14 connector on the power supply cable.

NOTICE

Allowed supply frequency: 50 – 60 Hz

Supply voltage: 100 - 240 VAC

Earth leakage: < 10 mA

Required material:

- Country-specific power cable

Procedure

1. Connect the power cable to the Control.
2. Connect the power cable to the power supply.

Connecting protective devices

If you want to connect external safety devices to slow down the arm and / or bring it to a standstill by means of category 1 or 2 stops (according to IEC 60204 1).

Please refer to Chapter 4.7 Installation of safety peripherals in this Manual.

MOUNTING AND INSTALLATION

WARNING

Risk of injuries

Connecting external devices with a separate power supply may jeopardize the system safety functionality.

Risk of severe injury, such as crushing, tearing of the skin, and puncturing from the Arm and end effectors.

- Ensure that the voltages in the connected devices are either SELV or suitably isolated to system-connected signals.

NOTICE

Material damage

Connecting external devices with a separate power supply may cause system damage if not adhering to electrical ratings.

- Voltages in connected devices must either be SELV or suitably isolated to system-connected signals.

NOTICE

Material damage to cables

Improper handling of cables leads to damage to cables.

- Do not bend, fold, or roll the connection cable.
- Lay the cable such that it will not be overstressed.

NOTICE

Material damage of Arm or end effectors

Unsafe connection or disconnection of live cables or end effectors during operation leads to equipment damage.

- Do not connect or disconnect cables when Franka Research 3 is connected to the power supply.
- Do not connect or disconnect end effectors when Franka Research 3 is connected to the power supply.

Using the supplied emergency stop device

Required material

- Supplied emergency stop device or customer-provided protective device (not included in scope of delivery)
- When using a customer-provided protective device: an additional connecting cable (not included in scope of delivery)

Procedure

1. connect the supplied emergency stop device to the X3 terminal.

or

Connect the customer-specific configured matching connector to the X3 connection and the protective devices to be connected (not included in the scope of delivery). Connect the protective device to the X3 connector.

MOUNTING AND INSTALLATION

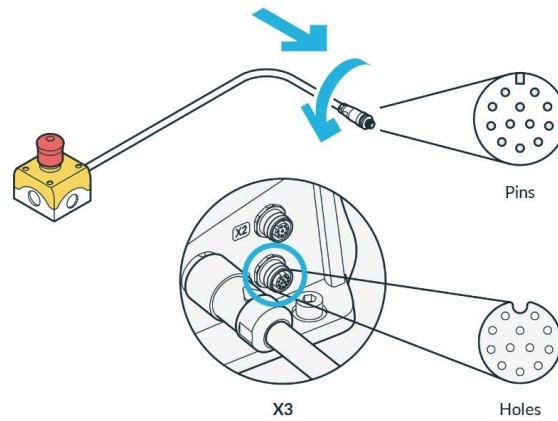


Fig. 7.25: Connection of the protective device (here emergency stop command device)

2. The plug itself is pulled into the connector port by turning the movable front part of the connector.
3. Turn hand-tight

For more information on safe input, please refer to Chapter 4.10 Safe input in this Manual.

For more information on safe input, please refer to Chapter 4.10 Safe functions in this Manual.

For more information, please refer to Chapter 4.7 Installation of safety peripherals in this Manual.

NOTICE

Safety devices need to be checked for proper functioning before initial operation and at regular intervals.

7.7 Mounting End Effectors

WARNING

Falling and/or flying tools from the end effectors

Tools that remain inserted in the end effector may turn into projectiles during later motions of the Arm and lead to injuries.

- Do not leave any tools inside the robot.

CAUTION

Sharp-edged, pointed designs and moving parts

Attached end effectors may cause injuries to hands, fingers, upper body, and head.

- Always wear personal protective equipment (e.g., safety goggles).
- The integrator needs to perform a risk assessment of any attached end effector.
- Do not stand in the maximum workspace during operation.

MOUNTING AND INSTALLATION

NOTICE

Connecting external devices with a separate power supply may jeopardize the system's safety if it does not adhere to electrical ratings.

In addition, the voltages in the connected devices must either be SELV or suitably isolated from system-connected signals.

The arm has a flange for the mechanical connection of an end effector.

The information on the end effector flange can be found in Chapter 7.3.1 Arm in this Manual.

NOTICE

Note that the pilot grip has a 45° twist to the end effector flange.

Interface X6 - End effector

The X6 electrical connection on the end effector flange can be used to supply power to the end effector if required and to communicate with the controller via the can bus. Note that the X6 connection has been configured specifically for Franka Hand and other devices may not be compatible with this connection. If an end effector is to be coupled that cannot be directly connected to such a port, external wiring for power supply and control of the end effector can be designed and implemented accordingly.

For detailed description of the wiring of interface X6 please refer to Chapter 7.6 Wiring and electrical installation in this Manual.

CAUTION

Opening and closing end effectors

Malfunctions of the Control may lead to unexpected opening and closing of end effectors.

- Always wear personal protective equipment (e.g., safety goggles).
- The integrator must perform a risk assessment on any attached end effector.
- Do not stand in the maximum workspace during operation.

CAUTION

Moving Arm

Risk of severe injury, such as crushing, tearing of the skin, and puncturing

- Always wear personal protective equipment (e.g., safety goggles).
- The integrator must perform a risk assessment on any attached end effector.
- Do not stand in the maximum workspace during operation.

NOTICE

After connecting an end effector, a risk assessment needs to be performed. The risk assessment depends on the end effector and includes but is not limited to the following:

- Sharp-edged or pointed end effectors
- Movement or rotation of sharp rotating end effectors

MOUNTING AND INSTALLATION

- An unexpectedly moving Arm leads to the end effector impacting or crushing a human

An additional risk assessment is required for risks of a possible end effector failure. The failure risk assessment depends on the end effector and includes but is not limited to the following:

- Malfunction of the protective stop signal leading to not stopping of the end effector opening/closing
- Evaluation of power-loss of the end effector and its functions
- Control failures leading to unexpectedly opening/closing of the end effector



For information on mounting and de-mounting the end effector, please see the respective end effector manual.

7.8 Practical Tips for Usage and Positioning of Franka Research 3

7.8.1 Energy consumption

For standard operation, Franka Research 3 requires an average electrical power of 140–350 W. Temporarily, electrical power of up to 600 W can be drawn from the power supply.

NOTICE

In case of unexpected power loss, Franka Research 3 attempts a Cat. 1 Stop. If the stored energy does not suffice, a Cat. 0 Stop will be performed.

After an emergency stop, the Arm could have lost calibration or been damaged. If malfunctions are detected at the next start-up, the user will be informed and should follow the instructions in Desk.

7.8.2 ESD limits

NOTICE

It is required to connect functional earth to meet stated EMC levels.

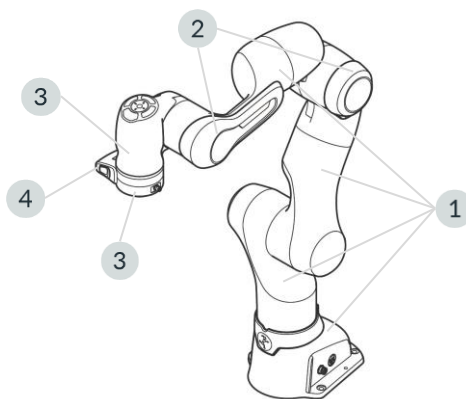


Fig. 7.26: Measuring points ESD measurement

The values in the table were measured at a temperature of 24.2 °C and relative humidity of 44 %.

	Resistance to ground R_G [Ohm]	Surface voltage [V]	Distance to ESDS [mm]
Coating robot (1)	$15,5 \times 10^9$	13	0

MOUNTING AND INSTALLATION

Forearm bumper (2)	$20,3 \times 10^9$	30	0
Wrist shell incl. Bumper (3)	$24,3 \times 10^9$	850	25
Flange (4)	$50,0 \times 10^3$	0	0
Pilot-Grip (5)	$25,7 \times 10^9$	279	25
Franka Hand (without fingertip) (6)	$38,1 \times 10^9$	615	25

The specified distances correspond to the DIN EN 61340-5-1 standard. Above 125 V to 2.000 V require 25mm spacing.

NOTICE

It is recommended to check the distances according to the application requirement and the DIN EN 61340-5-1 standard.

7.8.3 Design of the workspace

⚠ WARNING

Unexpectedly moving Arm

Risk of severe injury, such as crushing fingers, hands, upper body, head.

- Do not have sharp edges in the maximum workspace.
- Do not keep any pointed objects in the maximum workspace.
- Install the Arm in an ergonomic teaching position.



Fig. 7.27: Design of the workspace

When planning the installation, ensure that there is sufficient free workspace around the robot.

NOTICE

For recovery procedures in the event of a failure, it may be necessary to move the joints to the reference positions of the arm. It is therefore recommended to consider the position shown below in the setup planning so that the robot can reach the reference position if needed.

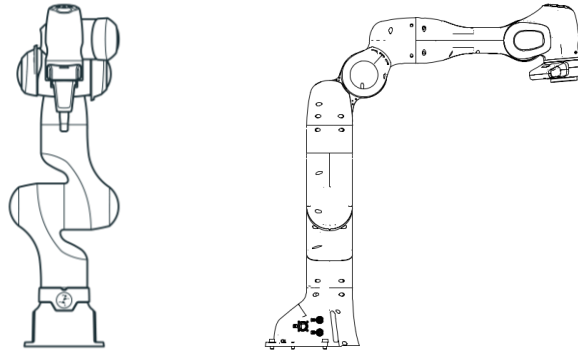


Fig. 7.28: Joint reference position of the arm

7.8.4 Personal safety and ergonomics

Free space for retracting

NOTICE

The following information on how to use and place the Arm is practical tips and may not be exhaustive when it comes to a specific application. They do not replace hazard and risk assessment but may suggest layout options.

Humans instinctively retract from unexpected motions. Therefore, the area where the operator or other persons stand should allow for sufficient space to retract or recoil.

Furthermore, ensure that this space is free of obstacles (e.g., cables, objects) to prevent people from tripping over them and harming themselves.

Greatest possible distance to the Arm

⚠ WARNING

Moving Arm

Risk of being trapped by the Arm.

- Keep the Arm at the greatest possible distance at any given time to allow the operator to react and retract.
- Do not operate the Arm while embracing it.
- Do not place your head or other body parts between or underneath segments of the Arm.
- Do not place body parts (especially hands, fingers) between the Arm, the end effector, or stationary objects.
- In case of acute mortal danger:
 1. Press the Emergency Stop Device to stop the robot from operating.
 2. Pull or push the Arm out of the dangerous position manually.

MOUNTING AND INSTALLATION

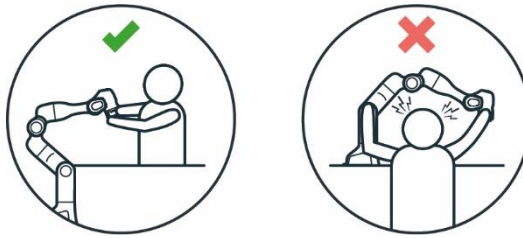


Fig. 7.29: Distance to arm bumping the head

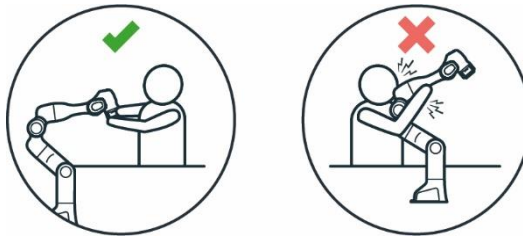


Fig. 7.30: Distance to arm bumping



Fig. 7.31: Distance to arm Squeezing the hand

Eye protection

WARNING

Unexpected movement of Arm and leaking oil

Contact with leaking oil may cause eye or skin irritation.

The use of various applications, the handled end effectors, and surrounding objects may lead to crushing, tearing of the skin, and puncturing.

- Always wear protective goggles.

Clothing and jewelry

CAUTION

Loose clothing or jewelry getting caught in Arm

Caught clothing or jewelry may cause imbalance and the risk of falling personnel.

- Do not wear loose clothing or clothing with ribbons
- Do not wear loose jewelry, e.g., necklaces or bracelets.

USAGE

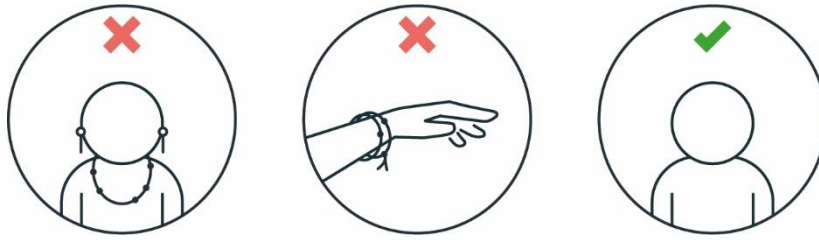


Fig. 7.32: Protective equipment does not wear jewellery

Further information

Corrosion might appear on visible surfaces. This does not affect the functionality of the robot.

NOTICE

The following information is practical and may not be exhaustive when it comes to preventing rust. In the event of rust, Franka Robotics assumes no warranty or liability, since rust does not impair the function.

- Ensure compliance with humidity and temperature ranges during use, transport, and storage.
- Store the robot in moisture-reducing materials, for example with drybags.
- Only work with clean and dry hands especially during handling, installation, and when teaching a task.

8 USAGE

8.1 Switching On

! DANGER

Short circuit due to developed condensation when the device is transported from colder to warmer and more humid surroundings

Risk of life-threatening injuries due to electrical shock.

- Leave devices to acclimatize after transport.
- Do not switch on wet devices.

Precondition

- Cables need to be plugged in correctly.
- The external power supply must be connected.
- Leave the maximum workspace.

Procedure

1. Switch on Control.

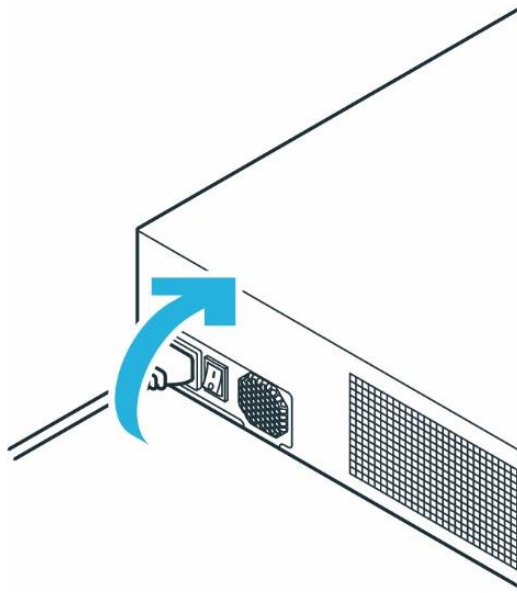


Fig. 8.1: Switching on Control

Control is now switched on.

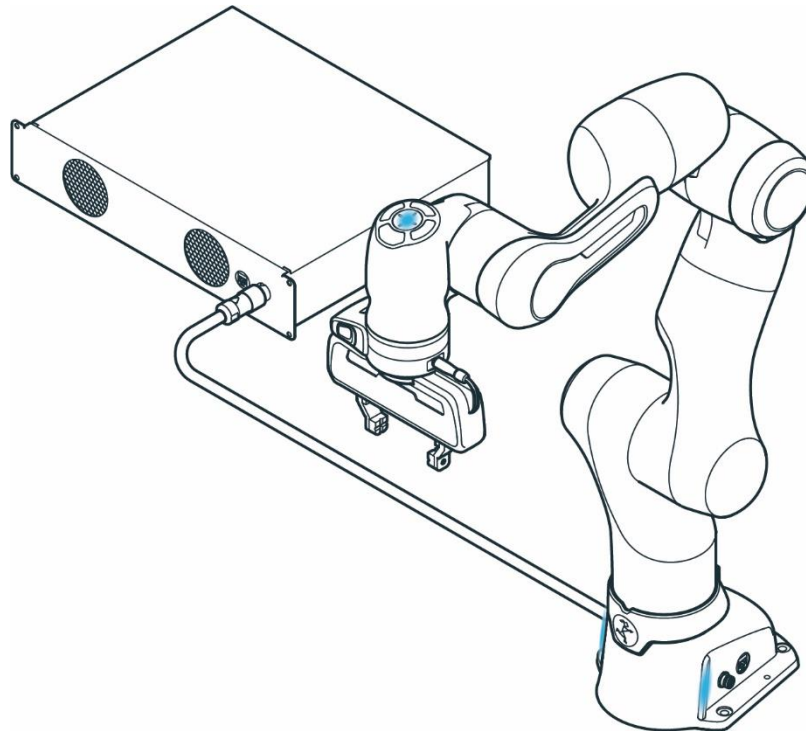


Fig. 8.2: Blue status lights on Arm

You can observe the following sequence:

- The cooling system will be activated and will be visible and audible.
- Booting up may take approximately 1 minute.
- The status lights on the Pilot and both sides of the base will start to flash.
- After the boot-up is complete, the status light will be continuously blue, indicating that the robot is in a stopped state in Execution mode.
If the status light flashes red, there exists a safety error. Then check whether the emergency stop is enabled or whether the Emergency Stop device has been correctly connected to input X3.

USAGE

- The fail-safe locking system is active. The joints are still mechanically locked. For information on unlocking the fail-safe locking system, see section Pre-guiding.
- The fail-safe locking system is active. The joints are still mechanically locked. For information on unlocking the fail-safe locking system, see section Pre-guiding in the event of a malfunction, ***please refer to Chapter 10 Maintenance and Disposal in this Manual.***

8.2 Safety Relevant Tests of Franka Research 3

8.2.1 Self-test of the robotic system

The self-tests of Control are executed while the system is running. The Arm is power cycled once to execute the self-tests of the Arm.

WARNING

Risk of injury due to falling objects

During the power cycle of the Arm, the power to the end effector is removed. Objects may fall from the end effector which could lead to injury.

- Remove all objects from the end effector.
- Move out of the hazard zone.

NOTICE

Every 24 hours, the user must initiate a safety diagnostic to detect potentially dangerous failures during operation. In the sidebar, the system will warn the user 2 hours before the time is exceeded.

If the time is exceeded, the robot will stop all operations and request the user to initiate the self-test. For this purpose, a message appears from which you can start the self-test.

You can also start the self-test manually at any time. To do this, proceed according to the following steps

1. Start Franka UI on your Interface device.
2. Go to "Settings".
3. Change to "Dashboard"
4. Click on Button "EXECUTE" next to the display of the countdown to the self-test

WARNING

Falling objects from end effectors due to a power supply cut off when SEEPO configuration is active

Risk of severe injury, such as crushing, tearing of the skin, and puncturing from the Arm and end effectors.

- Always wear personal protective equipment (e.g., safety shoes).
- Use the appropriate type of end effector to prevent objects from falling.
- Consider the shape, texture, and weight of the grasped objects in the risk assessment according to 10218-2. Using lightweight and/or round objects may reduce risks significantly.

8.2.2 Regular testing of safety functions

The function of some safety functions must be tested regularly. This applies to the following safety functions:

USAGE

Safety function	Test
Emergency Stop Device	<ol style="list-style-type: none">1. Press the Emergency Stop Device while the robot is not active.2. Check whether the brakes are locked.
Guiding Enabling Device	<ol style="list-style-type: none">1. Release the Pilot-Grip's Enabling Button while guiding. <i>The robot must stop.</i>2. Fully press the Pilot-Grip's Enabling Button while guiding. <i>The robot must stop.</i>
External Enabling Device	<ol style="list-style-type: none">1. Release the enabling button of the External Enabling Device while testing a task. <i>The robot must stop.</i>2. Fully press the enabling button of the External Enabling Device while testing a task. <i>The robot must stop.</i>
Any switch connected to X3.2 or X3.3	<ol style="list-style-type: none">1. Activate switch.2. Check if configured safety function is triggered accordingly.

NOTICE

- Activate the emergency stop system during start-up every 12 months.
- Reconnect the emergency stop system during start-up every 12 months.
- Check all safety installations, e.g., the emergency stop system, for function every 12 months.
- Check any additional safety measures that have been taken to secure safe operation.



For more information on the emergency stop system, ***please refer to Chapter 4.7 Installation of safety peripherals in this Manual.***

8.2.3 Testing the emergency stop

WARNING

Risk of severe injury due to a non-operating Emergency Stop Device

Using a non-operational Emergency Stop Device to stop an operation in case of an emergency may lead to staying trapped which leads to severe injuries, such as crushing, tearing of the skin, and puncturing from the Arm and end effectors.

- Store the Emergency Stop Device in a safe place.

NOTICE

Material damage

The end effector, workpieces, or the surroundings can be damaged when the device is stopped at an unfavorable point in the process.

- Only use the emergency stop in safety-critical situations.

USAGE

NOTICE

Any damage occurring to the Arm when the emergency stop is pressed will not cause harm to people, as the Arm will safely stop regardless of damage.

NOTICE

After an emergency stop, the Arm could have lost calibration or been damaged. If malfunctions are detected at the next start-up, the user will be informed.

NOTICE

Take other installed devices apart from Franka Research 3 into account that will be switched off by the emergency stop as well.

Precondition

- Franka Research 3 must be at standstill with no task running.
- The locking bolts of the fail-safe locking system must be open.
- The Arm may not move.

Procedure

1. Clear the space around the Arm to avoid any damage to grasped objects or surroundings.
2. Use the guiding to bring the Arm into a position that is free from obstacles, e.g., 200 mm above stationary objects.
3. Activate the emergency stop.

The Arm will slightly lower with a clicking noise when falling into the mechanical locking bolts.

NOTICE

For detailed instructions please refer to the Operating Manual, Chapter 2.1 Initial Connection in the respective Operating Manual corresponding to your system version (e.g., 5.6.0 or 5.8).

8.3 Switching Off and restarting

WARNING

Falling objects from end effectors due to a power supply cut off

Objects falling from the gripper lead to injuries to hands, fingers, feet, and toes.

- Always wear personal protective equipment (e.g., safety shoes).
- Use the appropriate type of grippers to prevent objects from falling.
- Consider the shape, texture, and weight of the grasped objects in the risk assessment according to 10218-2. Using lightweight and/or round objects may reduce risks significantly.

USAGE

Shutting down

NOTICE

The system is only completely shut down when the fans stop running.

Still running fans indicate that Franka Research 3 has not been completely shut down yet.

Repeat the Safety Instructions of Shutting down Franka Research 3.

SAFETY- INSTRUCTION

Procedure

1. Move out of the hazard zone.
2. In Franka UI, navigate in user menu to "Shutdown" and click on it.

The fail-safe locking system is activated.

Franka Research 3 will shut down.

Restarting

Turn off the system in Desk and wait for the fans to power down. Switch off the power switch on the Control. Wait one minute before Franka Research 3 again. To restart Franka Research 3, switch on the power switch on the back of the Control. Franka Research 3 will start to reboot again.

NOTICE

To avoid unintentional repowering of the system, secure the connecting cable in a safe place.

Disconnecting Franka Research 3 from the power supply

Procedure

1. Move out of the maximum workspace.
2. Go to Franka UI.
3. Select in User-Menu "Shut-down."

The system shuts down.

4. Switch off the power switch on the backside of the Control.
5. Pull the cable from the backside of the Control.

NOTICE

To avoid unintentional restarting of the system, stow the power cord in a safe place.

Franka Research 3 is disconnected from the power supply.

9 ROBOT LED INDICATOR SYSTEM

9.1 Overview of the status indicators

Status lights on both sides of the base take on the corresponding color, like a traffic light. The status lights will flash slowly during boot-up, when Franka Research 3 requires attention, or when the user enters values. The lights are lit in the corresponding color of the Franka Research 3 status during other processes. A circular status light in the middle of the Pilot-Disc indicates Franka Research 3's status as well.

When the operator is interacting with the Arm, the status light on the Pilot-Disc is switched off.

For more information on the color behavior, please refer to Chapter 8.1 Switching On in this Manual.

WARNING

Dangerous and uncontrolled motions of the Arm

Risk of severe injury, such as crushing, tearing of the skin, and puncturing from the Arm and end effectors.

- Ensure that the end effector and/or the object mass and the Center of Mass (CoM) are parameterized correctly.
- Keep out of the workspace during operation.

NOTICE

The motion speed in Teach or Hand-Guiding Mode is pre-set. The speed can be reduced according to the risk evaluation of the Arm within its application.

The robot uses six distinct LED colors to communicate its operational status. These lights can appear in two patterns:

- **Solid (Static)** – indicating a stable state
- **Blinking (Flashing)** – indicating a transitional or attention-required state

LED indicators are visible in three key locations:

- Desk Interface
- Base of the Robot
- Pilot-Disc

Each color and pattern combination provides critical information about the robot's current state and whether it is safe to interact with. These visual cues are designed to enhance operator awareness and safety.

9.2 LED Activation Behavior

- **Base LEDs:**
Always active. They continuously reflect the robot's operational status and are the primary reference for system state.
- **Pilot LEDs:**
Only active during programming or manual guidance. They provide feedback specific to user interaction during these modes.

Startup Checklist for Operators

When powering on the robot, always:

1. Verify that all indicator lights are functioning properly.
2. Ensure the LEDs are clearly visible from your working position.
3. Interpret the lights correctly to determine whether the robot is:
 - IDLE or in TEACH mode
 - Executing a Task
 - In an error or warning state
 - Awaiting user input

Important Safety Considerations

Whether it is safe to approach the robot in each state depends on the application-specific risk and hazard analysis and the configured safety scenarios. While LED indicators help identify the robot's state, they are not certified safety functions.

If in doubt, always use appropriate safety measures before approaching the robot:

- Press the Emergency Stop
- Activate a Protective Stop
- Confirm the robot is in a safe state via the user interface

9.3 Flashing Patterns

Pattern	Frequency	Meaning
Slow Flashing	~0.6 Hz (approximately 2 flashes every 3 seconds)	Indicates a transition between states or a request for user attention
Fast Flashing	~2 Hz (approximately 2 flashes per second)	Warns that motion is initiating, the robot is moving slowly, or the system is updating

Error and Communication Loss

If a visual indicator or its controlling device detects a loss of communication, this is signaled as an error with a solid red light.










9.4 LED Priority Logic

- The LED system always displays the most critical state.
- If multiple events occur simultaneously, the color with the highest significance is shown.
- Within the same priority level, only a one-color scheme is displayed at a time to avoid confusion.


9.5 LED Color Reference Table

Category	LED Color	LED Pattern	Status Meaning	User Action
Robot System Status	White	Static	System idle or in TEACH mode.	✅ Observe standard safety precautions before approaching.
	White	Slow Flashing	Booting or shutting down.	🚫 Do not interrupt. Wait until complete.
	White	Fast Flashing	System updating.	🚫 Do not unplug or interrupt. Wait until complete.
Brakes	Yellow	Static	Brakes locked/unlocked.	⚠️ Used during brake operations.
	Yellow	Slow Flashing	Waiting for the boot to finish.	⌚ Wait until solid yellow or next instruction.
Warnings	Yellow	Static	Warning state.	🚫 Do not approach. Check UI.
	Yellow	Slow Flashing	Warning: user interaction needed.	⚠️ Check UI and acknowledge warning.
Safety Errors	Red	Static	Severe errors (e.g., safety, system, communication).	🚫 Do not approach. Investigate via UI.
	Red	Slow Flashing	Safety violation or application error.	⚠️ Check UI. Approach only if safe and trained.

ROBOT LED INDICATOR SYSTEM

Category	LED Color	LED Pattern	Status Meaning	User Action
	Red	Fast Flashing	Error recovery in progress.	 Wait or reset via UI.
	Red	Slow Flashing	Input needed to recover from error. Error recovery possible with user input (e.g., joint limits exceeded during hand-guiding)	 Unlock joint or reset to resume operation.
Execution	Green	Static	Task executing autonomously.	 Do not approach. Robot is moving.
	Green	Fast Flashing	Execution will begin shortly (e.g., FCI countdown).	 Do not approach. Execution imminent.
	Green	Slow Flashing	Task active in assist mode.	 Approach with care. Follow safety protocols.
Execution Mode	Blue	Static	Execution phase ready. Brakes engaged.	 Approach with care. Robots may start moving.
	Blue	Slow Flashing	Brakes opening or collaborative mode (no active task).	 Robot may move slightly. Safe to approach.
	Blue	Slow Flashing	Execution interrupted. Awaiting feedback.	 Provide feedback to resume. Robot paused.
	Magenta	Static	Conflicting input detected (e.g., manual guidance vs. automation).	 Do not approach. Resolve input conflict.

MAINTENANCE AND DISPOSAL

Category	LED Color	LED Pattern	Status Meaning	User Action
Conflicts	Magenta	Slow Flashing	Input needed to resolve conflict.	 Provide input or guidance to continue.

10 MAINTENANCE AND DISPOSAL

10.1 Maintenance

The Arm and Control of the Franka Research 3 are designed for maintenance-free operation of approx. 20,000 hours under normal operating conditions. The normal operating conditions used as a basis were derived from various representative applications of the robot system and verified by analyses and tests. If a customer application deviates greatly from these underlying operating conditions, the duration may be extended or shortened under certain circumstances.

If the system approaches this operating duration, contacting Franka Robotics support (support@franke.de) is recommended. Evaluation of the system's log data by the support team will then indicate any necessary action.

NOTICE

If damage to the structure of the robot is detected during a visual inspection, the robot must be taken out of service regardless of the current operating time.

NOTICE

- Activate the emergency stop system during start-up every 12 months.
- Reconnect the emergency stop system during start-up every 12 months.
- Check all safety installations, e.g., the emergency stop system, for function every 12 months.
- Check any additional safety measures that have been taken to secure safe operation.

10.2 Cleaning

DANGER

Risk of electric shock

Improper use of liquid cleaning agents, as well as incorrectly disconnected devices from the power supply, can lead to fatal accidents.

- Do not clean devices that have not been safely disconnected from the power supply.
- Do not use liquid cleaning agents for cleaning the devices.

The following things need to be kept in mind while cleaning:

- Only qualified people may carry out cleaning.

MAINTENANCE AND DISPOSAL

- Cleaning is only permissible when Franka Research 3 is safely stopped and disconnected from the power supply.
- Switching off and disconnecting the device must be carried out by qualified people only.
- Do not use any liquids to clean the device.
- Do not use any cleaning chemicals.
- The components may only be cleaned with dry cloth.
- Moisture must not enter the device.
- Do not apply great force to the Arm. The parts to be cleaned must be supported manually to not overload and possibly damage the Arm.

NOTICE

Material damage to the devices

- Do not use liquid cleaning agents for cleaning the devices.

10.3 Disposal

Disposal

Disposal of Franka Research 3 must comply with the relevant country-specific laws, standards, and regulations.

Battery

The Control contains a coin cell battery. The cell battery must be disposed of separately according to the relevant country-specific laws, standards, and regulations.

To remove the battery, open the Control.

NOTICE

Opening the Control is only allowed for the purpose of removing the coin cell battery when disposing of it.

Return of waste of packaging

Please contact Franka Robotics to return any used packaging.

10.4 Mechanical exchange of Control

To replace the Control mechanically, proceed according to the following steps:

- Perform a backup, if the defective Control still allows it.
- Shut down the robot.
- Switch off the Control at the main switch
- Disconnect the Control from the mains supply
- Disconnect the plug connection C1 (connection cable to the robot) as well as the network connection C2 at the Control.
- Replace the defective Control with the replacement Control. Observe the required environmental conditions for the Control (see Chapter 10 "Assembly & Installation").
- Connect the robot's connection cable to the C1 plug connection on the Control
- Establish the network connection C2
- Connect the Control to the mains supply

TROUBLESHOOTING

- Connect your operating device to connector X5 on the robot base using an Ethernet cable
- Switch on the Control with the main switch.

11 TROUBLESHOOTING

Troubleshooting and Safety Error Reference

For a list of recoverable safety errors, *please refer to Chapter 4.10 – Safety Functions in this Manual.*

For troubleshooting the following issues, *please refer to Chapter Troubleshooting in the respective Operating Manual corresponding to your system version (e.g., 5.6.0 or 5.8)."*

- Loud clicking sound during shutdown
- Joint limit error
- Joint position error
- Failed to unlock joints
- Robot does not complete booting
- Desk continuously displays "Shutting down the system"
- Robot does not boot after turning on the Control
- Unexpected stop followed by restart

If recovery is unsuccessful, the issue may be due to hardware damage or a software malfunction. In such cases:

1. Shut down the system immediately.
2. Take the robot out of operation.
3. Contact your service partner or Franka customer support [Contact: support@franka.de](mailto:support@franka.de)

NOTICE

When contacting support, please provide the serial number of the Arm and the robot log files. You can download the log files via: Desk For detailed instructions, *see Chapter 13.2 – Log Download in the Operating Manual.*

12 TECHNICAL DATA

For the latest datasheet on Franka Research 3 (Document Number: **R02212**), please visit:

- Link <https://franka.de/documents>
- You will find the datasheet titled:
"FR3 Arm v1" available in both English and German.

NOTICE

The probability of failure per hour was evaluated at 40 °C. However, the safety rating holds for all functions within the temperature range, including the extended temperature range.

If the system probability of failure per hour values are used for calculation, take the temperature into account.

For more information on the probability of failure per hour, contact Franka Robotics at support@franka.de.

12.1 Ambient conditions for delivery and transport

The system can be stored and transported within the temperature range of -25 °C to +70 °C.

13 TRANSPORTATION AND HANDLING

WARNING

Heavy equipment

Due to the dead weight and partly due to the geometric design, lifting and handling the equipment may cause back injury and, if it falls, serious injury to fingers, hands, toes, and feet.

- Always wear personal protective equipment (e.g., safety shoes), when transporting, mounting, or demounting the equipment.
- Always lift the equipment with the help of a second person.
- The equipment must be placed on even surfaces to prevent it from tilting or sliding.
- Follow the existing company regulations on lifting loads and personal protective equipment.

NOTICE

Material damage to Arm, end effectors, and objects in the maximum workspace

Sensitive electromechanical components in the Arm and end effectors may be damaged if end effectors are connected to the Arm while bringing it into transport position.

- Dismount any end effector and attachments before bringing the Arm into transport position.

Do not leave loose objects in the maximum workspace.

DANGER

Material damage to Arm and Control

Mechanical shock may cause damage or loss of calibration of sensitive electromechanical components in the Arm and Control. Avoid shock.

- Do not set down the devices roughly.
- Always store and transport the devices in their original packing, even inside buildings.

13.1 Procedure

To ensure the safe and secure movement of the Franka Research 3 robot, users must follow a three-step procedure. This process is critical to protect the robot from damage during handling, transport, or storage.

Precondition

- The end effector and attachments must be removed from the Arm.
- The robot needs to move freely without being encumbered by obstacles to adopt the transport pose. In case of obstacles in the robot cell, consider moving the robot closer to the transport pose via hand-guiding.

TRANSPORTATION AND HANDLING

The procedure consists of the following steps:

13.1.1 Step 1. Transportation pose of the Arm

Before handling or transporting the robot, it must first be moved into a safe and compact configuration called the transport mode.

To do this, use the "Move to pack pose" function found in the Settings menu of the Franka UI. This ensures that all joints are correctly positioned to minimize the risk of damage during lifting or packaging.

For detailed instructions with images on how to prepare the robot for transport mode using the "Move to pack pose" function in Programming and Execution mode Settings in the Franka UI, *please refer to Chapter 13.5 Transportation pose of the Arm in the respective Operating Manual corresponding to your system version (e.g., 5.6.0 or 5.8).*

After that, follow the handling and lifting and re-packing the Arm instructions.

13.1.2 Step 2. Handling and lifting

Always lift the Arm in the positions intended for lifting (see graphic below) to not overstress the Arm joints during handling and lifting. In particular, the Arm may never be carried in the extended position with one person holding each end of the Arm.

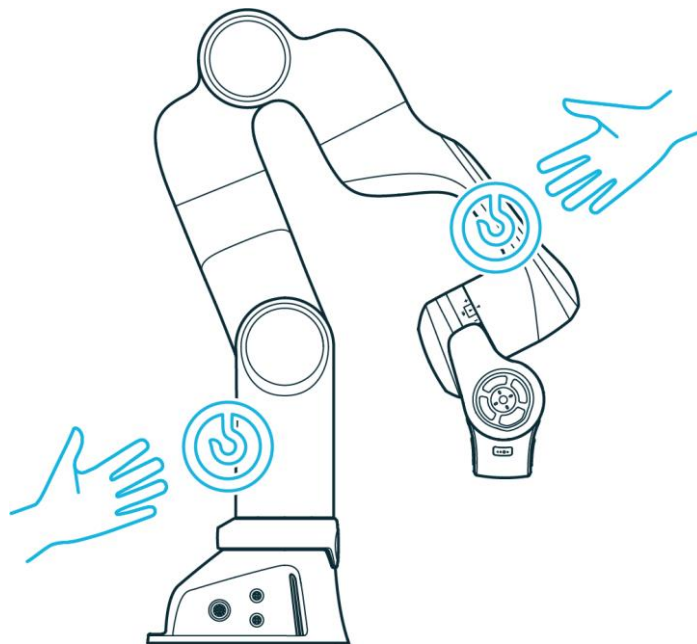


Figure 3: Lifting positions

NOTICE

Material damage to the Arm

Moving the Arm by force in a locked state will lead to a momentary slipping of internal parts which causes loss of calibration and damage to the Arm.

- Handle, lift, and transport the Arm only at the points indicated in this manual to avoid overstressing the joints of the Arm.
- The Arm is to be handled gently even when set-up and switched on or off.

TRANSPORTATION AND HANDLING

NOTICE

Do not step or lean on the Arm.

NOTICE

Pay attention to the weight.

The weight of the Arm is approx. 17,8 kg

WARNING

Heavy equipment

Due to the dead weight and geometric design, lifting and handling the equipment may cause back injury, and, if it falls, serious injury to fingers, hands, toes, and feet.

- Always wear personal protective equipment (e.g., safety shoes) when transporting, mounting, or demounting the equipment.
- The equipment must be placed on even surfaces to prevent it from tilting or sliding.
- Follow the company regulations on lifting loads and personal protective equipment.

WARNING

Damaged components

Electrical risks may lead to severe injury.

- Check if the packaging is in sound condition and fulfils its protective function.
- Check the cables, plugs, and mechanical housing for cracks and broken insulations.
- Do not use damaged cables, plugs, and mechanical housing for operation. In cases of doubt, contact Franka Robotics.

NOTICE

Material damage to Arm and Control

Mechanical shock may cause damage or loss of calibration of sensitive electromechanical components in the Arm and the Control.

- Avoid shock.
- Carefully set down the devices.
- Always store and transport the devices in their original packing, even for short-distance transport.

NOTICE

Material damage to Arm, end effectors, and objects in the maximum workspace

Sensitive electromechanical components in the Arm and end effectors may be damaged if end effectors are connected to the Arm while bringing it into transport position.

- Dismount any end effector before bringing the Arm into transport position.
- Do not leave loose objects in the maximum workspace.

13.1.3 Step 3. Re-packing the Arm

Precondition

- The robot needs to be in the transport pose.

Procedure

1. Open the box.
2. Grasp the Arm at the indicated lifting positions in twos and carefully put it inside the bottom protective layer.

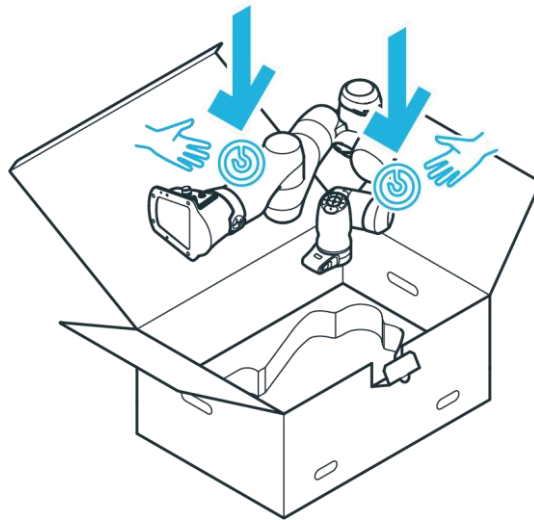


Figure 4: Lifting the Arm

3. Insert the middle protective layer.

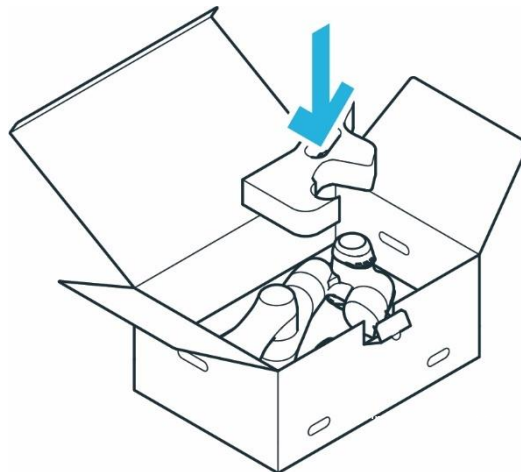


Figure 5: Packing the Arm

TRANSPORTATION AND HANDLING

4. Insert the top protective layer.

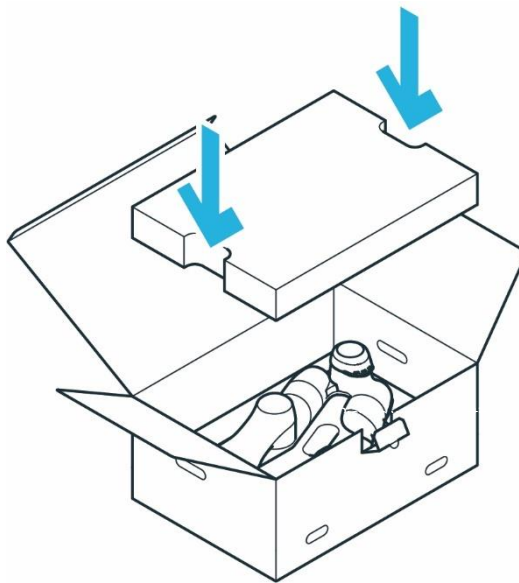


Figure 6: Closing the box

5. Close the foil coating.
6. Close the box.

14 APPENDIX

14.1 Stopping Times and Distances

The following diagram shows the stopping times and stopping distances of axes 1 to 4 according to the requirement from EN ISO 10218-1.

The following table illustrates the joint positions for the different extension states of the Franka Research 3.

	Extension 100%	Extension 66%	Extension 33%
Joint 1			
Joint 2			
Joint 3			
Joint 4			

Fig. 14.1: Illustration of extension states

14.2 Stop Category 1

The following table summarizes the maximum values of axes 1-4 for stop category 1.

Joint	max. stopping distance [rad]	max. stopping time [s]
1	0,803	0,735
2	0,798	0,682
3	0,684	0,534
4	0,455	0,418

14.2.1 Joint 1

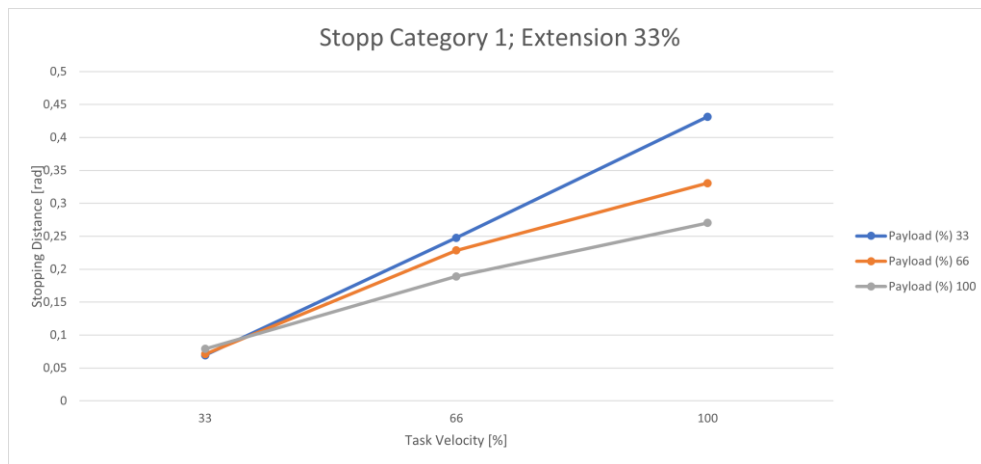


Fig. 14.2: Distance; Stopp Cat 1; Joint 1, 33%

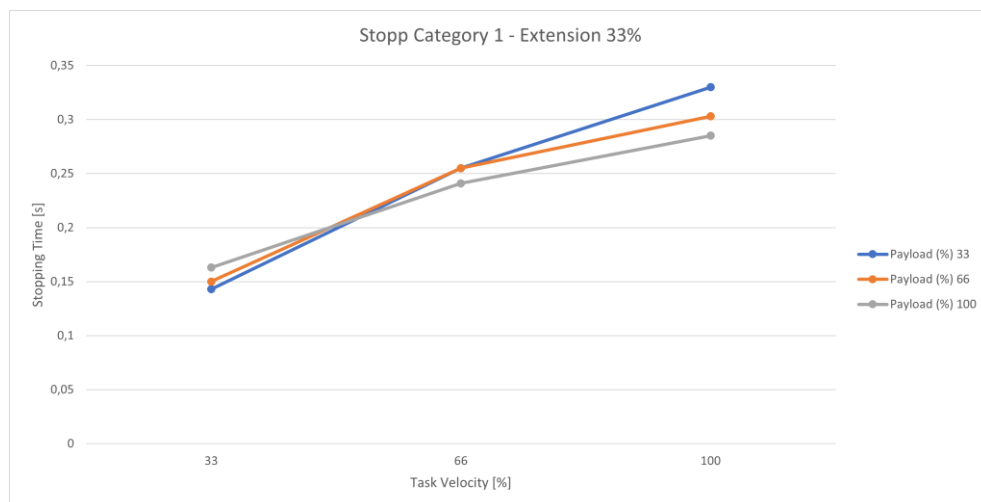


Fig. 14.3: Time; Stopp Cat 1; Joint 1 33%

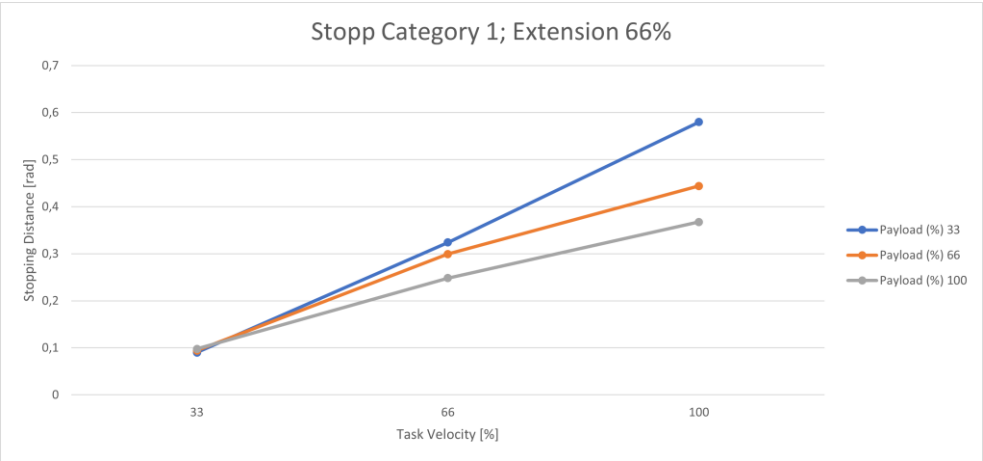


Fig. 14.4: Distance; Stopp Cat 1; Joint 1 66%

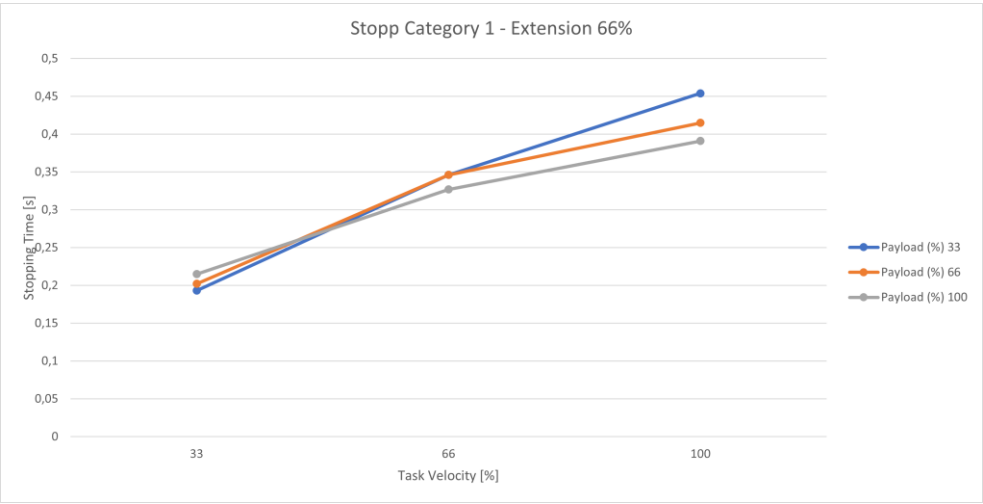


Fig. 14.5: Time; Stopp Cat 1; Joint 1 66%

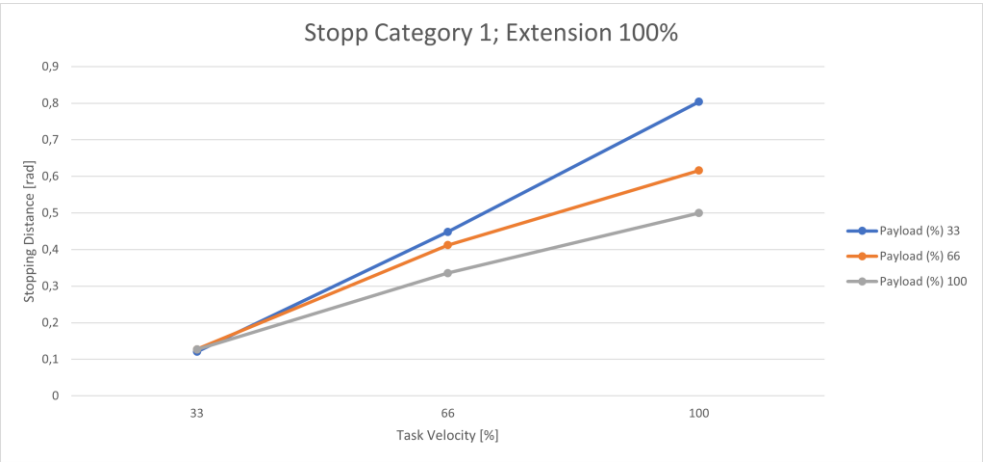


Fig. 14.6: Distance; Stopp Cat 1; Joint 1 100%

APPENDIX

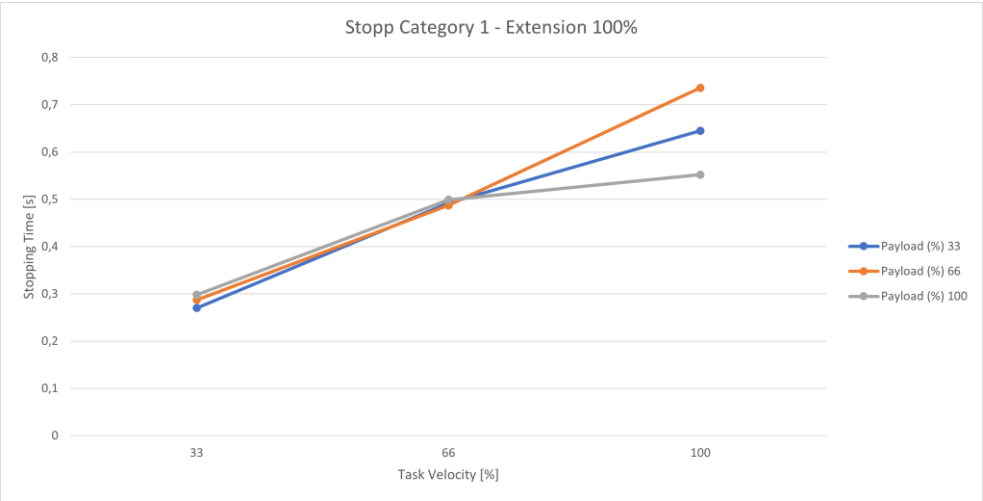


Fig. 14.7: Time; Stopp Cat 1; Joint 1 100%

14.2.2 Joint 2

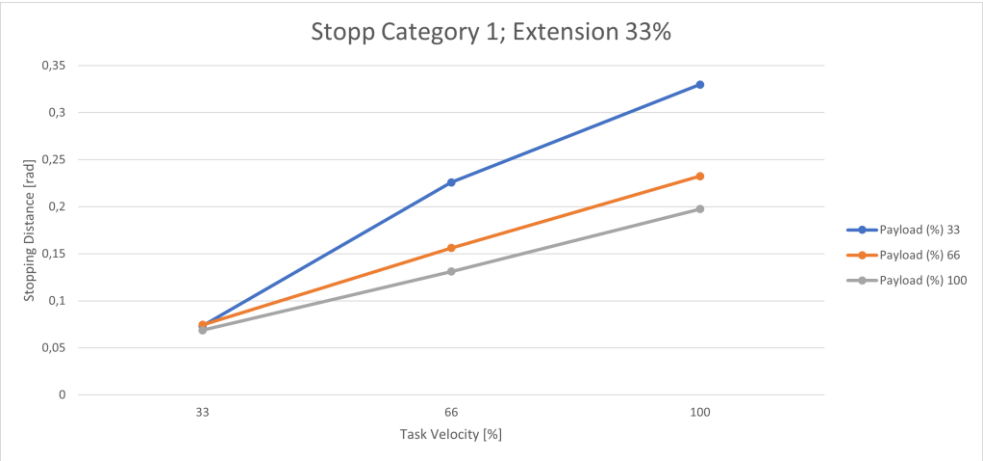


Fig. 14.8: Distance; Stopp Cat 1; Joint 2 33%

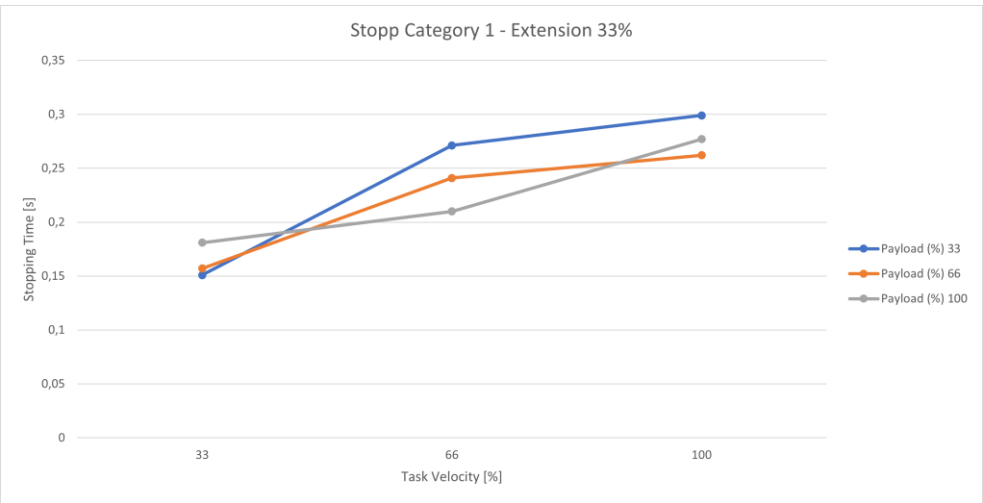


Fig. 14.9: Time; Stopp Cat 1; Joint 2 33%

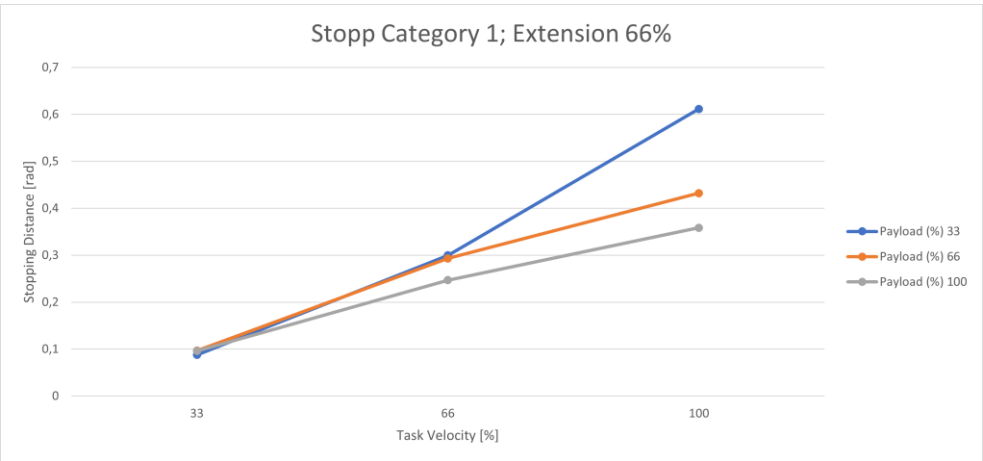


Fig. 14.10: Distance; Stopp Cat 1; Joint 2 66%

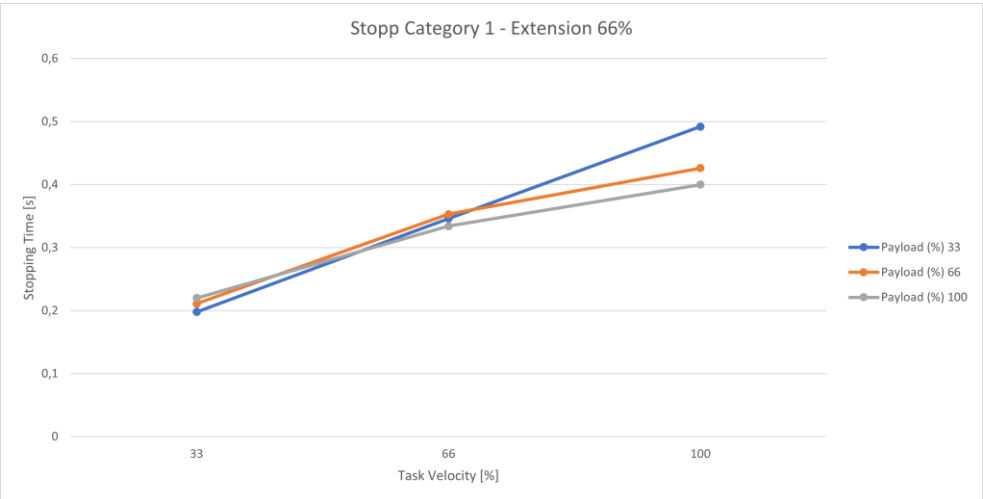


Fig. 14.11: Time; Stopp Cat 1; Joint 2 66%

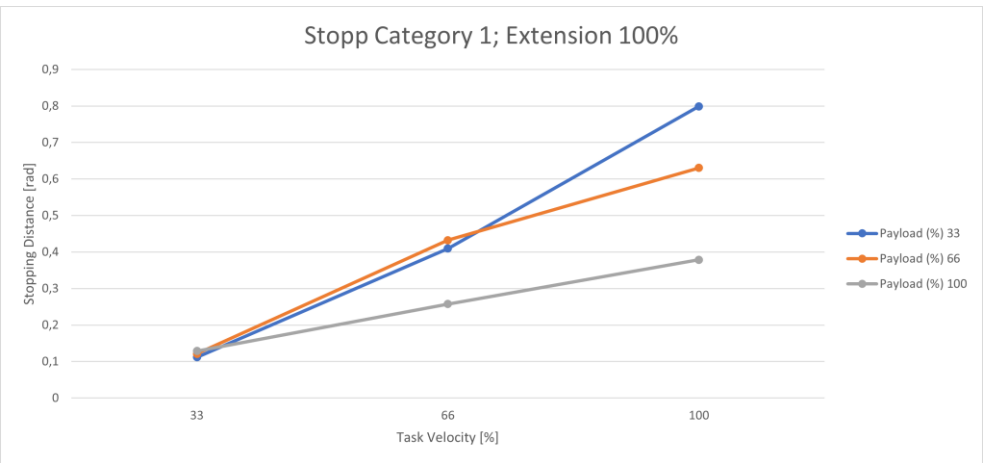


Fig. 14.12: Distance; Stopp Cat 1; Joint 2 100%

APPENDIX



Fig. 14.13: Time; Stopp Cat 1; Joint 2 100%

14.2.3 Joint 3

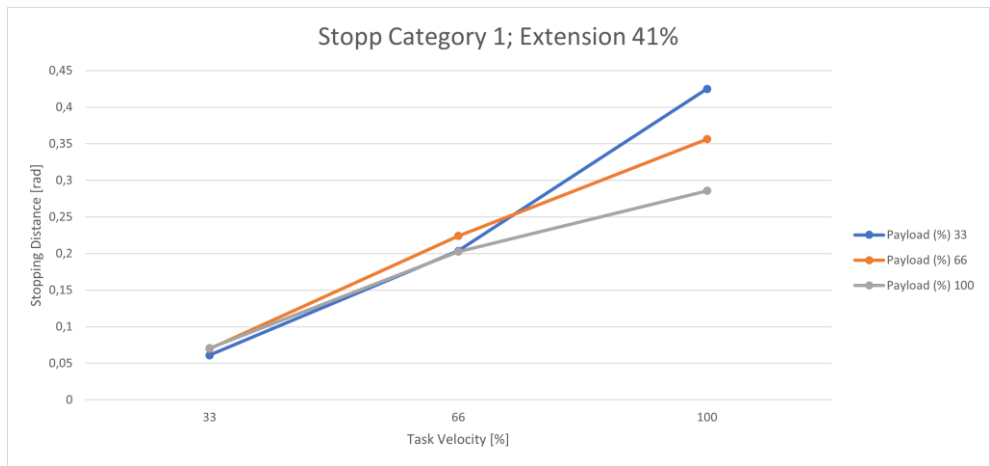


Fig. 14.14: Distance; Stopp Cat 1; Joint 3 33%

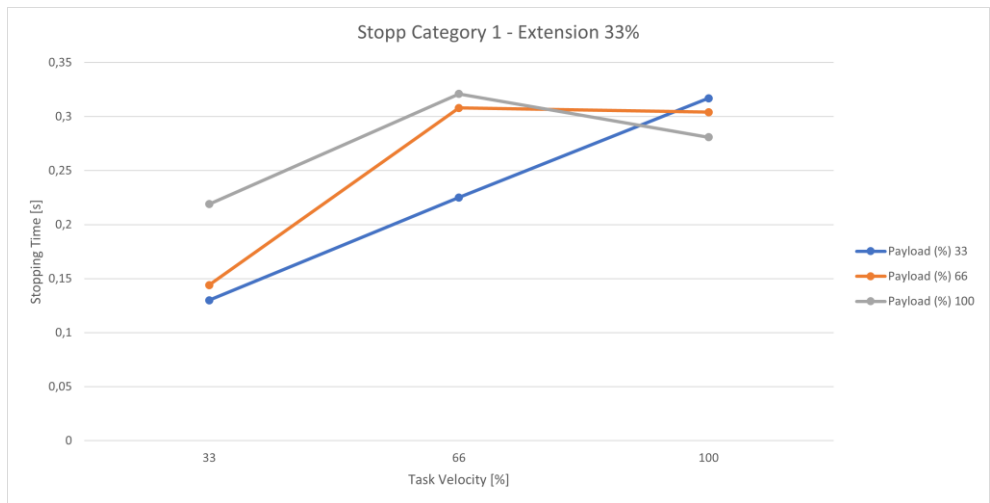


Fig. 14.15: Time; Stopp Cat 1; Joint 3 33%

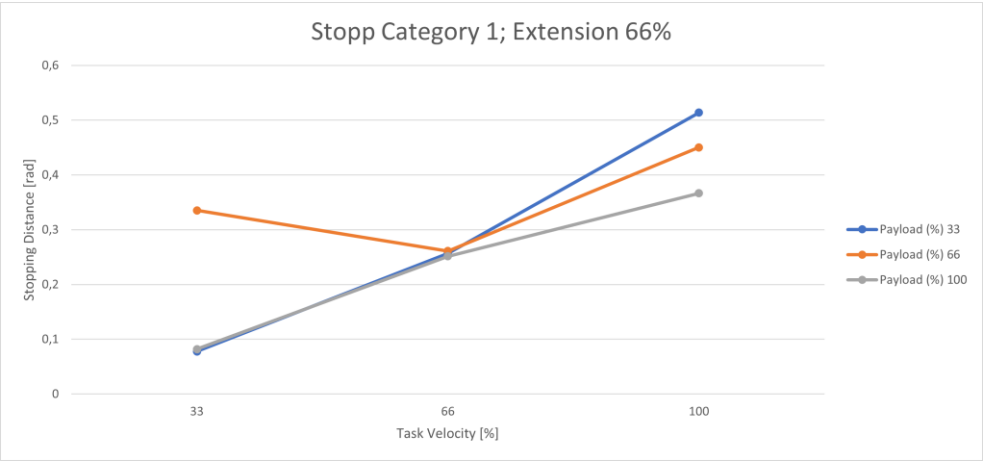


Fig. 14.16: Distance; Stopp Cat 1; Joint 3 66%

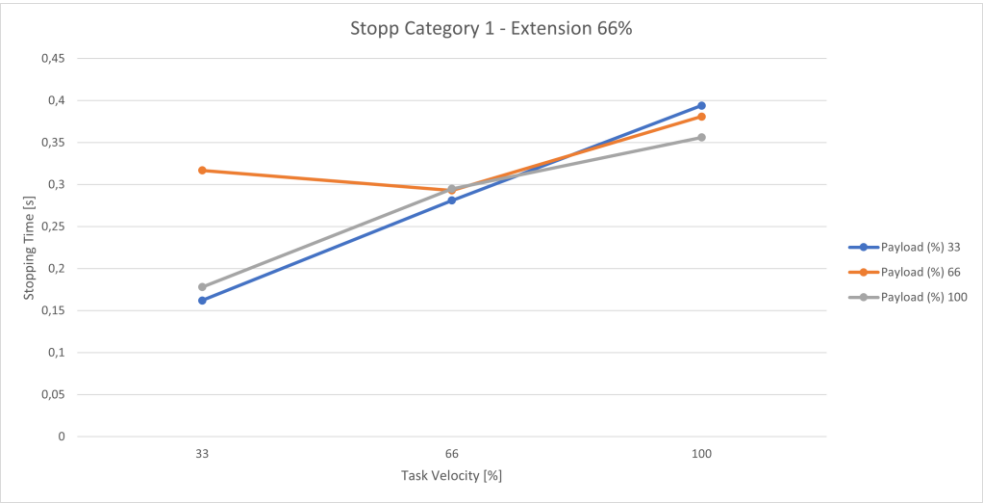


Fig. 14.17: Time; Stopp Cat 1; Joint 3 66%

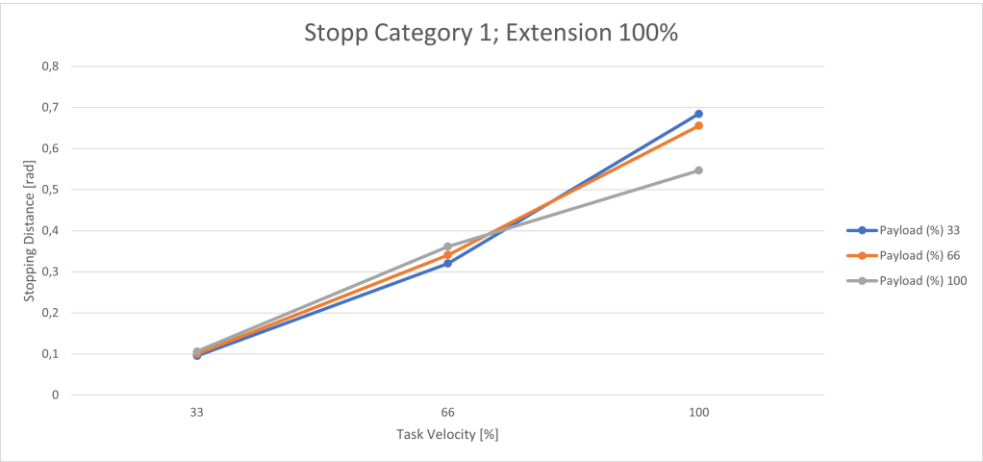


Fig. 14.18: Distance; Stopp Cat 1; Joint 3 100%

APPENDIX

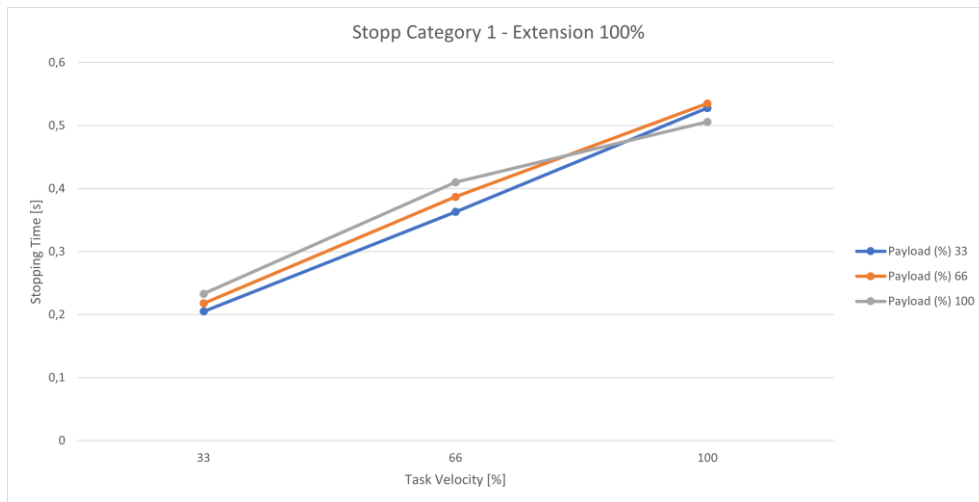


Fig. 14.19: Time; Stopp Cat 1; Joint 3 100%

14.2.4 Joint 4

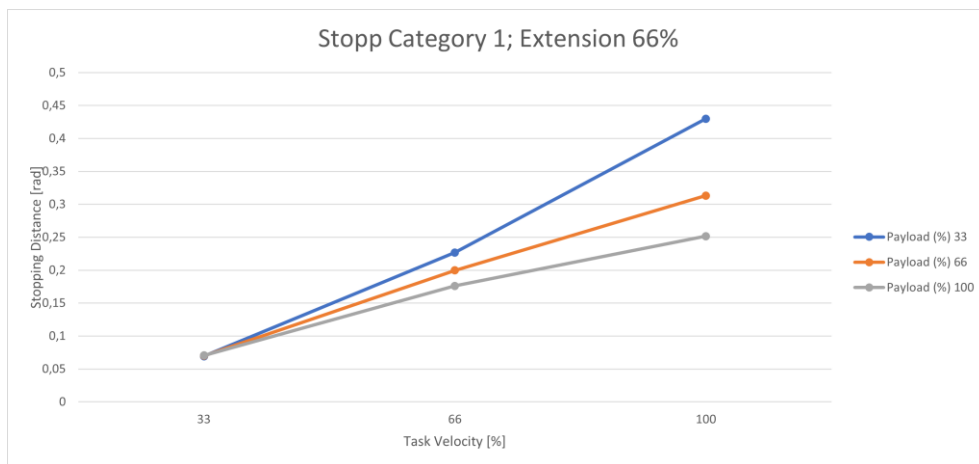


Fig. 14.20: Distance; Stopp Cat 1; Joint 4 66%

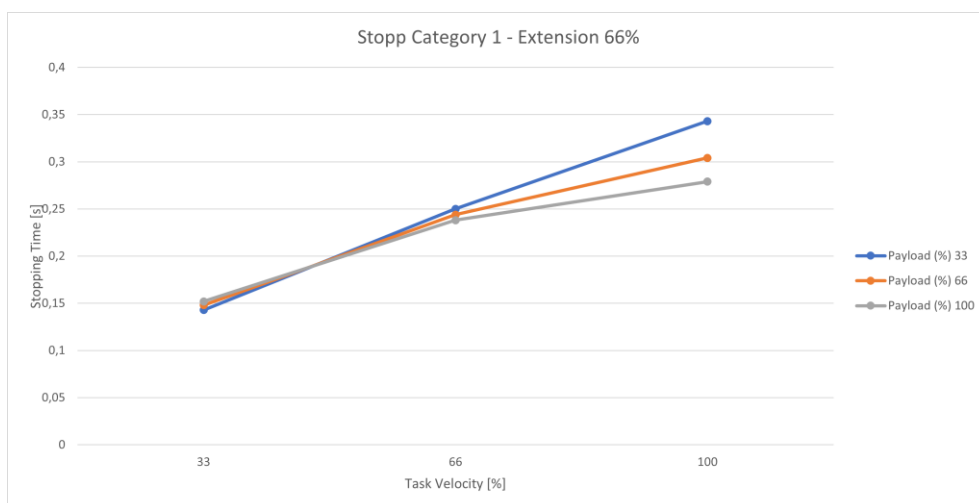


Fig. 14.21: Time; Stopp Cat 1; Joint 4 66%

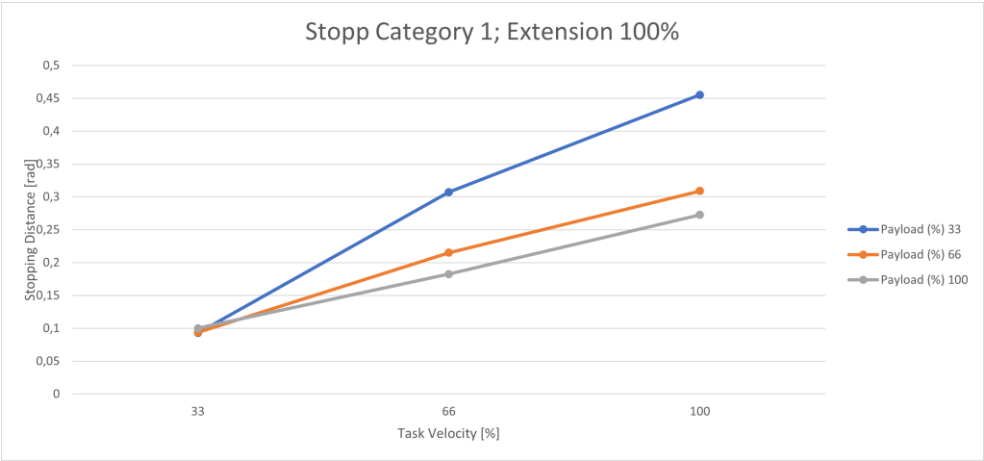


Fig. 14.22: Distance; Stopp Cat 1; Joint 4 100%



Fig. 14.23: Time; Stopp Cat 1; Joint 4 100%

14.3 Stop Category 2

The following table summarizes the maximum values of axes 1-4 for stop category 2.

Joint	max. stopping distance [rad]	max. stopping time [s]
1	0,806	0,743
2	0,805	0,684
3	0,695	0,542
4	0,455	0,417

14.3.1 Joint 1

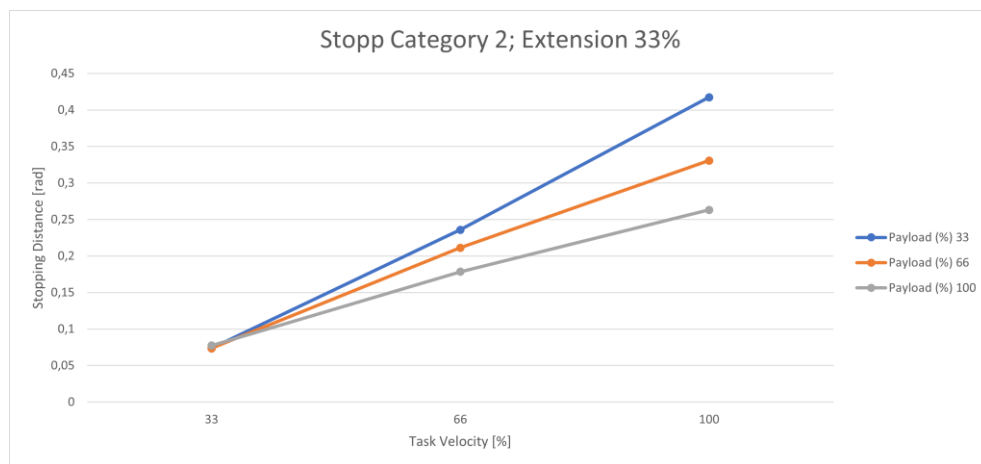


Fig. 14.24: Distance; Stopp Cat 2; Joint 1 33%

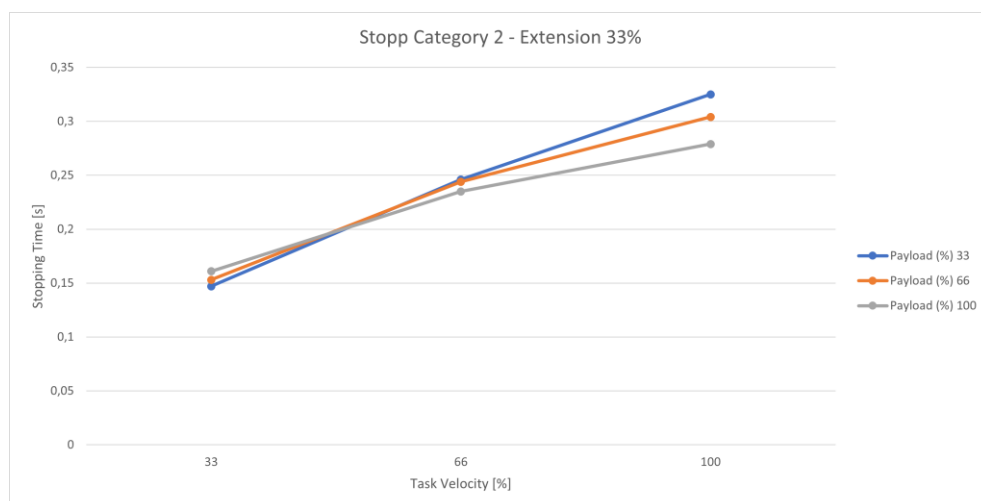


Fig. 14.25: Time; Stopp Cat 2; Joint 1 33%

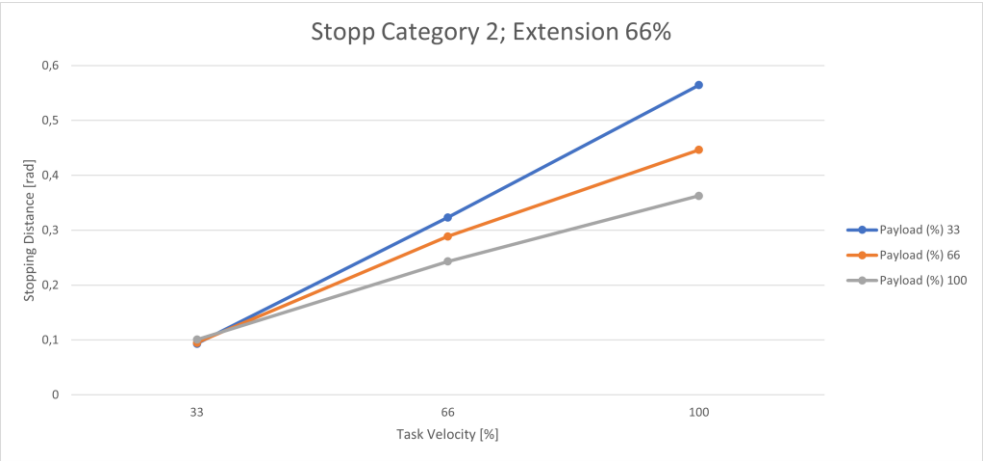


Fig. 14.26: Distance; Stopp Cat 2; Joint 1 66%

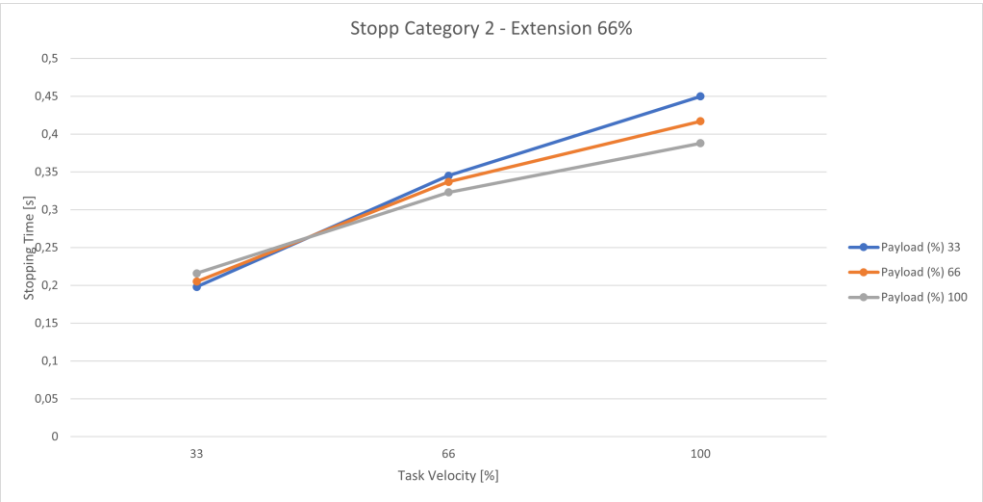


Fig. 14.27: Time; Stopp Cat 2; Joint 1 66%



Fig. 14.28: Distance; Stopp Cat 2; Joint 1 100%

APPENDIX

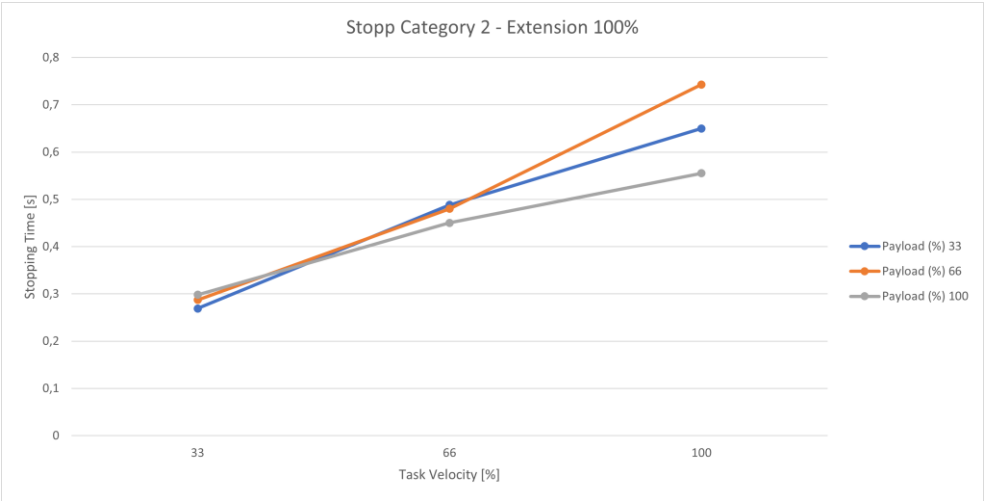


Fig. 14.29: Time; Stopp Cat 2; Joint 1 100%

14.3.2 Joint 2



Fig. 14.30: Distance; Stopp Cat 2; Joint 2 33%

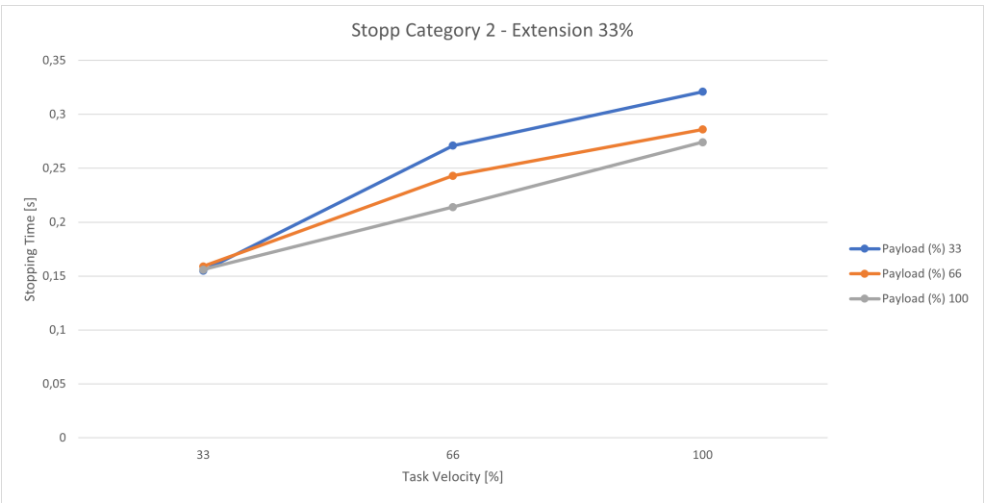


Fig. 14.31: Time; Stopp Cat 2; Joint 2 33%

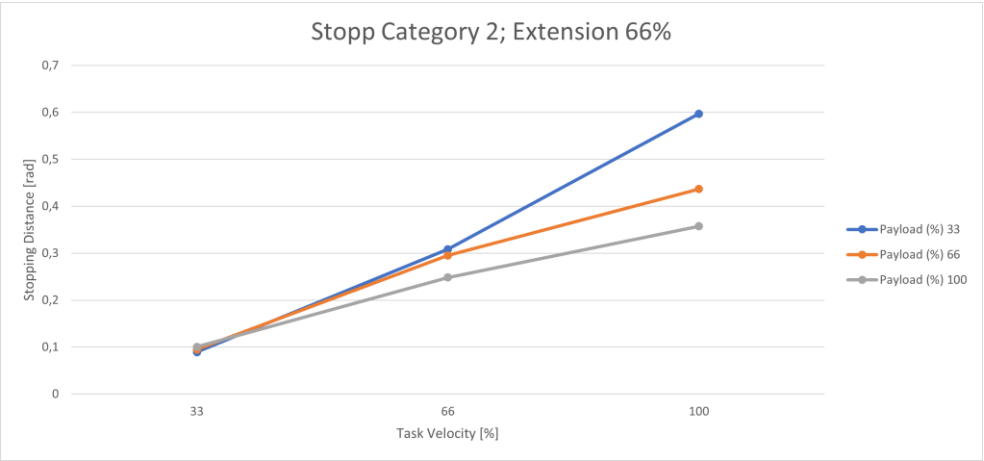


Fig. 14.32: Distance; Stopp Cat 2; Joint 2 66%

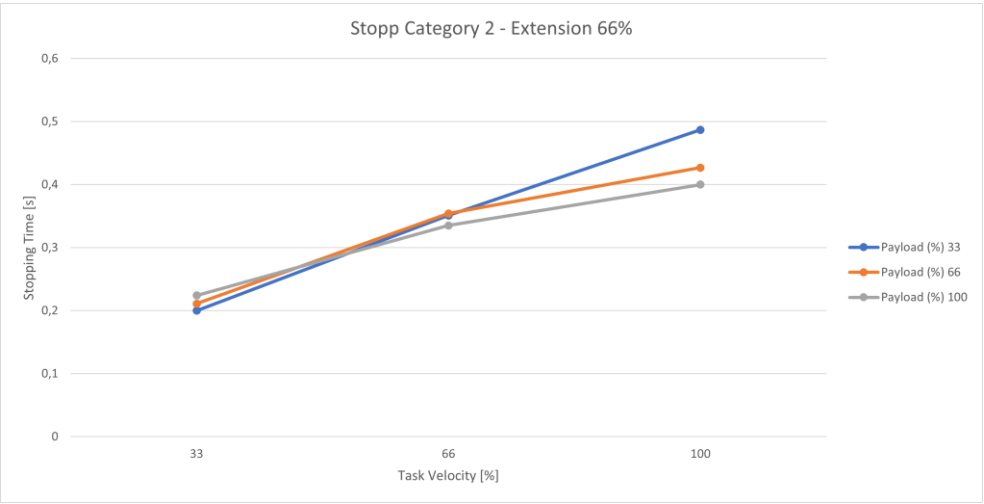


Fig. 14.33: Time; Stopp Cat 2; Joint 2 66%



Fig. 14.34: Distance; Stopp Cat 2; Joint 2 100%

APPENDIX



Fig. 14.35: Time; Stopp Cat 2; Joint 2 100%

14.3.3 Joint 3

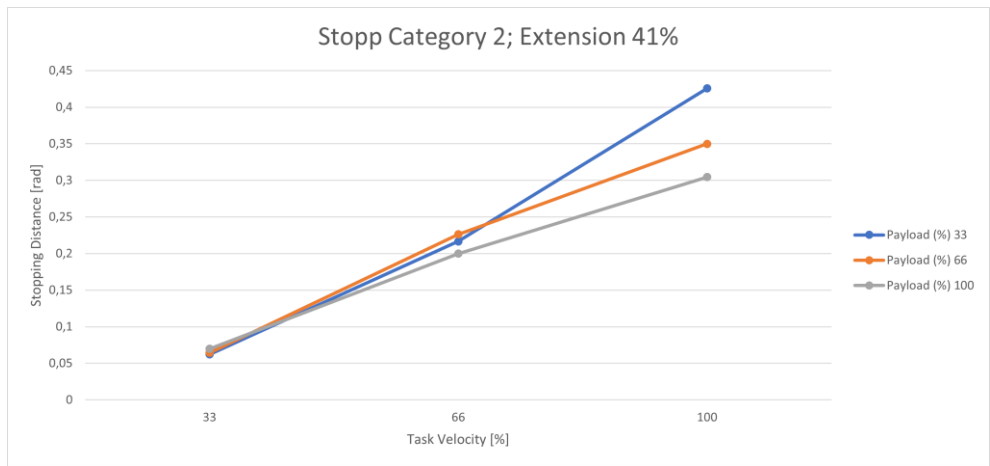


Fig. 14.36: Distance; Stopp Cat 2; Joint 3 41%

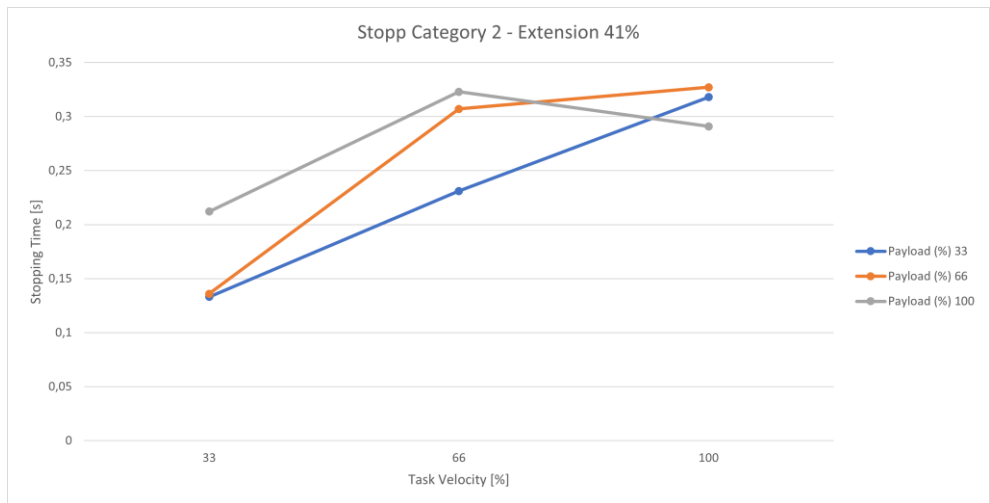


Fig. 14.37: Time; Stopp Cat 2; Joint 3 41%

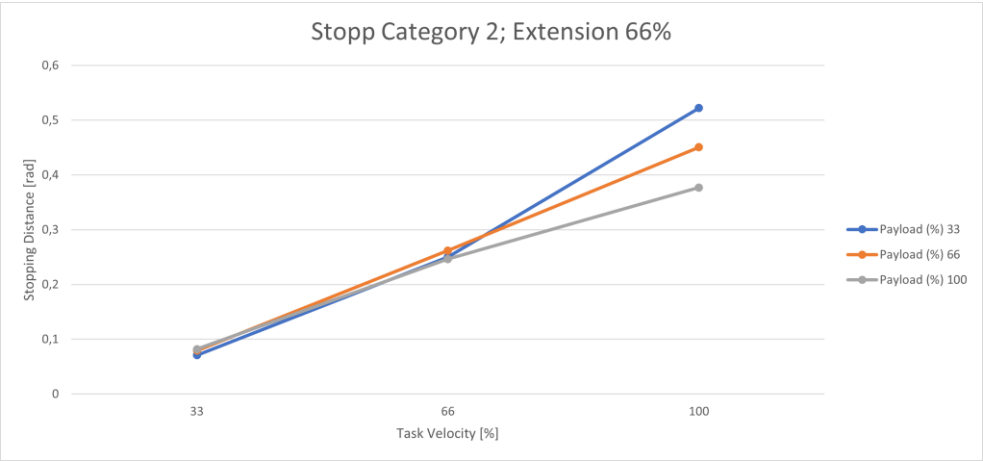


Fig. 14.38: Distance; Stopp Cat 2; Joint 3 66%

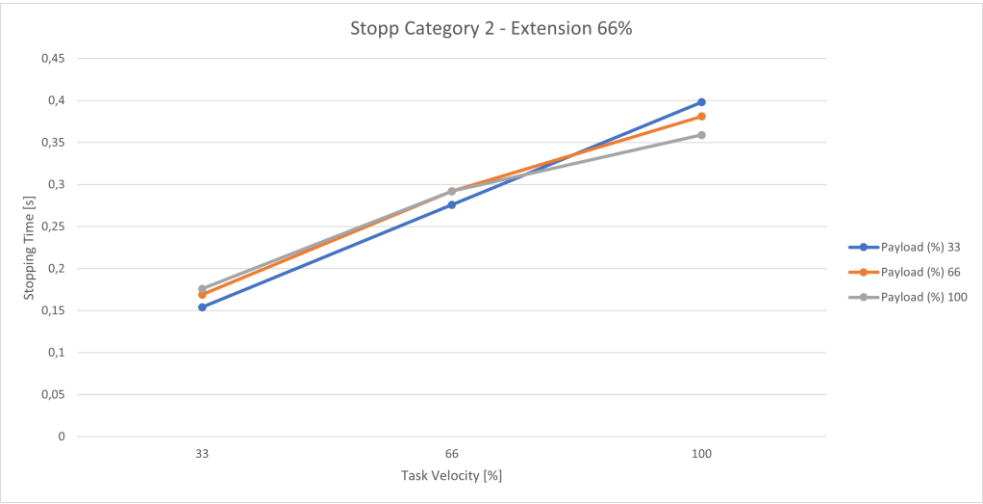


Fig. 14.39: Time; Stopp Cat 2; Joint 3 66%

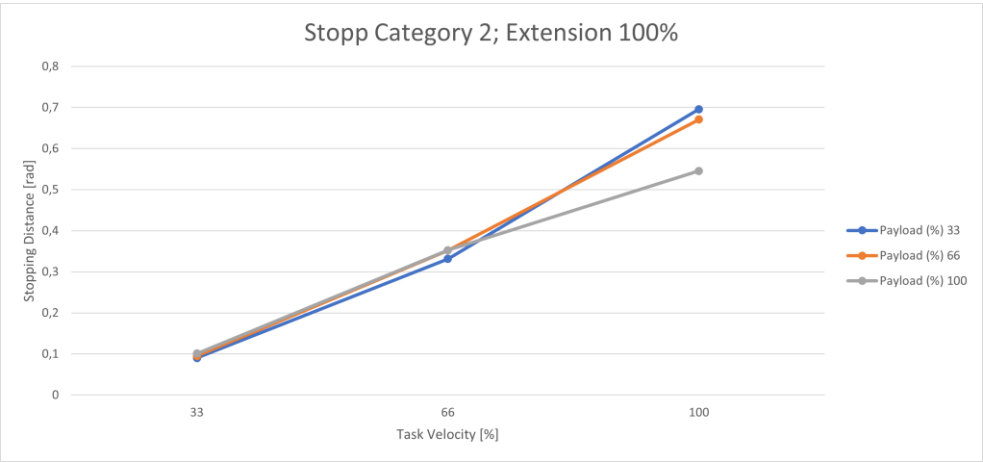


Fig. 14.40: Distance; Stopp Cat 2; Joint 3 100%

APPENDIX

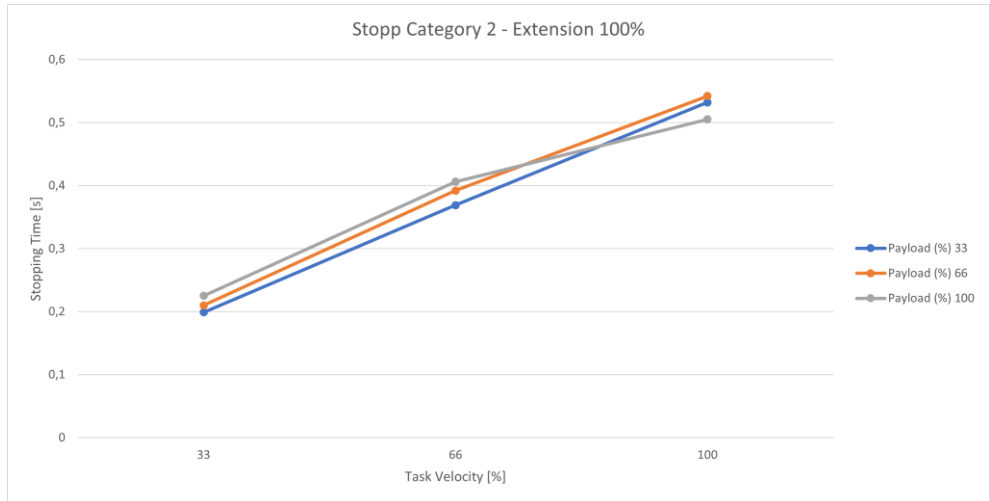


Fig. 14.41: Time; Stopp Cat 2; Joint 3 100%

14.3.4 Joint 4

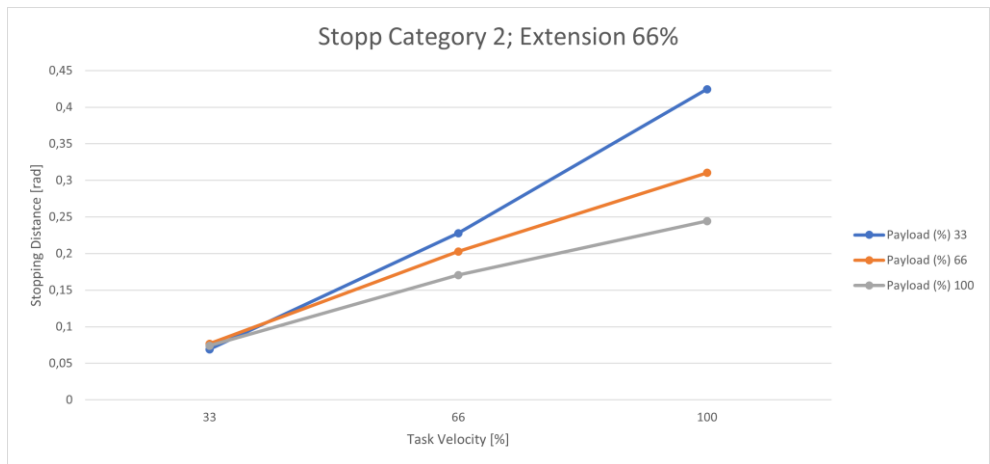


Fig. 14.42: Distance; Stopp Cat 2; Joint 4 66%

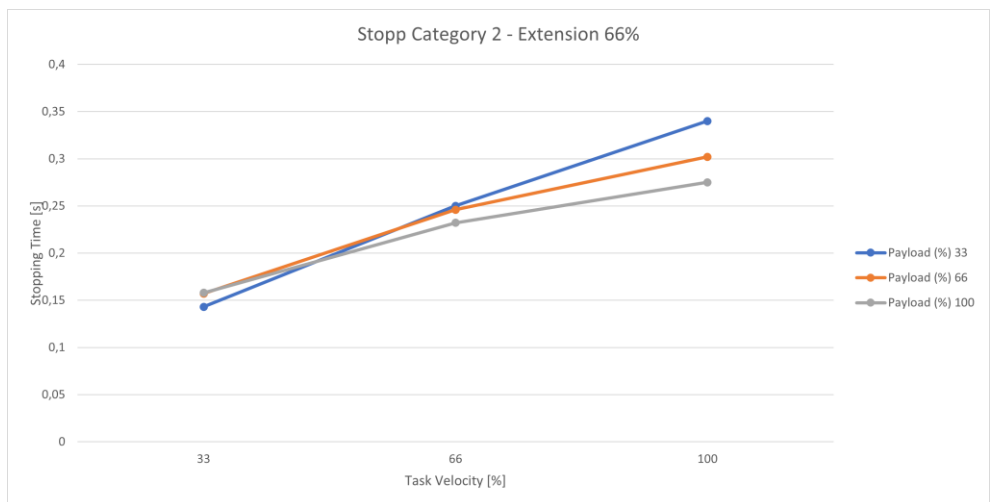


Fig. 14.43: Time; Stopp Cat 2; Joint 4 66%

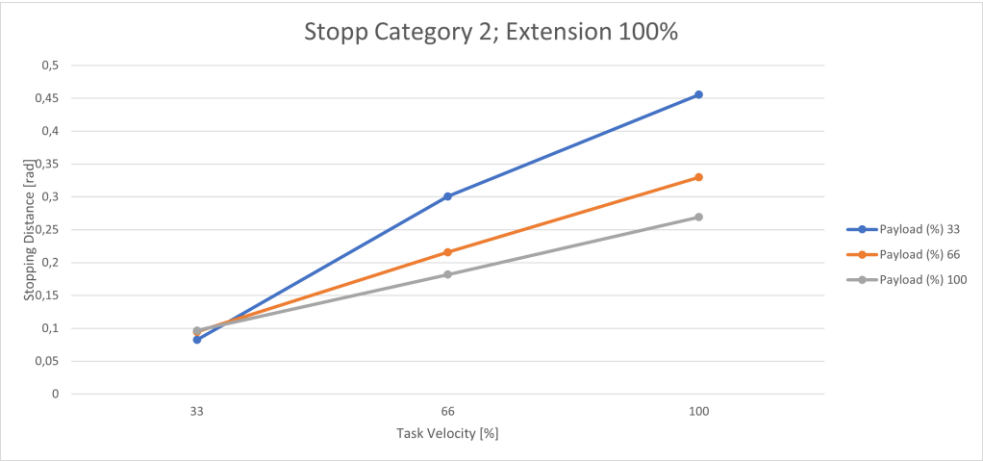


Fig. 14.44: Distance; Stopp Cat 2; Joint 4 100%



Fig. 14.45: Zeit; Stopp Cat 2; Joint 4 100%

14.4 Response times

Input	Reaction time motor	Reaction time Safe Output grip
User input	42 ms	48 ms
Safety function violation	36 ms	42 ms
	42 ms	48 ms
Central reaction – joint error	42 ms	48 ms
Central reaction – subordinate error	10 ms	–
Local reaction – joint error	10 ms	–
Local reaction – subordinate error	30 ms	36 ms
Reaction – Master Control error		

14.5 Safe Position Accuracy

If any safety function, that is based on safe position measurement is used, limited accuracy needs to be considered.

Main contributing factors to be considered include, but are not limited to structural elasticity, mechanical tolerances, assembly tolerances, signal processing, sensor accuracy.

The effect on the accuracy of these factors is strongly influenced by external forces, payload, dynamic state such as velocity and kinematic configuration (pose).

Safe position accuracy needs to be considered while parameterizing the safety functions to be used in an application. How much of a margin needs to be used is highly dependent on the aforementioned factors and therefore needs to be evaluated and validated individually.

15 GLOSSARY

Naming	Description
Administrator	<p>An administrator is a person authorized by the person responsible to access the robotic system and utilize the user interface Desk as follows:</p> <p>The administrator sets and changes roles, access rights, and passwords.</p> <p>The administrator sets and changes non-safety-related parameters of the system (e.g., when changing the end effector settings).</p> <p>The administrator programs and teaches the robot system.</p>
App	Apps are modular robot programs, each representing a partial step of a robot Task. They can be purchased in the Franka Store and be parametrized in Desk to form entire automation tasks.
Arm	The Arm is a tactile robotic arm with seven axes. It is part of Franka Research 3.
Axes	The Arm consists of seven consecutive axes. The motion is created in the axes.
Cartesian	The Cartesian space is the three-dimensional space where all axes (X, Y, and Z) are perpendicular.
Category 0 Stop	A Category 0 Stop is the stopping by immediate removal of power to the machine actuators (acc. to EN 60204:2019).
Category 1 Stop	A Category 1 Stop is a controlled stop with power available to the machine actuators to achieve the stop and then the removal of power when the stop is achieved (acc. to EN 60204:2019).
Category 2 Stop	A Category 2 Stop is a controlled stop with power available to the machine actuators (acc. to EN 60204:2019).
Center of Mass (CoM)	The center of mass is the center of gravity of an object. At this point, the gravitational force comes into effect.
Collaboration space	Space is accessible by both the operator and the robot during the execution of Tasks.
Connecting cable	The connecting cable connects the Arm with the Control.
Control	The Control is the main control unit and is part of Franka Research 3. The main control unit allows monitoring and control of the mechanical structure of the robot.
Desk	Desk is Franka Robotics web-based, intuitive, and graphical programming and user interface to exchange information and issue commands. It is part of the Franka UI.
Emergency Stop Device	<p>The Emergency Stop Device must be connected to the system to stop Franka Research 3 and execute a Cat. 1 Stop in case of an emergency. This will cause Franka Research 3 to brake at maximum capacity and the locking bolts to mechanically lock the Arm.</p> <p>The Emergency Stop Device is connected to the X3 connector at the Arm's base.</p>
Emergency Unlock Labels	The Emergency Unlock Labels can be found in three different areas on the Arm. They show where to use the Emergency Unlocking Tool to move the robot manually in case of an emergency.

GLOSSARY

Naming	Description
Emergency unlocking	Using the Emergency Unlocking Tool to unlock the fail-safe locking system to move the Arm manually is called emergency unlocking.
Emergency Unlocking Tool	The Emergency Unlocking Tool is a tool to manually unlock the fail-safe locking system in case of an emergency. Using the tool, the Arm can be moved even when the Arm has no power.
EMI Directive 2014/30/EU	The EMI Directive (2014/30/EC), referred to as EMI Directive, regulates the electromagnetic compatibility of devices within the European Economic Area, Switzerland, and Turkey.
Enabling Button	The Enabling Button is part of the Pilot-Grip and the Pilot. It allows the activation of the Arm's motion.
External Enabling Device	The External Enabling Device is connected to the X4 connector into the Arm's base. While half-pressing the External Enabling Device, it is possible to test and run automatic robot programs as long as the device remains half-pressed.
Fail-safe locking system	The fail-safe locking system locks all seven axes of the Arm. The Arm maintains its position, even when the power is turned off.
FCC rule 47 CFR part 15	FCC is the Federal Communications Commission. It is an independent US agency regulating radio, satellite, and cable communications. It regulates issues regarding the electromagnetic compatibility of devices.
Franka Robotics GmbH	Franka Robotics GmbH (abbreviated FR) is the company name. We have developed and now produce Franka Research 3.
Franka Research 3/ Franka Research 3 System	The Arm and Control system components compose the Franka Research 3 System, simply called Franka Research 3.
Franka Store	Franka Store is Franka Robotics online store, offering apps, bundles, and hardware for easy online ordering. It is part of Franka World: https://franka.world/
Franka UI	The Franka UI is the software framework for the web browser accessible user interface for Franka Research 3. It contains "Desk," "Watchman," and the "Settings" interface.
Franka World	Franka World is an online platform that interconnects customers, partners, and software and hardware developers, whose activity revolves around Franka Robotics' products and services. Franka World provides tools for managing Franka Research 3, access to an online store that features a continuously growing portfolio of software and hardware products, and the possibility of becoming part of an active and passionate community. Visit https://franka.world/ to make use of all the benefits.
Guiding / Hand-guiding	Guiding describes moving the robot by haptic interaction, e.g., to teach a new pose.
Guiding Button	The Guiding Button is located on the right of the Pilot-Grip. The Arm can be moved by pressing the Guiding Button and half-pressing the Enabling Button.

GLOSSARY

Naming	Description
Guiding Mode	Guiding modes ease guiding by locking or unlocking different directions or rotations in space, e.g., moving the Arm in three directions. One can switch between guiding modes using the Guiding-Mode Button on the Pilot-Grip or directly from Desk.
Hand/ Franka Hand	The Hand is an electrical two-finger parallel gripper and is optionally available. The Hand can be used for Franka Production3, Franka Research 3, and mountings according to the ISO flange design. The Hand is an end effector. It is not part of the certified machinery.
Integrator	<p>The integrator is responsible for assembling the partly completed machinery into the final machinery by combining the robot with other equipment or another machine, including additional robots to form a machine system.</p> <p>The integrator also conducts appropriate risk assessments to identify residual risks and to eliminate and minimize them according to ISO 12100.</p> <p>The integrator is responsible for the safety of the final application.</p>
Interaction	Franka Research 3 is designed to be easily programmed and operated, and to learn and relearn new Tasks quickly. When Franka Research 3 is in “monitored stop” mode or is being guided (teaching mode), Franka Research 3’s base is white to indicate that the Arm is ready for interaction.
Interface device	The interface device, a commercially available PC, tablet, or Notebook with a web browser, is connected to the Arm base via an Ethernet cable. The Franka UI can be accessed in a web browser via the interface device.
Low Voltage Directive 2014/35/EC	The Low Voltage Directive (2014/35/EC), hereinafter referred to as Low Voltage Directive (LVD), regulates the safety of electronically operated devices within the European Economic Area, Switzerland, and Turkey.
Machinery Directive (2006/42/EU)	The Machinery Directive (2006/42/EG), hereinafter referred to as Machinery Directive or MD, regulates a standardized level of protection to prevent accidents for machinery and partly completed machinery within the European Economic Area, Switzerland, and Turkey.
Maximum workspace	<p>Space that can be swept by the robot's moving parts plus the space that the end effector and the workpiece can sweep.</p> <p>For further information, see section Maximum-and safeguarded space in chapter Correct Installation.</p>
Operator	An operator is authorized to access Franka Research 3 and utilize the user interface Desk to use Franka Research 3 within the limits defined by the person responsible and the administrator. The operator is allowed to start, monitor, and stop the intended operation of Franka Research 3. Within Desk, the “operator” role can be assigned to users. Operators only have limited access to Desk.
Pilot	The Pilot is the user interface on the Arm for guiding and operating the Arm and/or Desk. It includes the Pilot-Grip and the Pilot-Disc.
Pilot-Disc	The Pilot-Disc is part of the Pilot and is used for interacting with the Arm and/or Desk.
Pilot-Grip	The Pilot-Grip is part of the Pilot and is used for manual guiding.
Pose	A pose is a combination of position and orientation in space.
Protective Measures	Protective measures achieve risk reduction according to 3.19 of ISO 12100. They are implemented and assessed by the following persons:

GLOSSARY

Naming	Description
	<p>Designer and/or Integrator (inherently safe design, safeguarding, and complementary protective measures, information for use)</p> <p>Responsible person/Integrator (organization: safe working procedures, supervision, permit-to-work systems; provision and use of additional safeguards; use of personal protective equipment; training)</p>
Responsible person	The responsible person is responsible for compliance with occupational health regulations and the operational safety ordinance. The person responsible for Franka Research 3 includes but is not limited to the entrepreneur, the director of the institute, the employer, or a delegate responsible for the use of Franka Research 3.
RoHS Directive 2011/65/EU	The RoHS Directive (2011/65/EU), hereinafter referred to as RoHS Directive, restricts the use of certain dangerous substances in electrical and electronic equipment within the European Economic Area, Switzerland, and Turkey.
Safeguarded space	<p>Perimeter safeguarding defines safeguarded space.</p> <p>For further information, see section Classification of spaces in chapter Correct Installation.</p>
Safety rules	A safety rule consists of a parameterized safety function, optional activation conditions, and a reaction to be carried out when the safety function is violated.
Safety scenarios	A set of safety rules defined in Watchman covering a specific risk situation, e.g., the "Test & Jog" scenario covering all rules for the Test & Jog mode.
Safety set-up	Definition of general safety-relevant settings such as safe input behavior or an end effector collision model.
SEEPO	Safety functions are "Safe End Effector Power Off." See chapter Safety functionalities for details.
Settings interface	The Settings interface is a web browser accessible user interface to set non-safety-related parameters of Franka Research 3, e.g., network settings, user roles, or passwords. It is part of the software framework Franka UI.
Single Point of Control (SPoC)	Single Point of Control (SPoC) is a functionality that allows only one user at a time to trigger critical actions, i.e., edit system settings and Tasks or trigger active robot actions like unlocking joints and running tasks.
SLD	Safety function "Safely limited distance." See chapter Safety functionalities for details.
SLP-C	Safety function "Safely limited Cartesian Position." See chapter Safety functionalities for details.
SLP-J	Safety function "Safely limited joint angle." See chapter Safety functionalities for details.
SLS-C	Safety function "Safely limited Cartesian speed." See chapter Safety functionalities for details.
SLS-J	Safety function "Safely limited joint speed." See chapter Safety functionalities for details.
SMSS	Safety function "Safely monitored standstill." See chapter Safety functionalities for details.
Stopping Distance	The stopping distance is the distance the Arm covers after a stop demand is received until the Arm comes to a full stop.

GLOSSARY

Naming	Description
Stopping Time	The stopping time is the time that passes after the stop demand is received, e.g., by the emergency device, until the Arm comes to a full stop.
Task	A Task in Desk represents an entire automation routine. A Task consists of one or several Apps.
Teaching	Teaching describes the process of parameterizing a Task and the contained Apps by manually guiding the robot or end effector. This includes, among others, the teaching of "poses" by guiding the robot to that poses.
Tracking Error	The actual motion of the Arm follows the target motion with a small deviation, a so-called tracking error.
Watchman	Watchman is a web browser accessible user interface to set and visually validate safety-related parameters of Franka Research 3, e.g., safety-related monitored speed or safety-related monitored spaces. It is part of the Franka UI.
Web browser	A software application runs on an interface device, serving as the Franka UI connection. A web browser provides the environment for Desk, Watchman, and the Settings interface. Examples include the following: Chrome, Edge, and Firefox.

16 INDEX

Additionally Equipment	41	Intended Use.....	16
Arm	12, 33, 41, 43, 52	Interaction	34
Manually Moving the Arm	23	Interface Device.....	32, 42
Unpacking the Equipment.....	44	Labeling on the Equipment	12
Cartesian.....	26	Functional earth label	13
Center of Mass (CoM).....	18, 77	Grasping position label	13
Certificates	10	Low Voltage Directive 2014/35/EC.....	110
Cleaning	82	Machinery Directive 2006/42/EC	16
Connecting cable.....	44, 61	Maintenance	81
Control	13, 17, 32, 41, 43, 53	Maximum space	48, 66, 71, 87
Desk.....	34, 62	Misuse	16, 19
Disposal.....	11, 81, 82, 83	Operator	26
Emergency.....	14, 17	Pilot.....	33, 34
Emergency Stop Device	14	Pilot-Disc.....	34
Emergency stop installation.....	22	Pilot-Grip.....	34
Emergency unlock label.....	12	Pilot-Mode	34
Emergency Unlocking	12	Ports	57
Emergency Unlocking Tool.....	23	C2.....	59
EMI Directive 2014/30/EU	109	X1.....	37
Enabling Button.....	74	X2.....	57
End effector	39, 65, 84	X3.....	58
External Enabling Device.....	14, 41, 58, 74	X4.....	58
Fail-safe Locking System	23	X5.....	59
FCC rule 47 CFR part 15	109	X6.....	59
Franka.....	7, 109	Practical Tips for Usage and Positioning.....	67
Franka Store.....	109	Protective Measures	17
Franka UI	7, 20	Responsible person.....	7
Gelenk-Referenz-Position des Armes.....	69	Restarting	75
Glossary	108	RoHS Directive 2011/65/EU	10
Guiding		Safety	15
Guiding Button.....	36	Installation of Safety Peripherals.....	22
Guiding-Mode Button.....	36	Safe inputs.....	26
Hand	34, 39, 42	Safe outputs.....	28
Handling.....	84, 85, 86	Safeguarded space	48
Installation.....	17, 40, 42, 43	Safety Functionalities	25
Preparing the Installation Site	51	SEEPO	27
Wiring and Electrical Installation	56	Self-test.....	73
Integrator.....	19	Single Point of Control (SPoC)	20

INDEX

Site of installation	49, 50	Category 2 stop.....	28
SLD.....	28	Stopping time.....	22
SLP-J	27	Stopping time and distances.....	89, 107
SLS-C	27	Switching Off.....	75
SLS-J	28	Switching On.....	71
SMSS.....	27	Task.....	74
Stopping distance.....	22	Transport	71
Stopping functions.....	28	Unpacking	44
Category 0 stop.....	28	User interface	34
Category 1 stop.....	28		

17 TABLE OF FIGURES

Fig. 3.1: Type label.....	12	Fig. 10.11: Lift out control	47
Fig. 3.2: Emergency unlock label.....	12	Fig. 10.12: Classification of spaces.....	48
Fig. 3.3: Hot Surface label	13	Fig. 10.13: Preparing the baseplate.....	52
Fig. 3.4: Functional earth label	13	Fig. 10.14: Drilling template	53
Fig. 3.5: Lifting position label.....	13	Fig. 10.15: Mounting of the arm.....	55
Fig. 3.6: Type label control.....	13	Fig. 10.16: Overview connection diagram.....	57
Fig. 3.7: Type label external device	14	Fig. 10.17: X2 interface	58
Fig. 3.8: Type label emergency stop device.....	14	Fig. 10.18: X3 - Safe Inputs.....	58
Fig. 4.1: Connection Emergency Stop Device.....	22	Fig. 10.19: X6 interface	59
Fig. 4.2: Emergency unlock	24	Fig. 10.20: Connection with the functional earth	60
Fig. 4.3: Pushing away manually	24	Fig. 10.21: Connection of connecting cable the arm	61
Fig. 6.1: Equipment overview	32	Fig. 10.22: connection of connecting cable to the	62
Fig. 6.2: Overview Arm.....	33	Fig. 10.23: Connection of external Enabling Device	62
Fig 6.3: Pilot	34	Fig. 10.24: Connection of operating device	63
Fig. 6.4: switching status Enabling Button.....	35	Fig. 10.25: Connection of the protective device	65
Fig. 6.5: Enabling Button	35	(here emergency stop command device).....	65
Fig. 6.6: Guiding Mode Button.....	36	Fig. 10.26: Measuring points ESD measurement	67
Fig. 6.7: Guiding Mode Button.....	36	Fig. 10.27: Design of the workspace.....	68
Fig. 6.8: Guiding Button.....	36	Fig. 10.28: Joint reference position of the arm ...	69
Fig. 6.9: Guiding Button.....	37	Fig. 10.29: Distance to arm bumping the head....	70
Fig. 6.10: Connection ports at robot base	37	Fig. 10.30: Distance to arm bumping.....	70
Fig. 6.11: End effector flange	39	Fig. 10.31: Distance to arm Squeezing the hand.	70
Fig. 6.12: Dimensions and connection ports of	39	Fig. 10.32: Protective equipment does not wear	71
control	39	jewellery.....	71
Fig. 6.13	40	Fig. 11.1: Switching on Control.....	72
Fig. 7.1: scope of delivery Arm.....	40	Fig. 11.2: Blue status lights on Arm.....	72
Fig. 7.2: scope of delivery Control.....	41	Figure 65: Lifting positions.....	85
Fig. 7.3: Scope of delivery devices	41	Figure 66: Lifting the Arm	87
Fig. 7.4: Scope of delivery Connecting cable	41	Figure 67: Packing the Arm.....	87
Fig. 10.1: Overview of interfaces	43	Figure 68: Closing the box	88
Fig. 10.2: Packing	44	Fig. 17.1: Illustration of extension states	89
Fig. 10.3: Unpack main carton.....	45	Fig. 17.2: Distance; Stopp Cat 1; Joint 1, 33%....	90
Fig. 10.4: Removal of individual cartons	45	Fig. 17.3: Time; Stopp Cat 1; Joint 1 33%.....	90
Fig. 10.5: Inner cartons.....	45	Fig. 17.4: Distance; Stopp Cat 1; Joint 1 66%....	91
Fig. 10.6: Open box from Arm.....	46	Fig. 17.5: Time; Stopp Cat 1; Joint 1 66%.....	91
Fig. 10.7: Unpack arm	46	Fig. 17.6: Distance; Stopp Cat 1; Joint 1 100% ...	91
Fig. 10.8: Lift out arm.....	46	Fig. 17.7: Time; Stopp Cat 1; Joint 1 100%	92
Fig. 10.9: Open box of control	47		
Fig. 10.10: Removal packaging.....	47		

TABLE OF FIGURES

Fig. 17.8: Distance; Stopp Cat 1; Joint 2 33%.....	92	Fig. 17.36: Distance; Stopp Cat 2; Joint 3 41% .	102
Fig. 17.9: Time; Stopp Cat 1; Joint 2 33%.....	92	Fig. 17.37: Time; Stopp Cat 2; Joint 3 41%	102
Fig. 17.10: Distance; Stopp Cat 1; Joint 2 66% ...	93	Fig. 17.38: Distance; Stopp Cat 2; Joint 3 66% .	103
Fig. 17.11: Time; Stopp Cat 1; Joint 2 66%	93	Fig. 17.39: Time; Stopp Cat 2; Joint 3 66%	103
Fig. 17.12: Distance; Stopp Cat 1; Joint 2 100%.	93	Fig. 17.40: Distance; Stopp Cat 2; Joint 3 100%	103
Fig. 17.13: Time; Stopp Cat 1; Joint 2 100%.....	94	Fig. 17.41: Time; Stopp Cat 2; Joint 3 100%.....	104
Fig. 17.14: Distance; Stopp Cat 1; Joint 3 33% ...	94	Fig. 17.42: Distance; Stopp Cat 2; Joint 4 66% .	104
Fig. 17.15: Time; Stopp Cat 1; Joint 3 33%	94	Fig. 17.43: Time; Stopp Cat 2; Joint 4 66%	104
Fig. 17.16: Distance; Stopp Cat 1; Joint 3 66% ...	95	Fig. 17.44: Distance; Stopp Cat 2; Joint 4 100%	105
Fig. 17.17: Time; Stopp Cat 1; Joint 3 66%	95	Fig. 17.45: Zeit; Stopp Cat 2; Joint 4 100%.....	105
Fig. 17.18: Distance; Stopp Cat 1; Joint 3 100%.	95		
Fig. 17.19: Time; Stopp Cat 1; Joint 3 100%.....	96		
Fig. 17.20: Distance; Stopp Cat 1; Joint 4 66% ...	96		
Fig. 17.21: Time; Stopp Cat 1; Joint 4 66%	96		
Fig. 17.22: Distance; Stopp Cat 1; Joint 4 100%.	97		
Fig. 17.23: Time; Stopp Cat 1; Joint 4 100%.....	97		
Fig. 17.24: Distance; Stopp Cat 2; Joint 1 33% ...	98		
Fig. 17.25: Time; Stopp Cat 2; Joint 1 33%	98		
Fig. 17.26: Distance; Stopp Cat 2; Joint 1 66% ...	99		
Fig. 17.27: Time; Stopp Cat 2; Joint 1 66%	99		
Fig. 17.28: Distance; Stopp Cat 2; Joint 1 100%.	99		
Fig. 17.29: Time; Stopp Cat 2; Joint 1 100%.....	100		
Fig. 17.30: Distance; Stopp Cat 2; Joint 2 33%.	100		
Fig. 17.31: Time; Stopp Cat 2; Joint 2 33%	100		
Fig. 17.32: Distance; Stopp Cat 2; Joint 2 66%.	101		
Fig. 17.33: Time; Stopp Cat 2; Joint 2 66%	101		
Fig. 17.34: Distance; Stopp Cat 2; Joint 2 100%	101		
Fig. 17.35: Time; Stopp Cat 2; Joint 2 100%.....	102		

Franka Robotics GmbH
Frei-Otto-Strasse 20
80797 Munich
Germany